GEBRAUCHSANLEITUNG
Originalversion
OPERATING MANUAL
MODE D'EMPLOI
MANUAL DE INSTRUCCIONES

TW 7200

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Operating Manual Page 25 ... 46

Important notes:

The operating manual is part of the product. Before initial operation, please carefully read and observe the operating manual and keep it. For safety reasons the product may only be used for the purposes described in these present operating manual. Please also consider the operating manuals for the devices to be connected.

All specifications in this operating manual are guidance values which are valid at the time of printing. However, for technical or commercial reasons or in the necessity to comply with the statutory stipulations of various countries, the manufacturer may perform additions to the product without changing the described properties. A potentially more recent version of this manual is available on our internet website. The German version is the original version and binding in all specifications!

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TABLE OF CONTENT

1	Technical Specifications of the titration sample changer TW 7200	27
1.1	Notes to the operating manual	27
1.2	Intended Use	27
1.3	Technical Specifications	28
1.4	Warning and safety information	31
1.4.1	Warning signs	31
1.4.2	Chemical and biological safety	32
1.4.3	Flammable liquids	32
2	Installation and Commissioning	33
2.1	Operating environment	33
2.2	Unpacking and setting up	33
2.3	Back panel of the titration sample changer TW 7200	34
2.4	System preparations	35
2.4.1	Installation of the hose bracket	35
2.4.2	Installation of the titration head	35
2.4.3	Adjusting the lifting height	35
2.4.4	Connecting a titration system or to a PC	35
2.4.5	Electrodes, titration tips, stirrers	35
2.4.6	Cleaning	35
3	Working with the titration sample changer TW 7200	36
3.1	Sample tray	36
3.1.1	General	36
3.1.2	Versions	37
3.1.3	Setting up, filling and removing the sample tray	37
3.2	Built-in magnetic stirrer and external rod stirrer	38
3.3	I/O connection of the sample changer TW 7200	39
3.3.1	General I/O description	39
3.3.2	Electrical properties of the I/O	39
3.3.3	Inputs of the I/O connector	39
3.3.4	Outputs of the I/O connection	39
3.3.5	The "Pump 1" and "Pump 2" connections	39
4	Connection to a titrator or a computer titration system	40
4.1	Interface connections	40
4.2	Installation in an Daisy Chain	40
4.3	Data transmission parameters	40
5	Communication with the TW 7200	41
5.1	RS command list of the sample changer TW 7200	41
5.2	Network communication	43
5.2.1	Communication via UDP	43
5.2.2	Communication via TCP	44
5.3	Extended RS command list - network	44
6	Operating display and error messages	45
6.1	Operating display	45
6.2	Errors and messages	45
6.3	Firmware update	46
7	Guarantee	46
8	Storage and transportation	46
9	Recycling and Disposal	46
10	EC – Declaration of Conformity	46

Copyright

© 2021, Xylem Analytics Germany GmbH

Reprinting - even as excerpts - is only allowed with the explicit written authorization.
Germany. Printed in Germany.

1 Technical Specifications of the titration sample changer TW 7200

1.1 Notes to the operating manual

The provided operating manual will allow you the proper and safe handling of the product. For maximum security, observe the safety and warning instructions in the operating manual!

-  Warning of a general danger:
Non-compliance results (can result) in injury or material damage.
-  Important information for device use.
-  Refers to another part of the operating manual.

The menu screens shown in this operating manual serve as an example and may differ from what you see!

1.2 Intended Use

The titration sample changer TW 7200 allows serial titrations with automatic sample change. It can be used for many samples in conjunction with SI Analytics® titration instruments.

The examples of possible use include:

- Acid and base determination in aqueous solutions such as p and m value,
- Acid and base determination in non-aqueous solutions,
- Acid numbers in oils
- Titration of strong and weak acids and bases
Redox titrations such as iodometry, manganometry, chromatometry, and COD determinations,
other mV titrations, e.g. chloride

All these titrations are performed in an open beaker or comparable vessel, which is placed in the sample tray.

With a titration head specially designed for this purpose, COD titrations can be carried out directly in the standardized reaction vessels according to DIN 38 409, Part 41.

The TW 7200 can be controlled externally via various interfaces (RS, USB or LAN) from a PC or a titrator.

The TW 7200 has an easily replaceable round sample tray. There are two types of sample trays:

In one version, all samples are only in one outer circle on the sample tray. In the other version, the sample tray has two circles, in each of which samples can be placed. By means of a horizontally movable axis, it is possible to operate the inner circle of the two-row sample tray. The TW 7200 has 2 stirrers, one each at the measuring position of the outer and inner circuit.

There are 11 different sample trays, each with a different number of positions, for different sample vessels. Each position is numbered and can be targeted by the control software.

 The TW 7200 is not intended for use with potentially biohazardous substances.

 **Do not use the device in hazardous locations!**

 **General:**

The safety guidelines that are applicable to the handling of chemicals have to be observed under all circumstances. This applies in particular to inflammable and/or etching liquids.

1.3 Technical Specifications

Translation of the legally binding German version

(Release: 4. May 2021)

EMC compatibility according to the Council Directive: 2014/30/EU;
 applied harmonized standards: EN 61326-1
 Low-voltage directive according to the Council Directive 2014/35/EU;
 testing basis EN 61 010-1: for laboratory equipment
 RoHS Council Directive 2011/65/EU
 FCC Part 15B and ICES 003

Country of origin: Germany, Made in Germany

Automatic Titration Sample Changer TW 7200 with exchangeable sample tray.

Sample trays are available for 12, 16, 18, 24, 25, 28, 30, 38 and 48 beakers or for 24 COD reaction vessels according to DIN 38 409, part 41.

Lifting height: 270 mm, limited to 130 mm at the factory on delivery

Beakers for sample trays with 16/25 samples	Order No. / TZ No.
150 ml with spout, low form	285216747 / TZ 1784
150 ml with spout, low form, Super Duty	285202008 / TZ 1788
250 ml with spout, high form	285201997 / TZ 1787
250 ml with spout, low form, Super Duty	285226140 / TZ 1780
Beakers for sample trays with 18/25 samples	Order No. / TZ No.
100 ml without spout, high form	285202030 / TZ 1765
Beakers for sample trays with 24/38 samples	Order No. / TZ No.
50 ml without spout, high form	1064720 / TZ 1783
50 ml, of plastic	285226150 / TZ 3973
75 ml without spout	285216756 / TZ 1786
Beakers for sample trays with 12 samples	Order No. / TZ No.
250 ml with spout, low form	1064589 / TZ 1785
600 ml without spout, high form	285226160 / TZ 1766
Sample bottles for sample trays with 16, 30 and 48 positions for sample preparation viscometry	Order No. / VZ No.
40 ml sample bottles with ND 24	285420430 / VZ 7088
100 ml laboratory bottle DURAN®	285420869 / VZ 7081

Power supply: by external multi-range power supply from 100 – 240 V, 50/60 Hz
 Input voltage: 12 Volt DC, 2500 mA
 Power consumption 30 W
 Corresponds to protection class III:
 Protection class for dust and humidity IP 50 according to DIN 40 050

 **Only use the power supply TZ 1853 or a power supply approved by the manufacturer!**

RS-232-C Interface:

separated galvanically through photocoupler, Daisy Chain function available

Data bits: static, **8** Bit
 Stop bit: adjustable, **1** or 2 Bit (default: 1 Bit)
 Parity: adjustable: even / odd / **none**
 Baud rate: adjustable: **4800**, 9600, 19200 (Default 4800 Baud)
 Address: adjustable, (0 to 15; default: 3)

RS-232-1 for connection of titrator, PC, burette,
 input Daisy Chain

RS-232-2 for connection of balances of the types Mettler, Sartorius, Kern, Ohaus, (for more, please
 contact us),
 Exit Daisy-Chain

USB Interface:

USB-type A to connect USB stick for software update
 USB-type B Mini-B for PC connection (virtual COM port)

Ethernet Interface: for connecting a local network (LAN)

Pump: 2 potential-free extra-low voltage switching outputs for controlling pumps and valves

Maximum cable length:

3 meters per cable

Stirrer: External connection:
 0V to 3.3V DC out, 400 mA
 Power supply for rod stirrer

Material:

Device: Housing; powder coated aluminum, stainless steel
 Trays: Stainless steel, powder coated

Dimensions:

Device: 14.3 x 62 x 50.5 cm (W x H x D)
 Trays: Diameter 45 cm, H = 6.5 cm (for beakers)
 H = 11 cm (for COD reaction vessels according to DIN 38 409, Part 41)

Weight:

Device: approx. 10.3 kg
 Trays: with 12 beakers 250 ml: approx. 2.7 kg
 with 16 beakers 150 ml: approx. 2.6 kg
 with 24 beakers 50 ml: approx. 2.4 kg
 with 24 COD reaction vessels: approx. 6.3 kg

Connections:

- Stirrer: Control output on the back of the sample changer;
Plug connection: 3 pole DIN socket
- Power supply: Low voltage input 12 V / – on the back of the sample changer;
Plug connection: Socket for low-voltage connection -jack plug-,
positive pole on pin contact, inner contact $\varnothing = 2.1$ mm, USA/Japan
- USB: 1 x USB-type- A and 1 x USB-type-Mini B
- Serial interfaces: 2 x Mini-DIN 4 pin socket
- Ethernet Interface: 1 x RJ 45 jack
- I/O Port: 1 x 9 pin D-Sub-socket
- Pump connections: 2 x 2 pole Hirschmann STASEI 2 connectors

Operation display: 2colored LED display (red/green).
Operating states, error states and operating faults are signaled by color signals
and flashing codes

Ambient conditions:

Do not use the device in hazardous locations!

- Climate: Ambient temperature: + 10 ... + 40 °C for operation and storage
Humidity according to EN 61 010, Part 1:
Max. relative humidity 80 % for temperatures up to 31 °C,
linear decrease down to 50 % relative humidity at a temperature of 40 °C
- Altitude: Device: No restrictions
Power supply: up to 5000 m
- Pollution degree: Pollution degree IP 20, indoor use only.

1.4 Warning and safety information

The device corresponds to protection class III.

It was manufactured and tested according to EN 61 010 - 1, Part 1 „**Protective Measures for electronic measurement devices**“and control device and has left the factory in an impeccable condition as concerns safety technology. In order to maintain this condition and to ensure safe operation, the user should observe the notes and warning information contained in the present operating instructions. Development and production is done within a system which meets the requirements laid down in the Norm DIN EN ISO 9001 standard.

⚠ For reasons of safety, the device has only be used for the range of application described in the present operating manual. Nonobservance of the intended proper use of the device may result in personal injury or damage to property.

⚠ For reasons of safety, the device and the power supply must be opened by authorised persons only; this means, for instance, that work on electrical equipment must only be performed by qualified specialists. **In case of nonobservance of these provisions the device and the power supply may constitute a danger: electrical accidents of persons or fire hazard!** Moreover, in the case of unauthorised intervention in the device or the power supply, as well as in the case of negligently or deliberately caused damage, the warranty will become void.

⚠ Prior to switching the device on it has to be ensured that the operating voltage matches the mains voltage. The operating voltage is indicated on the specification plate (underside of the device and backside of the power supply). **Nonobservance of this provision may result in damage to the device and the power supply, or in personal injury or damage to property!**

If the plate rotates, the titration head moves or the horizontal axis moves, the operating display will light up red. In this state, the sample changer must not be touched and no sample vessels must be removed or adjusted.

⚠ **If it has to be assumed that safe operation is impossible, the device has to be put out of operation and secured against inadvertent putting to operation.** In this case please switch the device off, pull plug of the mains cable out of the power supply, and remove the device from the place of work.

Examples for the assumption that a safe operation is no longer possible,

- if the package is damaged,
- if the device shows visible damages,
- if the power supply shows visible damages,
- if the device does not function properly,
- if liquid has penetrated into the casing,
- if the device has been altered technologically or if unauthorized personnel tried or succeeded to open the device as attempt to repair it.

In case that the user operates such a device, all thereof resulting risks are on the user!

⚠ The device has not be stored or operated in humid rooms.

1.4.1 Warning signs

Observe all labels and signs attached to the devices. Non-observance may result in personal injury or damage to the devices!

Danger of hand injuries!

1.4.2 Chemical and biological safety

 The TW 7200 is not intended for use with potentially biohazardous substances.

 **The relevant regulations regarding the handling of the substances used have to be observed:** The Decree on Hazardous Matters, the Chemicals Act, and the rules and information of the chemicals trade. On the part of the user it has to be ensured that the persons entrusted with the use of the unit are experts in the handling of substances used in the devices or that they are supervised by specialized persons, respectively.

 When using biohazardous substances, the regulations for handling the substances used must be observed. In such cases, the use is the sole responsibility of the user.

 For all work with chemicals: **Always wear protective glasses!** Please observe the memorandums of the employer's liability insurance associations and the safety data sheets of the manufacturers.

 Dispose of all used solutions in accordance with national regulations and laws. Select the type of protective equipment according to the concentration and quantity of the hazardous substance at the respective workplace.

1.4.3 Flammable liquids

When handling flammable liquids, make sure that there is no naked flame in the vicinity of the equipment. Adequate ventilation must be provided. Only small quantities of flammable liquids should be kept in the workplace.

 When working with liquids that do not correspond to common reagents, particular attention must be paid to the chemical resistance of the materials of the devices (vgl.  1.3 Technical Specifications).

2 Installation and Commissioning

2.1 Operating environment

i Observe the following points to ensure that the device functions properly and has a long service life

- The ambient temperature must be 10 to 40 °C.
- The relative humidity must be below 80 % (without condensation).
- The titration head of the sample changer must be able to move freely.
- Do not use or store the device in dusty, damp or wet places.
- Make sure that no liquids can get into the device and wipe off any liquids on the device immediately.
- Protect the device from vibrations, direct sunlight, corrosive gases and strong magnetic and/or electric fields.
- Use only original spare parts and accessories.

2.2 Unpacking and setting up

The device has been individually assembled for you, therefore there may be deviations with regard to the described scope of delivery and the accessories. Please refer to the enclosed packing list for the exact scope of delivery. If you have any questions, please contact us directly (for the service address, see backside of this operating manual).

The device and all accessories as well as the peripheral devices have been carefully checked at the factory for function and dimensional accuracy. Please ensure that even the small additional parts are removed from the packaging without leaving any residue.

The device can be set up on any flat surface. Do not push any objects under the device.

2.3 Back panel of the titration sample changer TW 7200

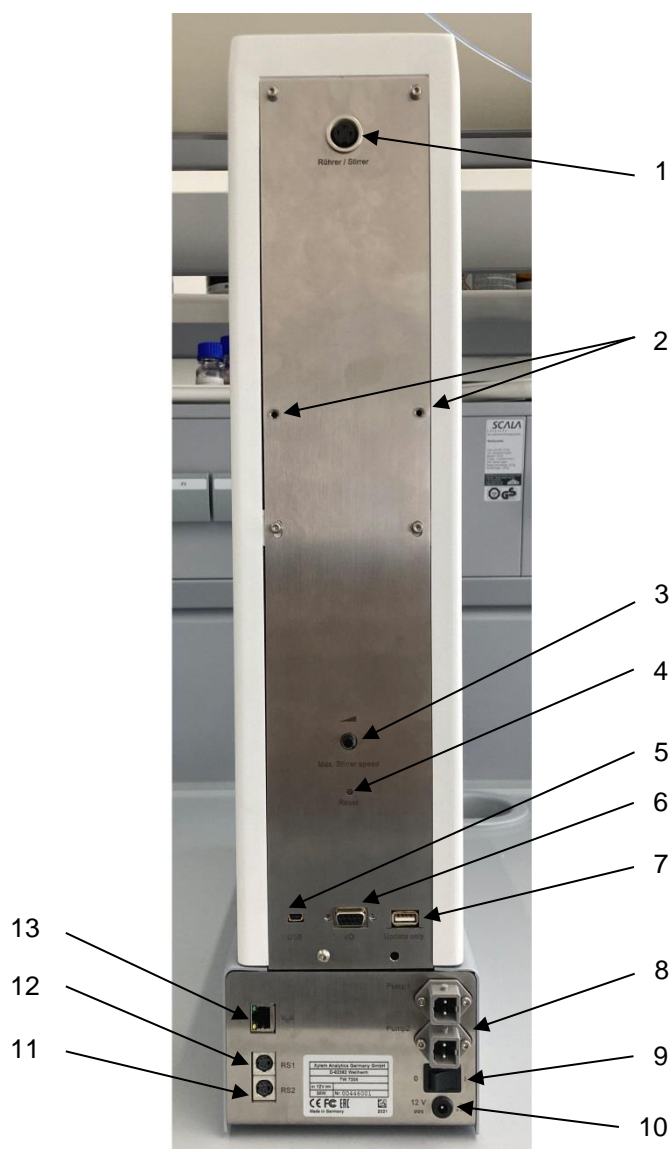
Fig. 1

The TW 7200 is equipped with the following connections (Fig. 1):

- 1) Connection for stirrer (3 pole DIN socket)
- 2) Fixing points for hose and cable holders
- 3) Setting the stirring speed of the built-in magnetic stirrers
- 4) Reset hole, for resetting the RS parameters
- 5) USB mini-B socket for connection to a PC
- 6) Connection for pump and valve modules, I/O ports, 9-pin D-submini female connector
- 7) USB-A socket for connecting a USB stick for software update
- 8) Two pump connections
- 9) Power switch
- 10) Power supply

Two RS-232 ports (Mini-DIN):

- 11) RS-2 port for data communication with other peripheral devices of a device chain
- 12) RS-1 port for communication with PC and data communication with titrators
- 13) Ethernet interface, for connecting a local network (LAN)

2.4 System preparations

2.4.1 Installation of the hose bracket

The hose holder is attached to the mounting holes provided for this purpose on the back of the TW 7200 using the knurled screws supplied.

When subsequently attaching the hoses and electrode cables, ensure that they are not kinked during the entire movement sequence.

2.4.2 Installation of the titration head

The sample changer is delivered from the factory without titration head.

When mounting or removing the titration head, make sure that the horizontal arm is retracted.

2.4.3 Adjusting the lifting height

The sample changer is factory-set to a stroke height of approx. 130 mm (half travel).

It is possible to change the stroke height by software commands.

2.4.4 Connecting a titration system or to a PC

The TW 7200 sample changer is connected to a titration system via the TZ 3094 data cable or directly to a PC via the TZ 3097 data cable or directly to a PC via the. For this purpose, the upper of the two 4-pole mini DIN sockets (data communication RS-232-C interface) can be used. In addition, a USB mini-B socket or the LAN connection are also available.

If the connection to a PC is made via the supplied USB cable (Type A (M) --- USB Type Mini-B (M), 2.0 m), a driver is required. This driver is located on the USB stick of the TitriSoft. If the PC is connected to the Internet, the driver will be installed automatically. Communication via the USB-RS interface is possible immediately. For further information see  section 4.

2.4.5 Electrodes, titration tips, stirrers

The required electrodes and titration tips are inserted into the corresponding bores (NS 14.5) of the titration head. A possibly used rod stirrer (TZ 1846 or TZ 1847) is also inserted into one of the bores. The electrical connection of the stirrer is established with a cable attached to the titration head to the socket "Stirrer" on the rear of the TW 7200.

The stirring speed can be adjusted by software commands. With the potentiometer on the back of the TW 7200 can be used to limit the maximum possible speed of the built-in magnetic stirrers. If the maximum possible stirring speed is to be adjustable by RS command, the potentiometer must be set to maximum level.

2.4.6 Cleaning

The TW 7200 is a maintenance-free device. For safe and precise operation, it is necessary to keep the device clean. A soft cloth moistened with water or a mild cleaning agent should be used for (external) cleaning.

 **Make sure that no liquid enters the device during cleaning!** If a titration or rinsing vessel overflows, or if small amounts of liquid spill over the TW 7200 sample changer, it must be switched off (then disconnect the power supply from the mains) and cleaned (dried). If larger quantities of liquid enter the device, it must be ensured that the device is completely dry and safety-checked before it is switched on again.

Decontamination

If the sample changer has been contaminated with substances that require special decontamination, it must be decontaminated in accordance with the relevant regulations.

Plug the low-voltage cable of the TZ 1853 power supply unit into the upper 12 V socket "in". Only then plug the power supply unit into the mains socket.

3 Working with the titration sample changer TW 7200

3.1 Sample tray

3.1.1 General

There are two types of sample trays for the TW 7200:

In the first version (Fig. 2), all samples are located only in an outer circle on the sample tray. At the measuring position there is a cup sensor, which detects in the outer circle whether a sample cup is set. The sample changer can be programmed so that the titration head only lowers when a sample vessel is at the titration position.

In the second version (Fig. 3), the sample tray has two circles into each of which samples can be placed. By means of a horizontally movable axis, it is possible to operate the inner circle of the double-row sample tray.

i The device automatically detects which type of sample tray is used.

i If a sample tray with 2 circles is used, it is not monitored whether a sample vessel is located at the measuring position. In this case, the operator must ensure that a sample vessel is always positioned at the measuring position.

Fig. 2

Fig. 3

3.1.2 Versions

Description	Ord. No.	TZ No.
for 12 sample vials, incl. 20 beakers 250 ml, low form	285228080	TZ 4012
for 16 sample vials without beakers. Suitable beakers are 150 ml TZ 1784, TZ 1788 or 250 ml TZ 1787, TZ 1780	285228090	TZ 4016
for 18 sample vials, without beakers. Suitable beakers are TZ 1765 (100 ml high form without spout)	285228100	TZ 4018
for 24 sample vials, without beakers. Compatible beakers/sample vials are 50 ml TZ 1783 / 75 ml TZ 1786	285228110	TZ 4024
double row for 25 sample vials, without sample vials. Suitable beakers are 150 ml TZ 1784, TZ 1788 or 250 ml TZ 1787, TZ 1780	285228120	TZ 4025
double row for 28 sample vials, without sample vials. Suitable beakers for this are TZ 1765 (100 ml high form without spout)	285228130	TZ 4028
for 30 sample vials, without sample vials. Suitable for 40 ml sample bottles with ND 24 VZ 7088	285228140	TZ 4030
double row for 38 sample vials, without beakers. Suitable beakers/sample vials are 50 ml TZ 1783 / 75 ml TZ 1786	285228150	TZ 4038
for 24 CSB vessels according to DIN 38 409, without CSB vessels	285228160	TZ 4040
double row for 48 sample vials, without sample vials. Suitable for 40 ml sample bottles with ND 24 VZ 7088	285228170	TZ 4050
for 12 sample vials, incl. 20 beakers 600 ml, high form	285228180	TZ 4053
for 16 sample vials, without sample bottles. Suitable for 100 ml laboratory bottles VZ 7081	285228190	TZ 4058

3.1.3 Setting up, filling and removing the sample tray

The TW 7200 automatically recognizes the sample tray size used by means of a code on the underside of the sample tray.

The sample trays are filled clockwise according to the numbers applied to the sample tray.

The connection to the sample changer is made by simply placing it on the drive cone provided. The sample tray must be turned until it locks into place.

The sample tray can be removed again by simply lifting it up.

3.2 Built-in magnetic stirrer and external rod stirrer

The TW 7200 has two built-in magnetic stirrers and the possibility of stirring directly in the solution from above by means of a connected rod stirrer (Fig. 4). In conjunction with the TitrISOFT PC software, these stirrers can be switched on and off via the interfaces.

Fig. 4

i When the sample tray is turned and the head is raised, the magnetic stirrers and, if used, the rod stirrer are switched off.

When the titration is started, the stirrer starts at the speed set in the method. When setting the stirring time, the time should not be less than 10 seconds.

When setting the stirring speed, make sure that the setting knob is set to full stirring speed, otherwise the maximum possible stirring speed is limited according to the level set on the setting knob. Fine adjustment of the desired stirring speed is then possible using the rotary knob on the backside of the TW 7200.

The speed of the magnetic stirrer can be preset and read out in revolutions/minute.

3.3 I/O connection of the sample changer TW 7200

3.3.1 General I/O description

A 9-pin plug connection is available on the TW 7200 sample changer for connecting an I/O box (Fig. 1, pos. 6). The devices available for this purpose e.g. the MP 25 pump or a peristaltic pump) can be connected via this I/O connection. The I/O connection is controlled via the TitrISOFT PC software. There are 4 outputs and 1 input available. The necessary commands can be taken from the command list (see section 5).

3.3.2 Electrical properties of the I/O

The inputs and outputs of the I/O connection are galvanically isolated from the titration sample changer electronics by means of TCMT 1600 optocouplers. No power consumers can be switched directly. The pump and valve units available from SI Analytics® can be connected directly to the I/O unit via a cable.

3.3.3 Inputs of the I/O connector

When using the input, the input must be connected to ground potential. The TCMT-1600 optocoupler has been used.

The pin assignment is shown as follows:

Input	Pin number
1	5
Common mass	6, 7, 8 and 9

The state of the input can be queried with the command "aaIP".

The response is: "aaIP0" or "aaIP1" depending on the state of the input.

3.3.4 Outputs of the I/O connection

The outputs of the I/O connector are assigned as follows:

Output	Pin number
1	1
2	2
3	3
4	4
Common mass	6, 7, 8 and 9

When using an output, the respective output is pulled to ground by the sample changer. Maximum current load of the TCMT 1600 optocoupler: 50 mA.

3.3.5 The "Pump 1" and "Pump 2" connections

These connections, located on the rear of the instrument, are used to control a connected pump (e.g. MP 25) or valve. The built-in relay can switch a voltage of 24 V and a current of 2 A. By means of 2 different commands, the two connections can be switched on and off. The connections can also be switched on time-controlled via a command.

4 Connection to a titrator or a computer titration system

4.1 Interface connections

For data communication with other devices, the sample changer TW 7200 has two serial interfaces (RS-232-C), a USB mini-B socket and a LAN connection. The connectors for these interfaces are located on the rear panel of the instrument (Fig. 1, pos. 5, 11, 12 and 13). The RS 1 socket handles the connection in the direction of the PC, while the RS 2 interface handles the connection to other devices, in the daisy chain.

The transmission parameters are set via software commands.

4.2 Installation in an Daisy Chain

The use of an instrument address makes it possible to operate several titration instruments in a chain (daisy chain). For this purpose, a connection is established from the PC to the first instrument in the chain (e.g. TITRONIC® 500) piston burette) using an RS-232-C cable. Another cable is used to establish a connection from the second RS-232-C interface of the burette to the upper RS-232-C interface of the sample changer. Additional devices can then be connected from the second (lower) RS-232-C interface of the sample changer (Fig. 1, pos. 11 and 12).

i It must be ensured that the devices in the chain have different addresses. The device address of the TW 7200 can be set via corresponding commands. The possible addresses range from 00 to 15. Address 3 is set at the factory for the sample changer.

i If the set parameters of the interfaces or the address of the sample changer are no longer known, they can be reset. To do this, press the reset key with a thin object while the sample changer is switched on. This sets the device address to address 3 and the parameters of all interfaces to the standard parameters.

4.3 Data transmission parameters

RS commands can be used to set the transmission speed and other parameters separately for each interface. Besides the transmission speed, the following parameters can be set:

Default parameter of the RS interface:

Data rate	= 4800 Baud
Number of data bits	= 8
Number of stop bits	= 1
Parity check	= even, odd, no

By RS command, the following different speeds can be set: **4800**, 9600, 19200, 28800, 38400 baud.

5 Communication with the TW 7200

5.1 RS command list of the sample changer TW 7200

The TW 7200 sample changer has a series of commands with the help of which it can be controlled by the TitriSoft titration software. If own application programs are to be created, the commands given below must be used.

All commands have the form: "address" (two digits); "command"; "CR"; "LF".

The address is used to address the correct device. Several devices, which have an addressable RS-232-C interface, can be connected to one PC. The addresses can be 00 to 15. In the following list the addresses are marked with "aa" in the general form.

Command	Description	Reply
aaGT	Query of the set sample tray gg = total position / zz = inner positions / cc = ID	aaGTgg;zz;cc
aaSCN	Automatic determination of the placed sample tray size	aaSCNgg;zz;cc
aaDV	Rotate the sample tray one position further	aaDV Y
aaDR	Turn the sample tray back by one position	aaDR Y
aaDPzz	Turn the sample tray to the specified position (zz = 1 to sample tray max.).	aaDP Y
aaPO	Query current sample tray position	aaPOzz
aaKPzzz	Move titration head to absolute head position in percent (zzz = 0 - 100) no cup check when moving the head	aaKP Y
aaGK	Query titration head absolute head position (zzz = 0 - 100)	aaGKzzz
aaKEA	Activate titration head end position up	aaKE Y
aaKEE	Activate titration head end position normal ("center")	aaKE Y
aaKH	Lift the titration head up to the set limit switch	aaKH Y
aaKR	Lower titration head, if no beaker available:	aaKR Y aaKR ERROR:NO BEAKER
aaKGzzz	Move titration head downwards relatively as a percentage of the total travel (end position top to bottom) (zzz = 1 - 100) if no beaker available:	aaKG Y aaKG ERROR:NO BEAKER
aaKUzzz	Move titration head upwards relatively in percent of total travel (end position up to down) (zzz = 1 - 100)	aaKU Y
aaQSz	Set stirring stage for rod stirrer and magnetic stirrer (z = 0 - 9)	aaQS Y
aaQA	Switch off the rod stirrer and magnetic stirrer	aaQA Y
aaQDzz	Preset magnetic stirrer speed (100 to 900 rpm)	aaQD Y
aaGQ	Read out magnetic stirrer speed actual value (zzz = rpm)	aaGQzzz
aaQRSz	Set stirring stage for rod stirrer (z = 0 - 9)	aaQRS Y

aaQRVzzzz	Preset voltage for rod stirrer (0mV = Off / 500-3300mV)	aaQRV Y
aaOE1..4	Switch on the specified outputs (1 to 4). Separate digits by ';' Ex.: aaOE1;3;4	aaOE Y
aaOA1..4	Switch off the specified outputs (1 to 4) Separate digits by ';' Ex.: aaOA1;2;3	aaOA Y
aaIP	Query of the input of the IO interface (z = 0 / 1)	aaIPz
aaRH	Query device identification	aaIdent: TW7200
aaVE	Query software version number	aaVersion: JJKWr
aaGS	Serial number query	aaGSzzzzzz
aaBA	Switch off Pump Unit #1 connection	aaBA Y
aaBE	Switch on Pump Unit #1 connection	aaBE Y
aaBS1..9	Switch on connection pump unit #1 for n-seconds (1-9)	aaBS Y
aaCA	Switch off connection pump unit #2	aaCA Y
aaCE	Switch on Pump Unit #2 connection	aaCE Y
aaCS1..9	Switch on connection pump unit #2 for n-seconds (1-9)	aaCS Y
aaRB	Check if beaker is present	aaRB Y
	if no beaker available:	aaERROR:NO BEAKER
aaSRS<par>	Parameterization of the serial interfaces	aaSRS Y
	<par>: interface;baudrate;data bits;stop bits;parity interface: 1=DC1 / 2=DC2 / 3 =DC1+DC2 / 4 =USB VCP Baud rate: 4800, 9600, 14400, 19200, 28800, 38400 Data bits: 8 (fixed) Parity: even, odd, no Stop bits: 1, 2	
	Example: aaSRS1;4800;8;1;no	
	The parameters are only accepted after a restart (on-off)!	
aaINIT	Re-initialization of all axes of the device Resets occurred errors first	aaINIT Y
aaSR	Stop all movements, switch off stirrers, pumps, outputs	aaSR Y
99AAzz	Numbering of all devices on the daisy chain	
aaAAzz	The first device in the chain gets the address zz. The next device gets the address zz+1.. Wenn zz =15, device gets the address zz+1. If zz =15, then the next device gets the address zz = 0	
	all devices respond with their new address	zzY
99AB..	send any command to all devices	
aaAB..	all devices respond to the command given after <99AB> Ex: 99ABVE returns all version numbers of all connected devices.	

5.2 Network communication

By default, address assignment via DHCP is activated on the TW 7200. This automatically assigns an IP address to the device in an IP network, provided that a DHCP server is active there.

The NWA command can be used to query the current connection parameters as well as to set new connection parameters. During the initial setup it is recommended to set up the address parameters via RS1 or USB. If a network connection with a known IP to the TW 7200 is established, the network parameters can also be changed directly via the network connection.

The UDP (User Datagram Protocol) and TCP (Transmission Control Protocol) protocols are available for direct communication with the device. The communication itself is ASCII-based with the identical commands from the RS command list. For this purpose, the command is written into the data field (payload) of the respective protocol.

The address of the device must still be prefixed despite IP communication. Thus, if required, further devices connected directly to the DC2 interface can be addressed via the TCP connection of the TW 7200. The control characters "CR"; "LF" at the end of the command are not necessary here, but are permitted.

5.2.1 Communication via UDP

The UDP protocol is mainly used to find and identify the device in the network. A forwarding of the UDP commands to DC2 does not take place, the TW 7200 answers the message directly. Even if a TCP connection to the device is already open, UDP requests are still answered. The following list shows the commands available via UDP:

Command	Description
aaRH	Query device identification
aaVE	Query software version
aaGS	Query serial number
aaGI	Query device information
aaNWA	Query / set network settings
aaBLINK	Identify device = LED 4x green flashing

Send message to the TW 7200

UDP port: 50.000 / Destination IP: Broadcast address (.255) or known IP of TW 7200

Command: From above list (e.g. 03RH)

Receive message from TW 7200

To receive the UDP response, the controlling device must provide a UDP server on port 50,000.

UDP-Broadcast

With the UDP broadcast a packet can be sent to a whole subnet. (e.g.: 192.168.0.1 to 192.168.0.254). However, the UDP Broadcast traffic does not traverse multiple networks; it is not forwarded outside the local subnet.

If the TW 7200 is in the same local subnet as the control unit (client), a command from the above list can be sent with the UDP broadcast address (.255), TW 7200 now only responds specifically to the control unit with a UDP packet with its IP address and the desired response of the command in the payload. Thus the initially unknown IP address of the TW 7200 can be determined quickly.

Example:

Action	UDP Payload	Source port	Destination port	Source IP	Destination IP
Control unit sends	03RH	optional	50.000	IP Control unit e.g. 192.168.0.10	xx.yy.zz.255 e.g. 192.168.0.255
TW 7200 answers	03Ident: TW7200	optional	50.000	IP TW7200 e.g. 192.168.0.72	IP control unit e.g. 192.168.0.10

To receive the answer of the TW 7200 the control device must listen on the UDP port 50.000
= provide UDP server

5.2.2 Communication via TCP

If the IP address of the device is known, a TCP connection to the TW 7200 (server) can be established on port 50,000. The device only allows a TCP connection to a control device (client). All commands from the RS command list are supported.

The TCP connection of the client must be persistent, i.e. it must remain open as long as communication is to take place. This enables the TW 7200 to send responses to the client independently via the open TCP connection. It is therefore similar to a connection via serial interface. In some terminal programs instead of a COM port a TCP port can be defined in the same way, which can be used to communicate with the same command list.

Likewise, incoming messages that do not correspond to the address of the TW 7200 are forwarded to the DC2 interface, and the responses from devices connected there are sent back via the TCP connection.

5.3 Extended RS command list - network

The commands listed separately here are also available via USB / RS.

Command	Action	Response from TW 7200
aaMAC	Query MAC address	03MACaa-bb-cc-dd-ee-ff z.B. 03MACD8-80-39-97-D3-D7
aaNWA	Query network parameters	aaNWA <ma>;<ip>;<mk>;<gw>;<dns>
aaNWA<ma>;<ip>;<mk>;<gw>;<dns>	Set network parameters	aaNWA Y
aaGI	Device information	aaGI<id>;0;<ser>;<rh>;<ve>;<ip>;<ma>

Parameter	Designation	Example	Comment
<ma>	DHCP / static IP	M = static IP A = DHCP an	
<ip>	IP address	192.168.0.21	
<mk>	Subnet mask	255.255.255.0	
<gw>	Gateway	192.168.0.1	
<dns>	DNS	0.0.0.0	Not required for TW7200 ; parameter can be omitted when setting (compatibility reasons)
<id>	Device type number	72 (solid	
<ser>	Serial number	12345	
<rh>	Device name	TW7200	
<ve>	Software version	2109	

Examples:

Command	Action	Response from TW 7200
03NWA	Parameter query	03NWA A;10.76.54.115;255.255.255.0;10.76.54.25;0.0.0.0
03GI	Device information	03GI 72;0;1234;TW7200;2106;10.76.54.115;A

Action	Command
03NWA	03NWAA
03GI	03NWAM;10.76.54.211;255.255.255.0;10.76.54.24;0.0.0.0 or: 03NWAM;10.76.54.211;255.255.255.0;10.76.54.24

6 Operating display and error messages

6.1 Operating display

The LED **lights up green** when the sample changer is switched on and ready for operation.

The LED **lights up red** when the sample changer is in motion.

The LED flashes green when error conditions are present.

The LED flashes red when serious operating faults have occurred.

i The respective blink code indicates the malfunction. A distinction is made between normal operation and firmware update (different flashing frequency):

2x Green flashing: Parameter reset by user (reset button) Restart required

3x Green flashing: No plate detected

2x Red flashing: Internal memory error

3x Red Flashing: Error horizontal drive

4x Red Flashing: Error vertical drive

6.2 Errors and messages

The RS commands respond with ERROR if the command could not be executed or there is an error. In this case, the command itself is repeated.

Example: Command "03DP3" not possible, because plate is still rotating

Response: TW7200: "03DP ERROR:BUSY"

Message	Problem/Action
aaBB ERROR:Command or aaERROR:Command	Command not available, syntax or value incorrect
aaBB ERROR:BUSY	Device moves axis and is not ready
aaBB ERROR:KEIN BECHER	NO VIAL No sample vial present with KR,KG or RB command
aaBB ERROR:< number code>	1: Memory error 20: Error vertical drive 30: Error horizontal drive 40-42: Error sample tray drive 43: No sample tray detected / set 100: Other error

aa = Address / BB = Command repetition

6.3 Firmware update

A firmware update can be performed via the USB type A connection using a USB stick.

1. Copy the firmware file (e.g. " TW720x_FW_*.bin ") directly - without subdirectory (root) - to a FAT formatted USB stick. The file must not be renamed. There must not be another update file in the root directory on the USB stick!
2. Switch off the TW 7200 and connect the USB stick to the device.
3. Switch on the TW 7200. After the USB stick has been recognized, the loading of the update file is signaled by the status LED. This flashes red & green alternately, interrupted by repeated red phases.
4. continuous, slow flashing (1Hz) after approx. 10 seconds signals the successful update.

In the event of an error, the status LED flashes quickly red (5Hz). The update must be performed again. To do this, briefly switch off the power supply and restart the TW 7200.

5. After the update has been completed, switch off the device and remove the USB stick.

7 Guarantee

We provide guarantee for the device described for two years from the date of purchase. This guarantee covers manufacturing faults being discovered within the mentioned period of two years. Claim under guarantee covers only the restoration of functionality, not any further claim for damages or financial loss. Improper handling/use or illegitimate opening of the device results in loss of the guarantee rights. The guarantee does not cover wear parts, as lobes, cylinders, valves and pipes including the thread connections and the titration tips. The breach of glass parts is also excluded. To ascertain the guarantee liability, please return the instrument and proof of purchase together with the date of purchase freight paid or prepaid.

8 Storage and transportation

If the device has to be stored over some time, or to be transported, the use of the original packing will be the best protection. However, in many cases this packing will not be available anymore, so that one will have to compose an equivalent packaging system. Sealing the devices in a foil is hereby recommended. The devices should be stored in a room with a temperature between + 10 and + 40°C, and the (relative) humidity of the air should not exceed 70 %.

9 Recycling and Disposal

Please observe the applicable local or national regulations concerning the disposal of "waste electrical and electronic equipment".

The devices and their packaging have been manufactured as far as possible from materials that can be disposed of in an environmentally friendly manner and recycled properly. If you have any questions about disposal, please contact our service department (see back of these operating manual).

10 EC – Declaration of Conformity

The corresponding declaration of conformity of the device can be found on our homepage. It will also be made available to you on request.

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SI Analytics is a trademark of Xylem Inc. or one of its subsidiaries.
Xylem Analytics Germany Sales GmbH & Co.KG
© 2021 Xylem, Inc. Version 211112 M 825 318 6