INSTALLATIONSANLEITUNG

Originalversion

INSTALLATION GUIDE

WinVisco

Software zum AVS[®] 370
Software for the AVS[®] 370

SI Analytics

a **xylem** brand

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Important notes:

The installation guide does not replace the operating manual on the attached CD. Before initial operation, please carefully read and observe the installation guide and keep it. Please also consider the operating manuals for the devices to be connected.

All specifications in this installation guide are guidance values which are valid at the time of printing. However, for technical or commercial reasons or in the necessity to comply with the statutory stipulations of various countries, the manufacturer may perform additions to the product without changing the described properties. A potentially more recent version of this installation guide is available on our internet website. The German version is the original version and binding in all specifications!

The menu screens serve as an example and may differ from what you see!

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TABLE OF CONTENT

1	Properties of the WinVisco Software	13
1.1	Intended use.....	13
1.2	Warning and safety notes	13
1.3	System Requirements	14
2	Installation.....	15

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Deutschland, Printed in Germany.

1 Properties of the WinVisco Software

1.1 Intended use

The WinVisco Software serves for the control of the measurement device AVS® 370. The AVS® 370 carries out automatic measurements with SI Analytics® capillary viscometers and determines absolute and relative kinematic viscosities. The operation of the WinVisco Software takes place with a personal computer (PC). The software controls the measurement sequence automatically. The determined measurement values are included in further calculations and are stored in a database. Different documentation possibilities visualize the respective data. User profiles must be created for working and operating the software. The operating elements are only released through logging in of a user.

The WinVisco software can control up to 8 ViscoPump modules. The modules can be operated in parallel. A flexible and fast viscosity determination with different configurations and samples is possible hereby.

In addition, the WinVisco software can control up to 8 small devices (burettes or titrators).

Devices of the company Schott® Geräte, Schott® Instruments, SI Analytics® and Xylem Analytics Germany are exclusively supported as measurement devices. Measurement devices of third parties are not supported.

 The software WinVisco has predecessor versions until version 3.83.00. These versions are not compatible with the present version 4.x.x.x and above.

1.2 Warning and safety notes

 **The connected devices are controlled via the WinVisco software!**
Please also observe the operating manuals for the devices to be connected.

 **When changing the work mode (sucking or pressing) of the ViscoPump** (see  section 5.3.2) **the following has to be considered:** With a wrong operating type, the sample liquid can escape from the viscometer and e.g. reach the surrounding thermostat bath or the ViscoPump. **The hose connection of the system must be adapted in particular!**

1.3 System Requirements

The product is subject to various software and hardware requirements. The minimum requirements must absolutely be fulfilled for a smooth operation.

i Contact your IT administrator prior to installation, if you are not clear about the specification of your PC system.

Minimum requirements

Operating system (OS)	from Microsoft Windows7 SP1 \ 8.x \ 10
Processor (CPU):	1 GHz
Memory (RAM):	1 GB (with 32-Bit), 2 GB (with 64-Bit)
Hard drive (HDD):	up to 2.5 GB of free disk space*
Graphics (GPU):	128 MB DirectX 9-capable graphics card with at least 1024x768 pixel resolution
Connections:	at least one RS-232 Type C interface (COM)

* WinVisco requires free storage space as a pure software product. The storage requirement increases if components necessary for the operation are missing on the target system. The setup recognizes missing software components, notifies this and offers the possibility to install it as well.

Recommended requirements

Operating system (OS)	Microsoft Windows 10
Processor (CPU):	Intel Core 2 @ 2 GHz AMD Athlon X2 @ 2GHz
Memory (RAM):	4 GB
Hard drive (HDD):	up to 2.5 GB of free disk space*
Graphics (GPU):	256 MB DirectX 10-capable graphics card with at least 1280x1024 pixel resolution
Connections:	at least one RS-232 Type C interface (COM)

Further requirements

Microsoft Access Database Engine 2010	(can be installed automatically through the setup)
Microsoft .NET Framework 4.6.2	(can be installed automatically through the setup)

2 Installation

WinVisco is made available of different data carriers.

The setup starts automatically when installing a CD.

When installing a USB stick or other sources, the setup must be started manually.

i The setup is exclusively provided in English.

If required components should be missing, the installation routine starts with the following window:

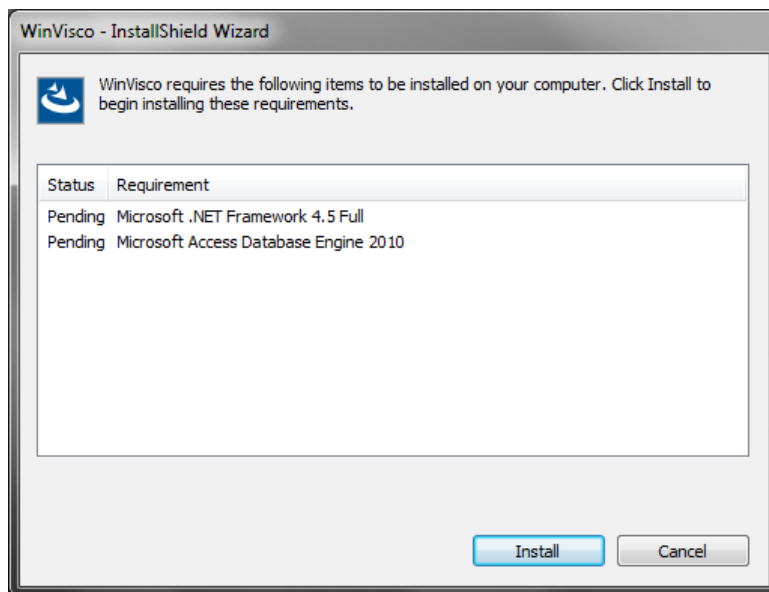
Fig. 1 Installation of necessary software components

By means of clicking on “Install”, the necessary components are installed.

After the successful installation of the software components (or if these were already present), the installation routine reports with the following window:

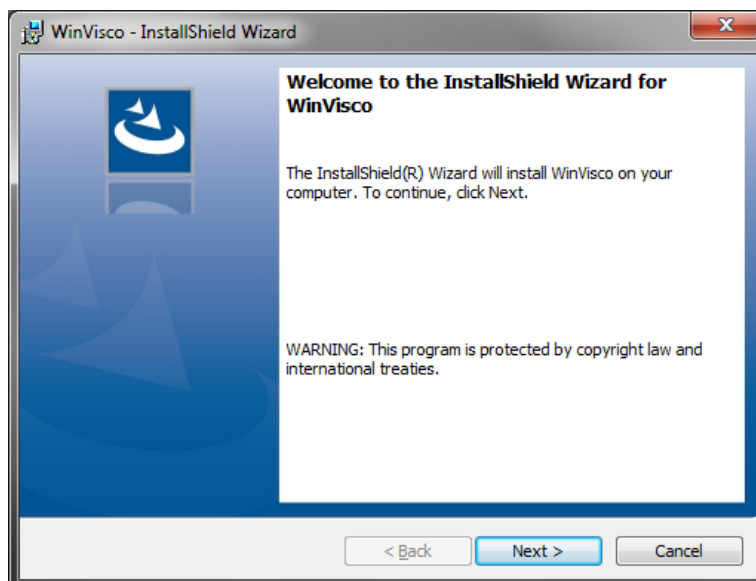
Fig. 2 Start of the installation routine

Via “**Next**” you are forwarded to the license agreements:

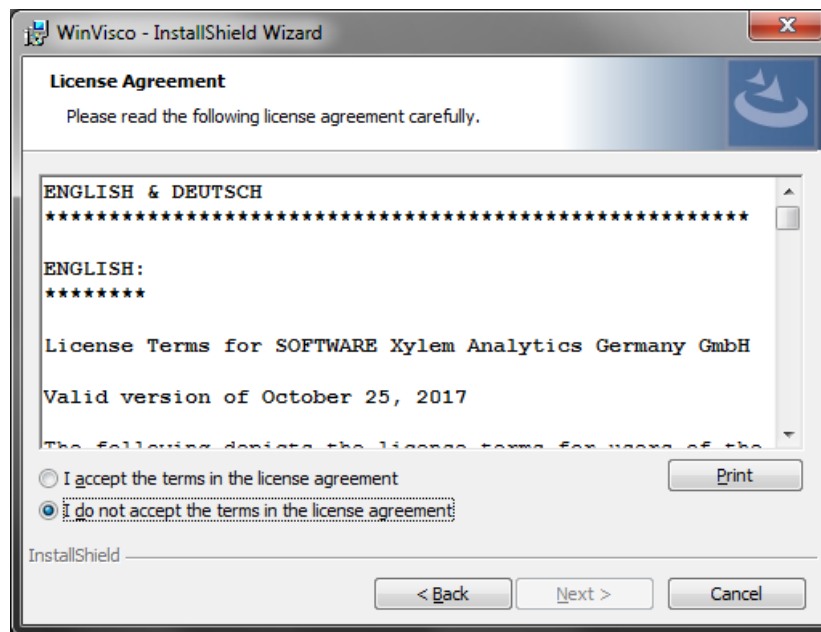
Fig. 3 License agreements prior to product installation

With agreement and “**Next**”, the installation routine is ready for the actual product installation:

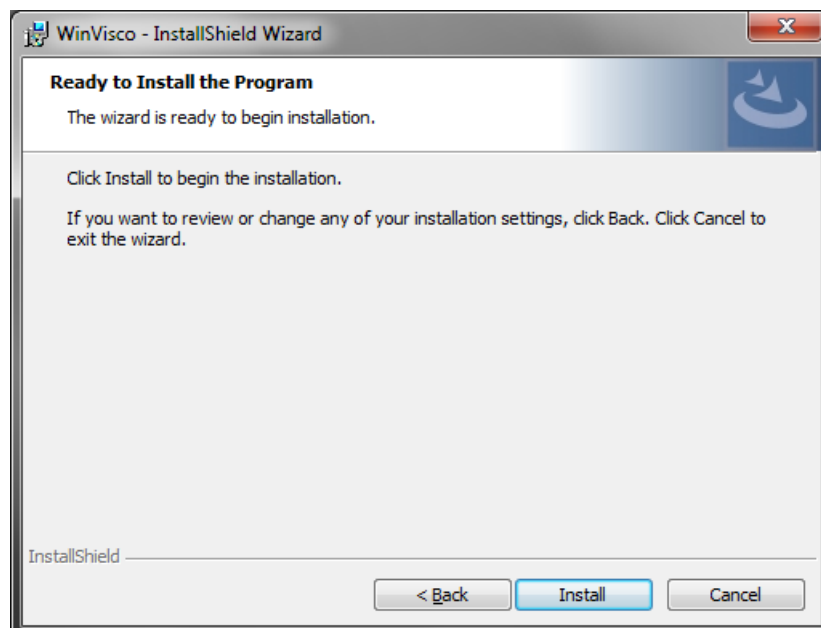
Fig. 4 Start of the software installation

WinVisco is installed on your system by means of “**Install**”.

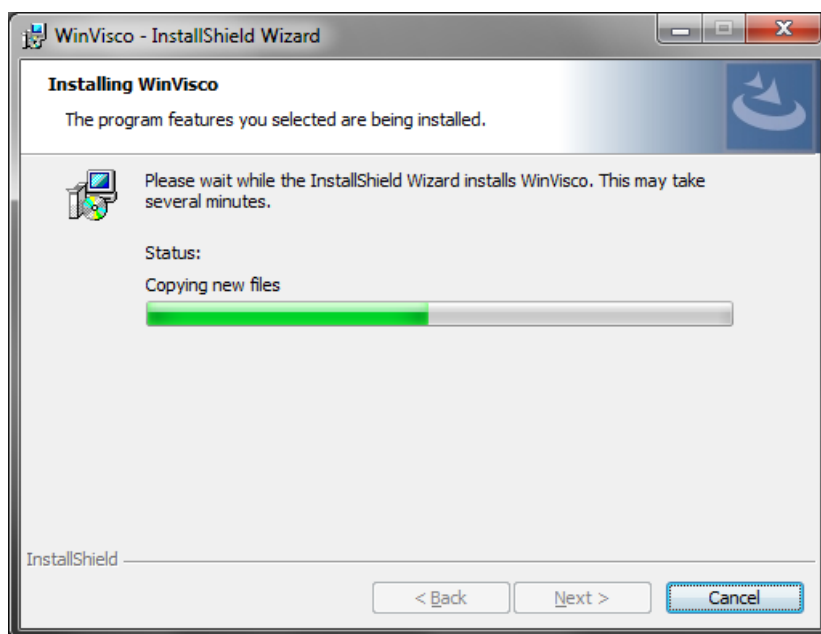
Fig. 5 Product installation

The successful installation is confirmed to you in the last window:

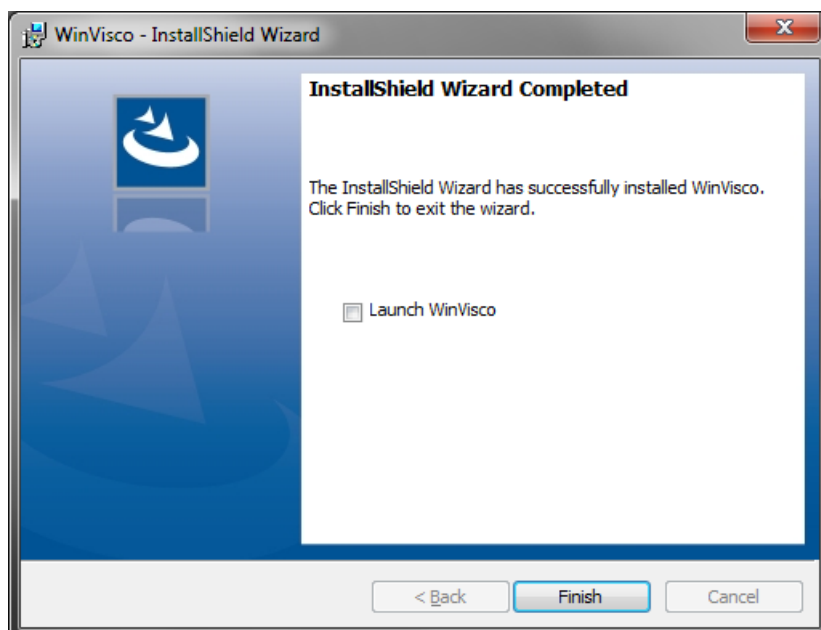
Fig. 6 Confirmation of the successful installation

i If you desire an automatic WinVisco start, activate the tick at “**Launch WinVisco**”.

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Tel. +49(0)6131.894.5042

E-Mail: Service-Instruments.si-analytics@xylem.com

SI Analytics is a trademark of Xylem Inc. or one of its subsidiaries.
© 2025 Xylem, Inc. Version 250918 **M 827 210 0**

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Xylem Analytics Germany GmbH
Xylem Analytics Germany Sales GmbH & Co.KG