SOFTWARE UPDATE

Originalversion

SOFTWARE UPDATE

AVS[®] Pro

Update (Version 22/04) zur Software
Update (Version 22/04) for Software

SI Analytics

a **xylem** brand

Update zur Software AVS® Pro Seite 3 ... 30

Update for Software AVS® Pro..... Page 31 ... 58

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The screenshot shows the 'Messparameter' dialog box with several sections:

- Parameterermittlung:** Radio buttons for 'Einzeleingabe', 'Gemeinsame Eingabe' (selected), and 'Blindwert ermitteln'.
- Beginn des Messvorgangs:** Radio buttons for 'Spülvorgang' (selected) and 'Proben transfer'. Below is a field 'Nur Proben ab Nr.' with the value '1' and a 'verarbeiten' button.
- Soll - Temperaturen:** Two input fields for 'Temperatur 1' and 'Temperatur 2', both set to '25,00'.
- Probenzuordnung:** A field 'Anzahl Proben' with the value '1' and a checkbox 'Indiv. Zuordnung'.
- Probenträgergröße:** A field 'Probenträgergröße' with the value '16'.
- Sicherheitsentleerung in...:** A list of radio buttons for 'Viskosimeter 1' through 'Viskosimeter 8' and 'Abfallstation'. 'Viskosimeter 1' is selected.

 At the bottom, there are three buttons: '< Zurück', 'Starte Messungen', and 'Messungen verzögert starten...' (highlighted with a red circle).

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- A time selection area with 'Um' followed by a time picker set to '04:00' and 'Uhr'.
- A date selection area with 'am' followed by a date picker set to 'Mittwoch, 21. Januar 2015'.
- The text 'die Messungen starten.' followed by an 'OK' button.
- The text '(Wartezeit 14,55 Stunden)' followed by an 'Abbrechen' button.

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Copyright

© 2025, Xylem Analytics Germany GmbH

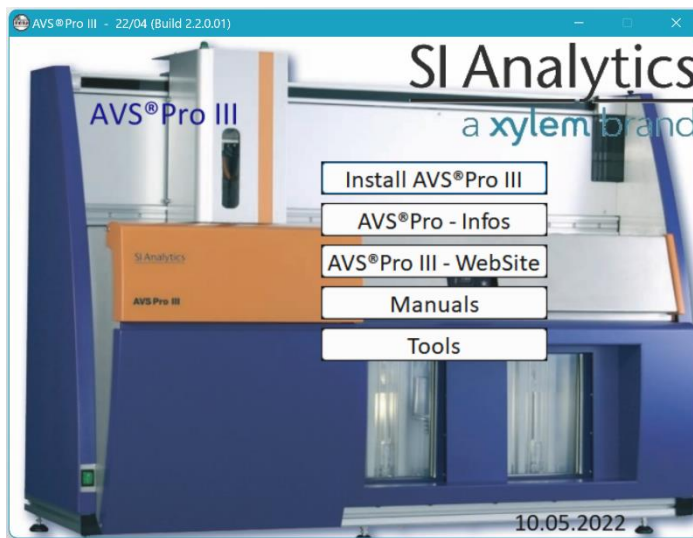
Reprinting - even as excerpts - is only allowed with the explicit written authorization.
Germany. Printed in Germany.

1 Installation of the software


Important notes

- The installation program can be used for a new installation, as well as for the software upgrade of an existing AVS® Pro system.
- If the installation program finds an existing AVS® Pro software on the PC, the data and settings of the existing version are automatically transferred. If you don't want the automatic transfer, uninstall your existing version completely before installing the new version!
- If the new software should be installed on a new PC, the data and settings can be taken over on the new PC semi automatically in the new software installation. In this case find a detailed description on the installation CD (\Tools\Backup_Restore_Migration\).

1. Insert the installation CD-ROM. The following window appears

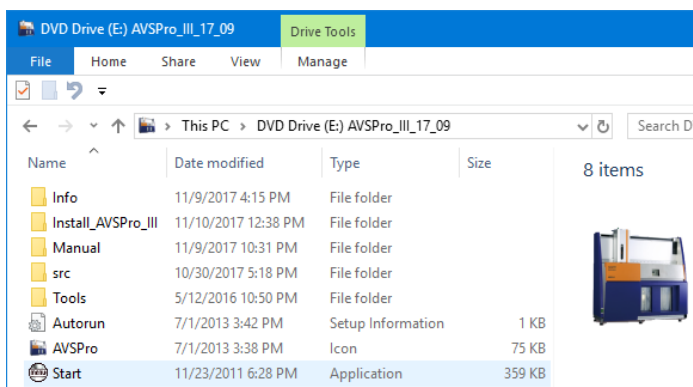
If this window does not appear, you must open the CD drive with a double click and run the file "Start.exe".


2. Click on «**Install AVS® Pro III**» to start the software installation.

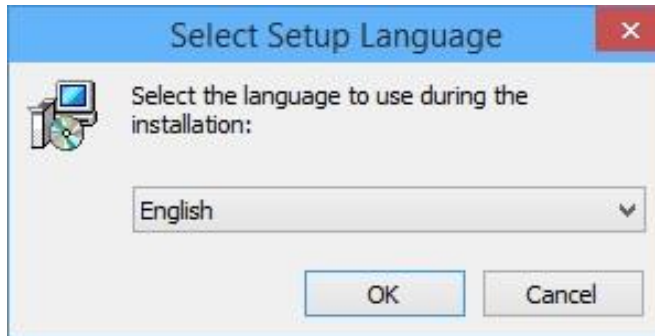In the next window you can choose between English and German.

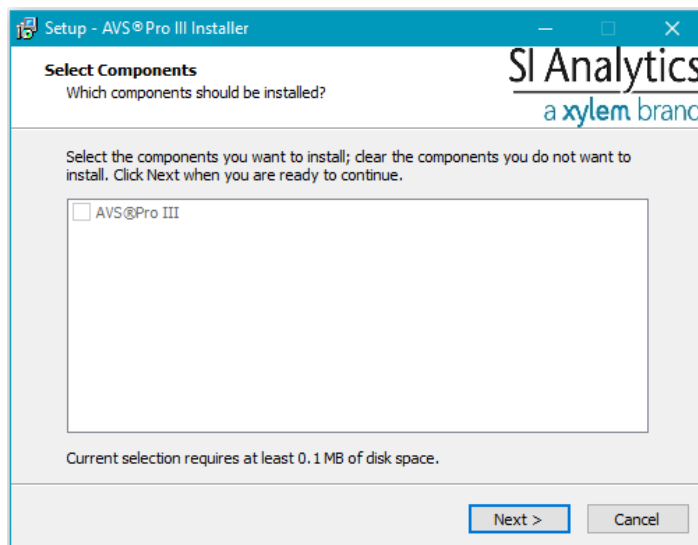
Select the desired language, then click «**OK**».


Confirm with «**Next**».

3. If an older installation of the AVS® Pro software or its data and settings is found on the PC, an automatic backup starts.

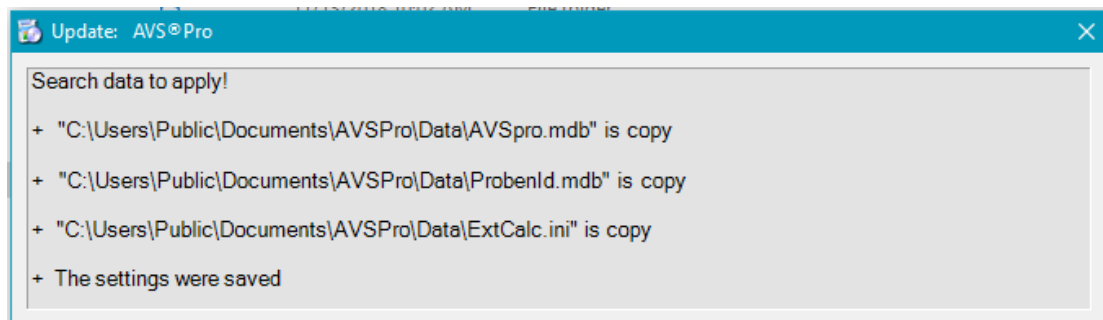
The short result of this search will be displayed:

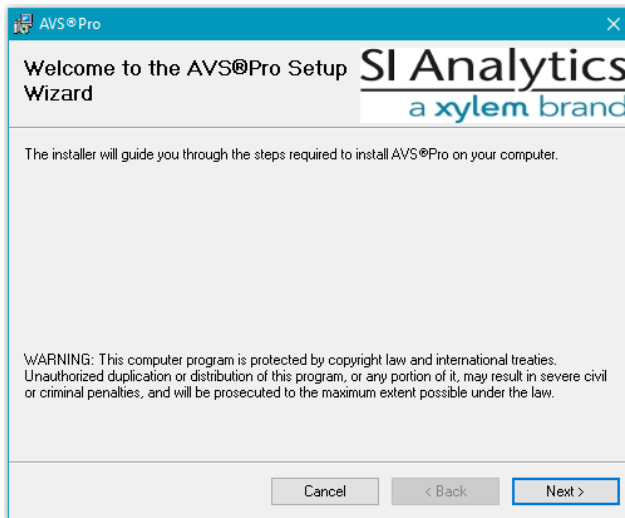
4. Now, the actual Setup Wizard must be started by clicking „**Next**“:


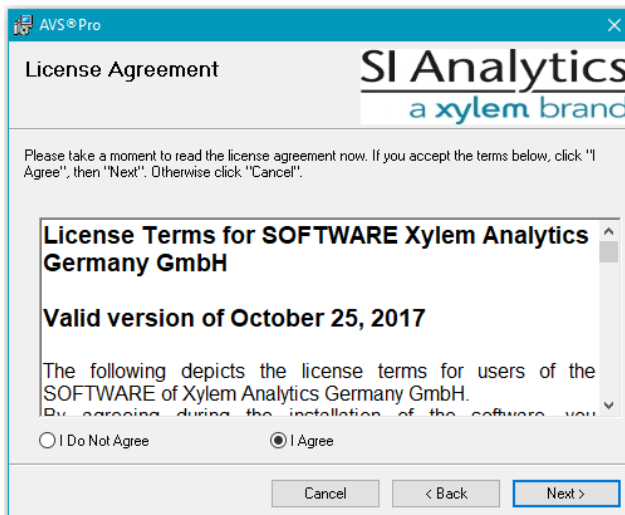
5. Then you have to agree to the license terms:

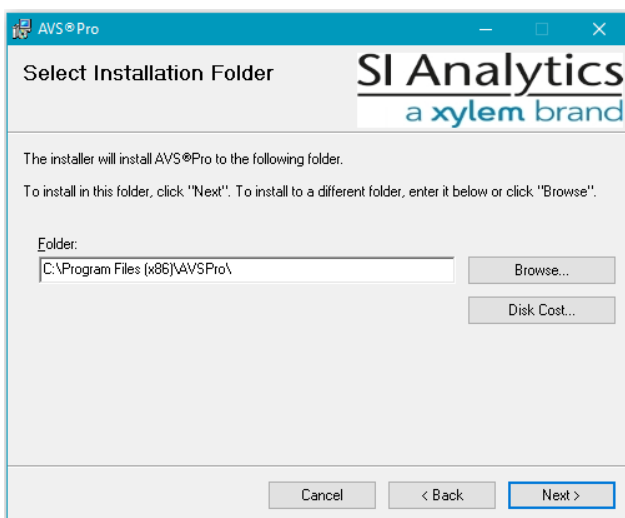
6. It is recommended accepting the installation folder as it is proposed here:

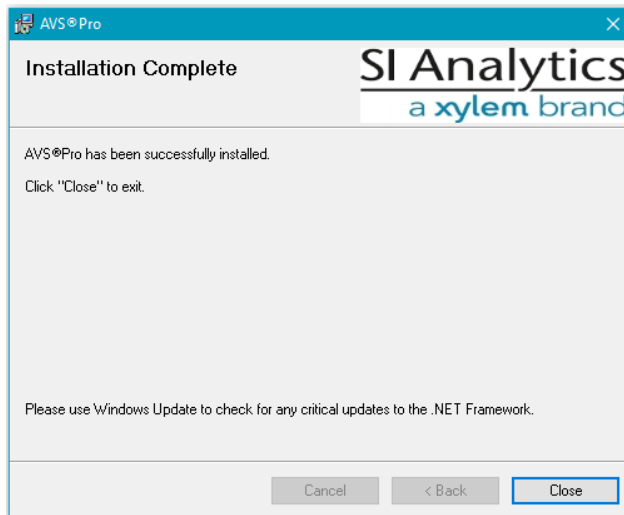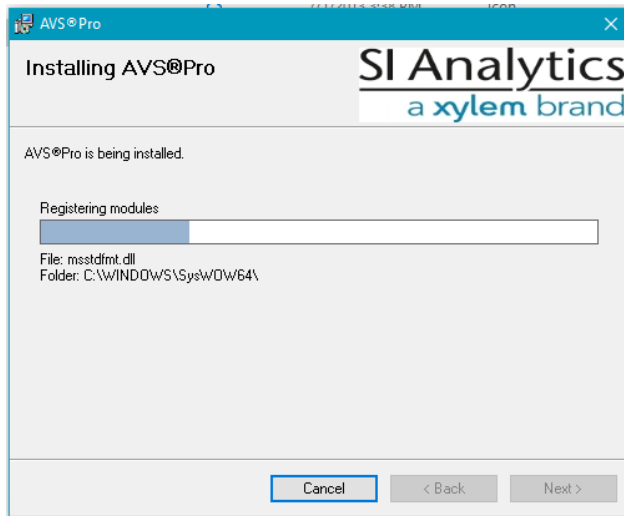7. The installation process runs automatically and the data & settings saved at the beginning are transferred:

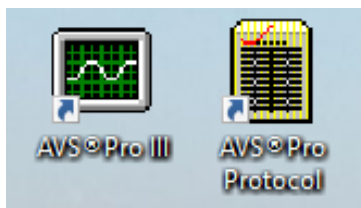8. If the installation has been completed successfully you find two icons on the desktop:


9. The icon „AVS® Pro III“ can be used to start the software.
The icon „AVS® Pro Protocol“ can be used to start the protocol.

2 First launch of the new AVS® Pro software

Important notes


- From version 13/47 of the AVS® Pro PC software on, a separate license code is required for the software.
- The purchased license code also applies to all future versions of the AVS® Pro PC software whose file version number has the same major number (e.g. "2.") hat. Only if the file version no. starts with "3." a new license code is obligatory.
- If you have upgraded from an already licensed version (for example: 13/47), the license information has been automatically transferred and will in general not have to be entered again into the system.

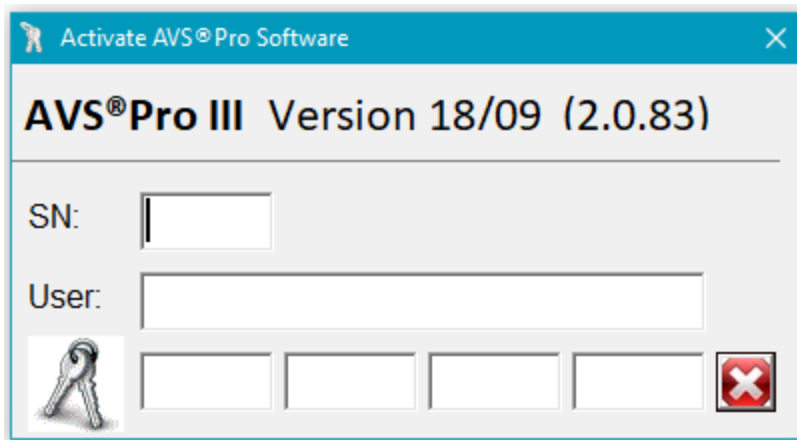1. If no license information can be found, the following window appears at your first start of the software:


Enter here the Serial number, the User information and the license code.
You will find this information in the license document.

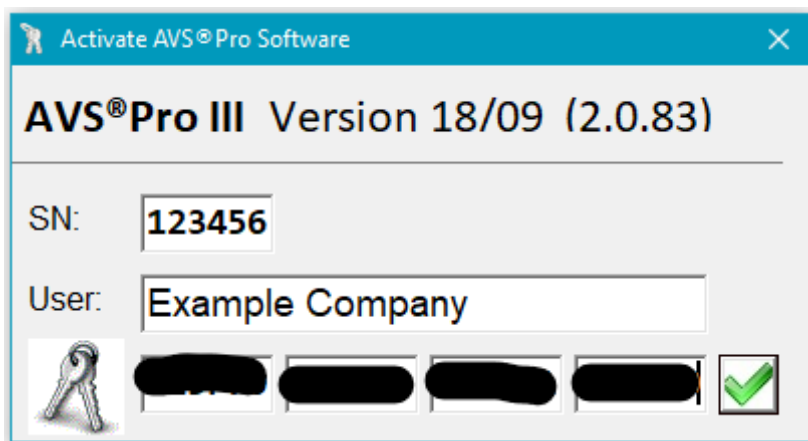2. After entering the correct data, a green checkmark will appear instead of the red cross.


Click on the green checkmark to permanently take over the license information!

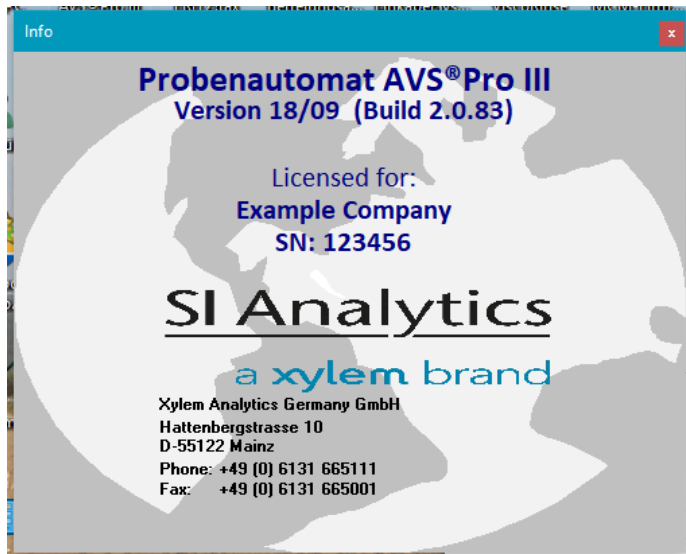3. No license scanning is performed from the second launch of the software onward. Instead, the Info window comes up with information of the licensee and the serial number immediately

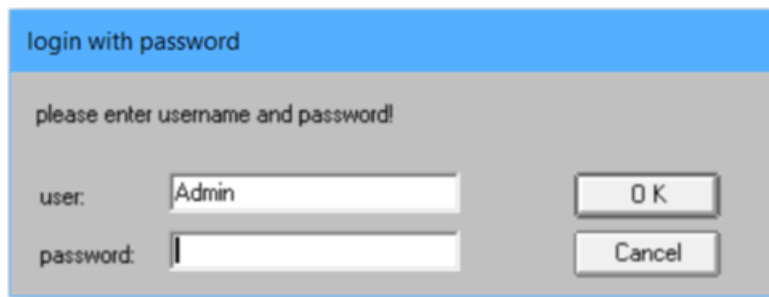
4. Then, the password login will appear:


5. By default, the administrator has the password **"Admin"**.
- If you have upgraded from a previous version, and have already changed the administrator password, please enter this one. As an "Administrator", you have all access rights!
 - There is also a default user **"User"** with password **"u"**, which has only restricted access rights.
 - The password can be switched off for more software launches in the menu «**master data**» → «**user**». Then the admin account is automatically used.
Refer to the chapter "User administration" for more information.

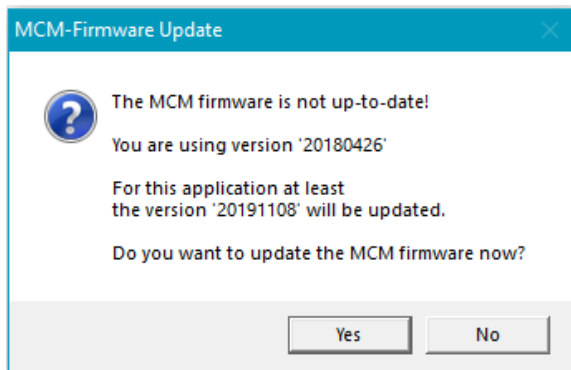6. Starting with version 19/09, the check is performed every Wednesday when the application is started, whether a firmware update for the “MCM step 420 module” should be carried out. If so, the following message appears:

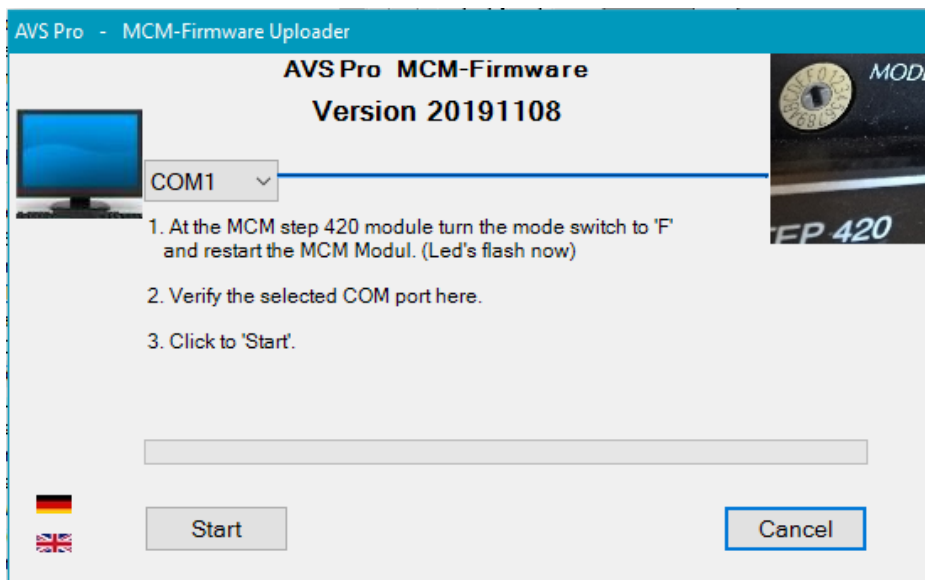With a click on «Yes» the MCM-Firmware Uploader is started and the AVS® Pro application closed:

If «No» was clicked here, the firmware update can also be started at any time via to start the info dialog:

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- User management is only available if you are logged in as administrator
- The administrator password can be changed in the user management.
- In the user management users with restricted rights can be created, changed (including password) or deleted.

Overview of how user management works:


	User password: switched on		User password: switched off
Program start	User name and password are asked for. Last user name is already filled in.		No user and no password query Registration is done with user "Admin"
Activate Password query	--		Admin password is queried
Deactivate Password query	Admin password is queried		--
Menu „master data => user“	active		inactive
Logged in as:	Admin	User	(Admin)
User management	• Create a user • Set passwords • Change username or password • Delete user	Access blocked	<u>Not</u> available
Missing menus in the main window			file => new login
Locked menus in the main window		processing => methods... parameters thermostatic extra => options... Profile-Management... use password master data => viscometer user	
Locked menu in the measurement window		processing => ViscoPump parameters... extra => show temperature warnings rinsing ViscoPump active mode blow out capillary Speed mode	
Locked menu options in the window "devices connected"		file => search extra => dosing parameters	

4 Profile management


See profile management in the main menu **«Extras»** → **«Profile Management »**

- If the settings of the software must be modified for different applications, you can save these changes with the profile management. You can then quickly and easily switch back and forth between the different settings.
- The respective profile stores all settings relevant to the measuring sequence such as rack size, solvent, rinse mode, sample label, ViscoPump/dosing parameter etc.

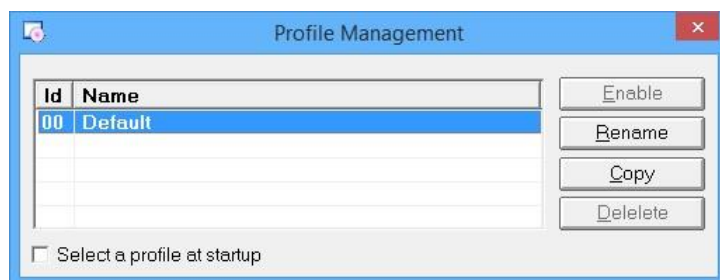The default profile is called "Default" and can be renamed at any time.

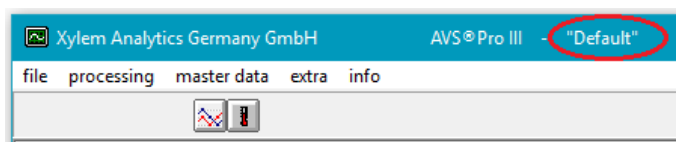The name of the currently used profile is displayed in the header bar of the program window:

Optionally, you can select the profile to use to at each program start. For this, insert the tick at **«Select a profile at startup»**.

If this tick is not set, the program automatically starts with the profile used last

Create a new profile

- Copy an existing profile, such as the "Default"-profile.
- Give the new profile a meaningful name.
- Select the new profile and activate it by using the button "Enable"
- Make all relevant changes to the settings for this new profile.

Change profile

- You can change the profile without restarting the program.

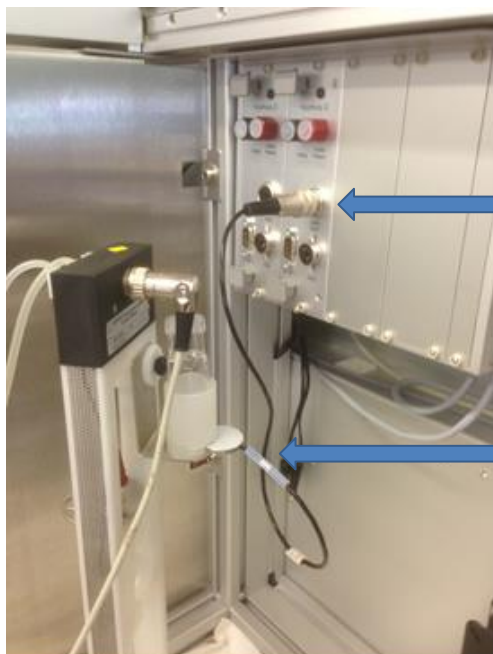
Delete profile

- The current profile cannot be deleted.
- To delete a profile you need to activate a different profile.
- You cannot delete the default profile (ID 00).

5 Monitoring the security bottle by ViscoPump

Important note


- If your AVS® Pro system is equipped with ViscoPumps type II (VZ 8511/ VZ 8561) or type III (VZ 8561), safety bottles can be monitored directly by the respective ViscoPump.
1. Connect the plug of the overflow sensor (VZ 8552) with the socket of the corresponding ViscoPump:


Plug of the overflow sensor

The sensor has to be adjusted (trigger point) with 3 - 5 ml of the used solvent.

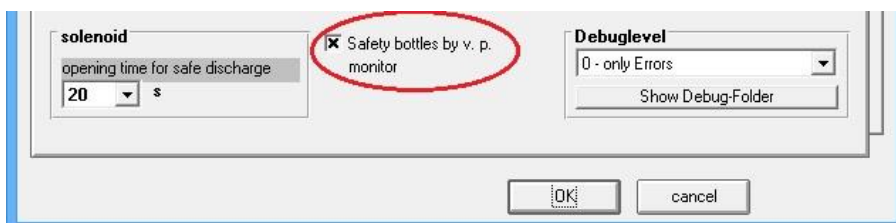2. Main menu: **«Extras» → «Options» → «Common»**
Select the option **«Safety bottles by v.p monitor»**

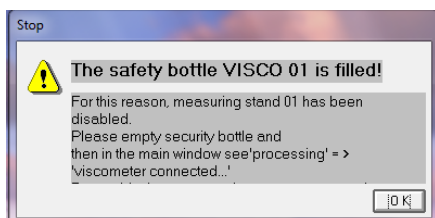
3. If an amount of solvent is in the safety bottle and the sensor has been you will get this message:

Empty the safety bottle and re-enable the ViscoPump affected.
(Menu: **«Processing» → «Viscometer connected ...»**).

6 New Options «Common»

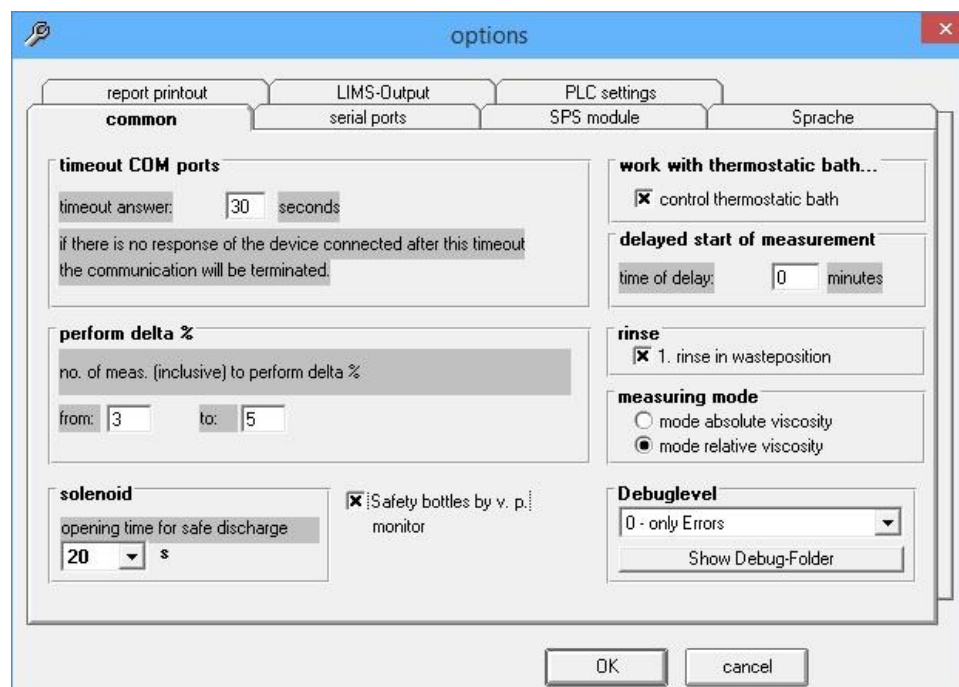Main menu «Extras» → «Options...» → «Common»

«1.Rinse in waste position»

- from the version 10/48 up.

«Opening time for safe discharge»

- since version 12/50 programmable.

«Debug level»

- In consultation with the support team a debug level can be used for problem analysis in case of an error. When a debug level is > 0 the communication between PC and machine is logged in text files. The log files can be accessed via the button "Show Debug-Folder".

«Safety bottles by v.p. monitor»

- is available since version 13/05, for more details refer to chapter «Monitoring the security bottle».

7 New Options «PLC settings»

Main menu «Extra» → «Options...» → «PLC settings»

«Rinsing next sample»

- Values for rinsing with the next sample can be pre-defined, which differ from the values used for the actual measurement.
- When the values entered here meet the defined input of 0.00, the same values are used for rinsing the sample as for the measurement of the sample.
In this case the values from main menu will apply:

«Processing» → «Viscometer connected» → «Extra» → «Dosing Parameters».

«Rinsing: solvent 1», «Rinsing: solvent 2»

- Specifications for the rinsing with external fluids.

«Vacuum drying times» (Only for AVS® Pro III!)

- When using the rinse mode “Vacuum drying” the drying time for the metering module determines, how long the dispensing module has to stay in the dry port. The drying time for the viscometer determines the minimum pumping time of ViscoPump when rinsing with solvent. It only works if “Vacuum drying” is specified as next step of rinsing.

«Flap switch check» and «Check vacuum»

- These options specify, whether the sample chamber door safety switch and the vacuum of the waste system should be checked when the program starts. (**This only applies to AVS® Pro I with VZ 71xx vacuum pumps**)

«Immersion depth»

- Here, the respective Immersion depth of the sample needle is specified. For the vials VZ 7080 and VZ 7081 a depth of 125 mm should be set.

«Dosing unit»

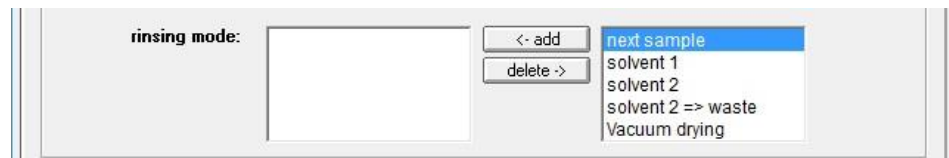
- For the quantity data to be taken into account correctly, if using a micro-dosing unit VZ 7032, please change dosing unit settings here.

«Use dosing prefill option» (Only for AVS® Pro II + III!)

- Pre-filling the dosing unit with the sample while the previous rinsing is still running (Requires at least MCM version 20180426).


8 New settings «methods»

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
«Rinsing mode»

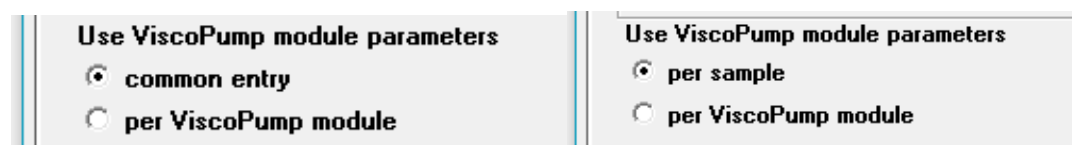- «**Solvent 2**» → «**Waste**»
In this rinse only the dosing unit is rinse with external solvent 2 and the solvent is then emptied into the waste position.
- «**Vacuum Drying**»
This rinsing step can be used after washing with a solvent for intensive drying of the dosing unit. The dispensing needle is inserted into the dry port for the time defined in the options.


«Use ViscoPump parameters»

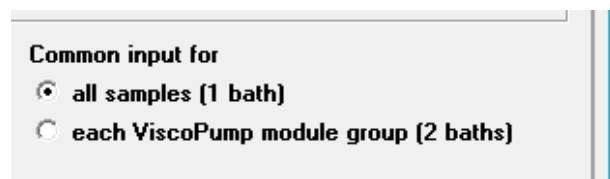- Both «**Common entry**», as well as «**Individual entry**», allows the possibility that the ViscoPump parameters are used per ViscoPump.
The corresponding parameters can be in the menu in the dialog «**Processing**» → «**Parameters of the VisoPump modules...**».

«Common input»

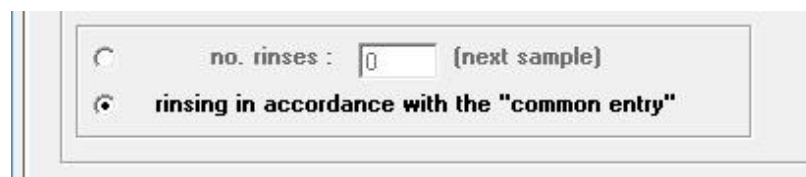- The «**Common input**» can take place separately for two groups of ViscoPumps.
That is important, e.g., if two thermostatic baths with different temperatures are used


«Rinsing determination of t_0 -flowtime»

- On the « **t_0 -flowtime** » tab is the option to adapt the rinse mode, as programmed in the «**Common entry**».
(This option is only available with the AVS® Pro III!)

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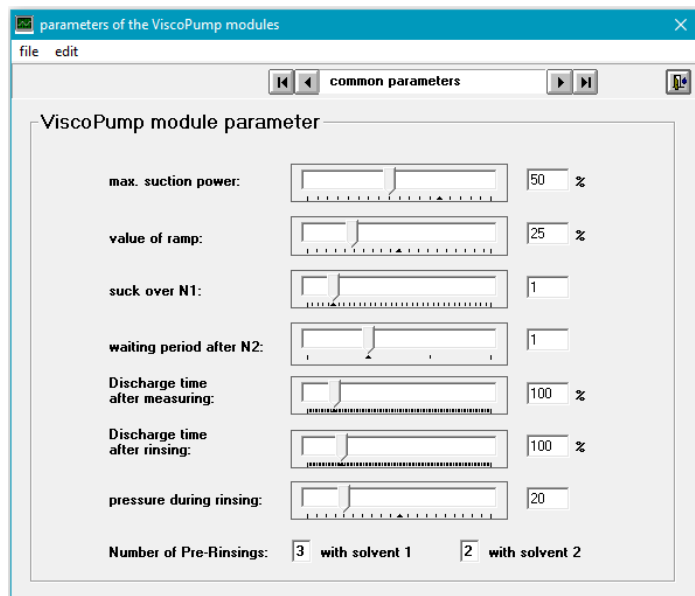«Discharge times»

- from vers. 13/03 an different times for the discharging can be set according to a measurement or after rinsing.

«Parameters for»

- So far, values per sample, could be set here for the «**common entry**» und die «**determination of t_0 -flowtime**».
- As of version 13/26 there is the possibility of defining your own always for each ViscoPump. In the dialog «**selection of methods**» you can choose whether to use the values for ViscoPump or the values per sample. Another option is to leave it at «**common entry**».

«Number of Pre-Rinsing»

- This option is available from vers. 14/12 on.
- Pre-rinsing refers to the pumping of the solvent into the ViscoPump in a washing cycle before it is sucked again.

10 Starting measurement up to 3 days delayed

As of version 13/40, there is the possibility of starting of measurement up to 3 days delayed

Thus a series of measurements can be prepared e.g. on Friday with an automated start on Monday morning.

1. After activating the button «**delayed measurement start**» the following dialog appears:

Set the desired time to start here and then click «**OK**».
Press «**Cancel**» to return to the measurement parameters.

2. Before the actual start of measurement, this dialog will be shown:

3. Waiting can be canceled however at any time with the button «**Cancel**» which brings you back to the measurement parameters.

Important note

- The delayed measurement option exists regardless of the setting in «**delayed start of measurement**» in the common options. The adjustable time is max. 300 minutes (5 hours) and applies with each starting is of the measurements.

11 New settings «measuring parameters»

In the measurement parameters there are 3 new setting options:

«Emptying viscometer»

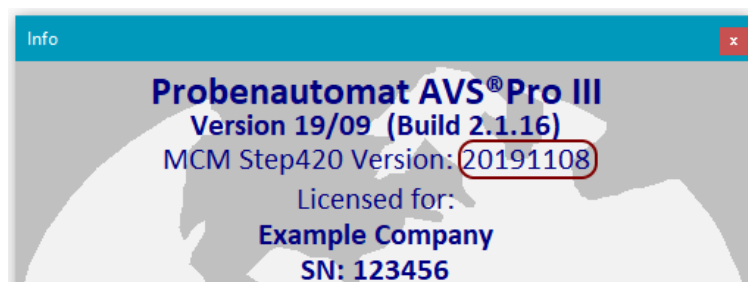
- The setting «**before each sample**» is used to set an additional action to empty the viscometer.
- For the first sample of each viscometer, this additional action, is suppressed because a safety drain has just been carried out before.

«Needle park position» (Only for AVS® Pro II and III!)

- This setting can be used to specify for safety reasons that the dosing unit is parked with the needle in the waste position after the completion of a series of measurements. This is a safety precaution.
- This setting is only available if the MCM version 20171006 or higher has been installed.

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
- With this setting a solvent rinse is initiated for each viscometer involved in the measurement series after the last sample.
- This setting option is only available if:
 - The «Common entry» method is active.
 - The «Common entry» method contains at least one solvent rinse.
 - At least the MCM version 20191108 is installed.

12 New menu «Extras» → «Service»

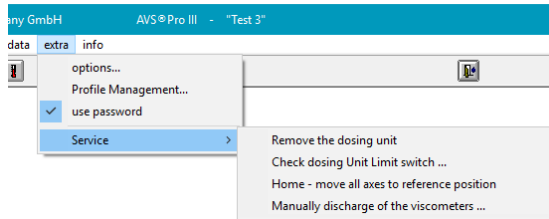There are 4 new menu items in the main menu under «**Extras**» → «**Service**»:




Important note:

- Applies only to AVS®Pro II and III!
- These menu items are available only if you are logged in as administrator


«Remove the dosing unit »

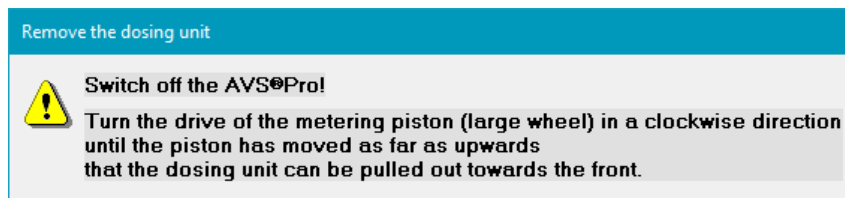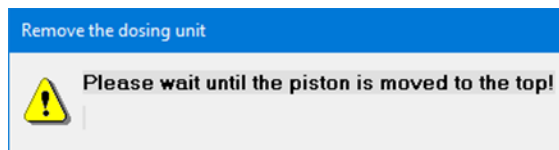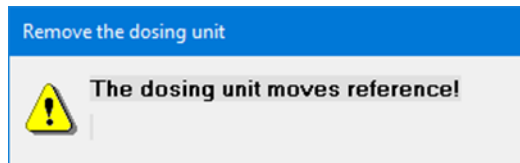- This menu item supports the removal or change of the dosing unit and replaces the additional software AVS®Pro RDU (Remove Dosing Unit) which used to be necessary previously.
- After selecting this menu item, the following windows are displayed:


«Check dosing Unit Limit switch ...»




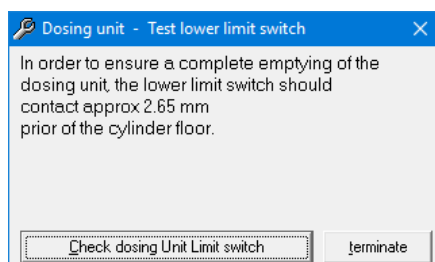
- This menu item is only available if the MCM version 20170919 or higher has been installed.
- This menu item can be used to check the optimal setting of the lower limit switch of the dosing unit. A complete emptying of the dosing unit can only be ensured if the lower limit switch of the dosing unit is optimally adjusted.
- After selecting this menu item, the following dialog appears:

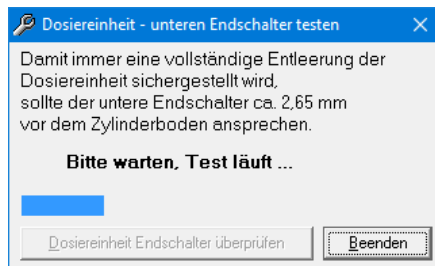- After you click on the „Check dosing Unit Limit switch“ button, you will see the following:

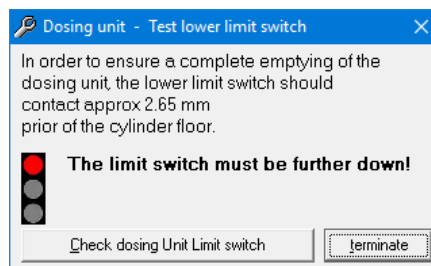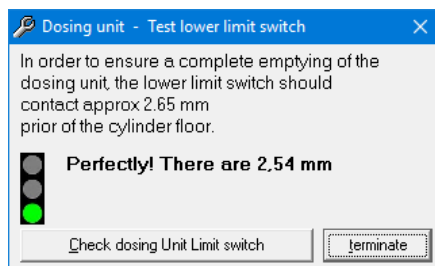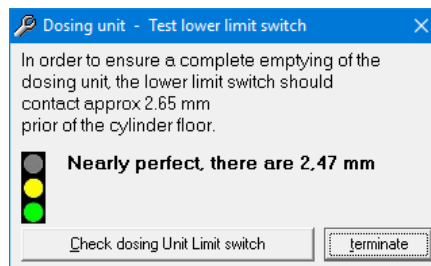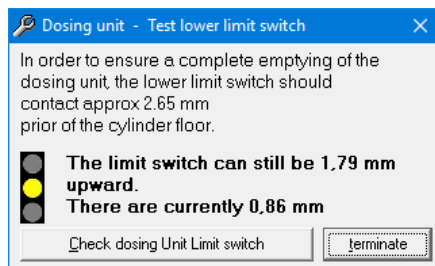
- This is how the test result could be displayed:


«Home – move all axes to reference position »

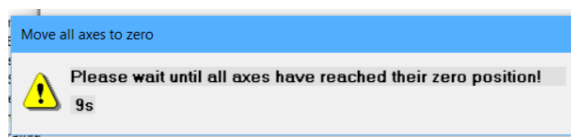- With this menu item, all axes, except of the dosing piston, can be referenced and moved back to their starting position.

«Manually discharge of the viscometers ...»

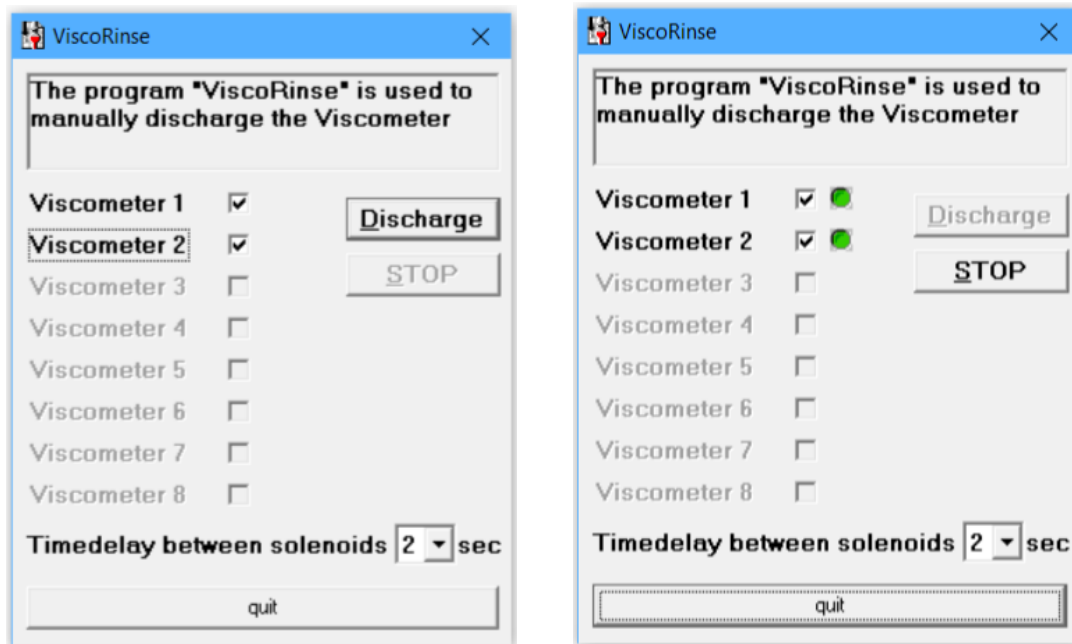
- This menu item starts the ViscoRinse program which allows a selection of viscometers to be manually discharged

- With the option-checkbox near the texts “Viscometer 1-8” you can preselect which viscometers should be rinsed.
- These will be rinsed one-by-one after clicking the button “Discharge”.
- You can freely choose the “Time delay between solenoids” from 1 to 9 seconds.
- The solenoids remain open until the button “Stop” is selected which will close it.


13 New in the «Measurement» window

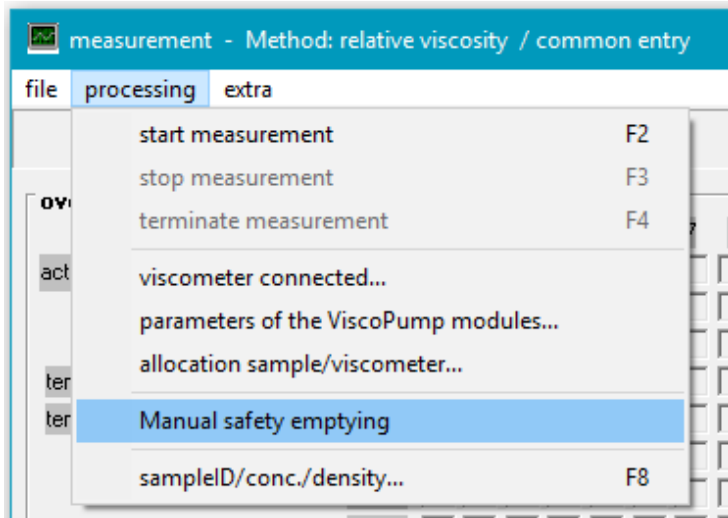The subject line of the «Measurement» window shows the current method.



New menu option «Manual safety emptying»

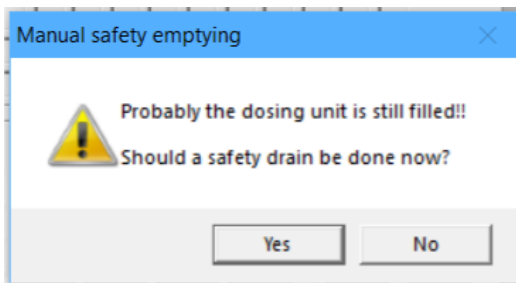- If the dosing unit is still filled after the cancellation of a series of measurements, a safety discharging can be started manually with this menu item.
- This menu item is automatically activated when closing the “measurement” window if it is recognized that the dosing unit may still be filled.
In this case the following security query is shown first:

14 New calculations “IV Billmeyer” and “SV value”

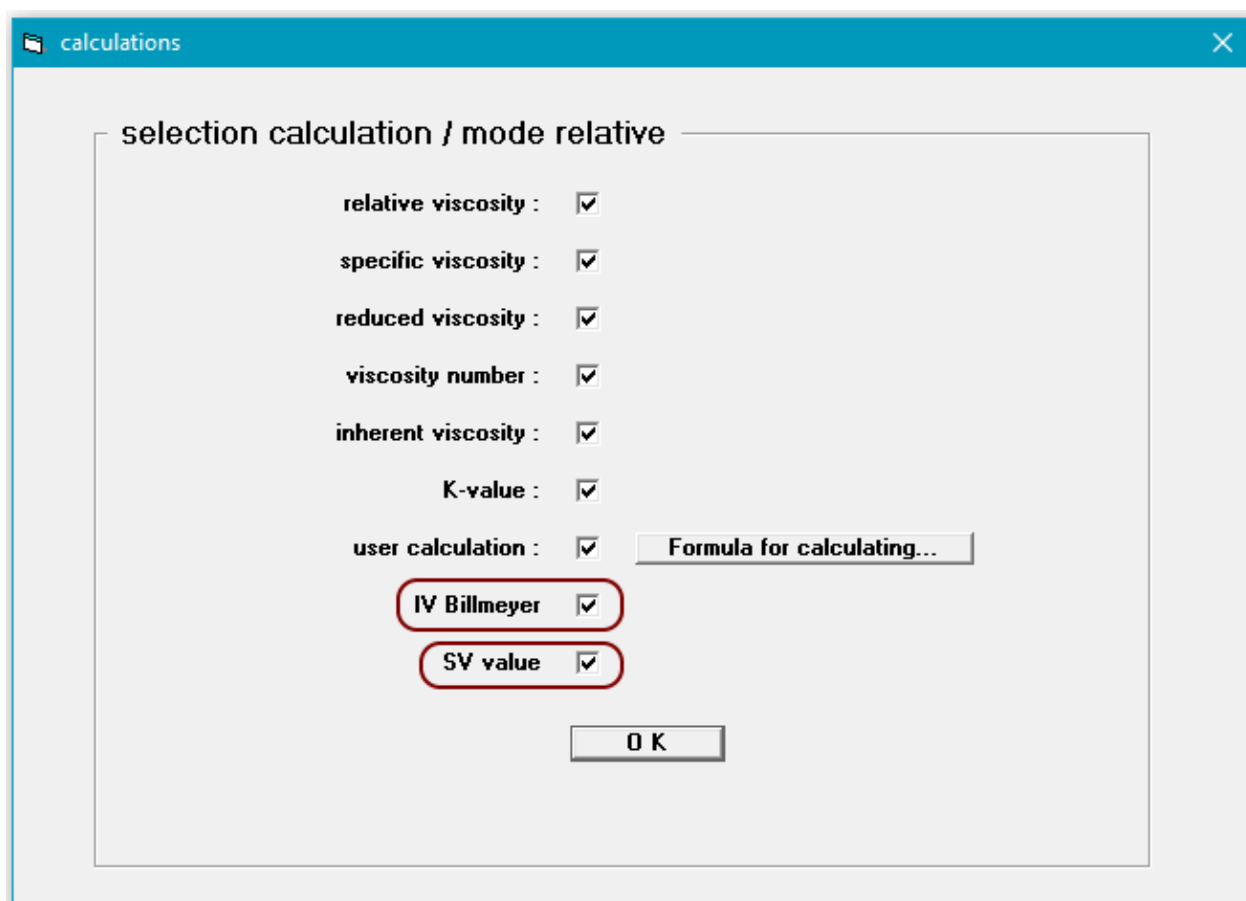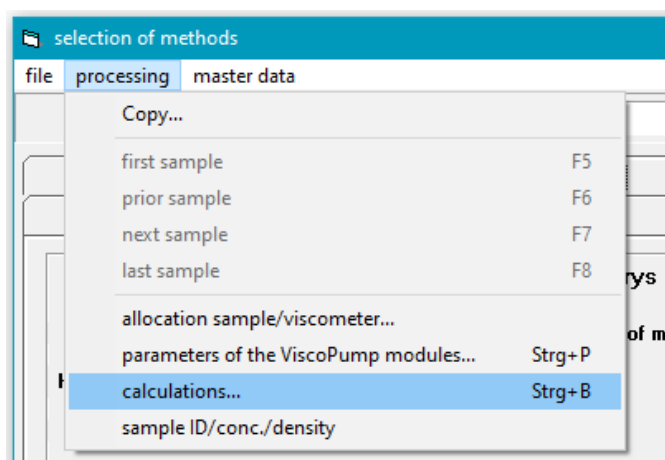
For the possible calculations in «mode relative», two additional calculations are available from version 19/09 onwards:


The results are also output in the measurement reports.

15 Prioritization of repositioned samples

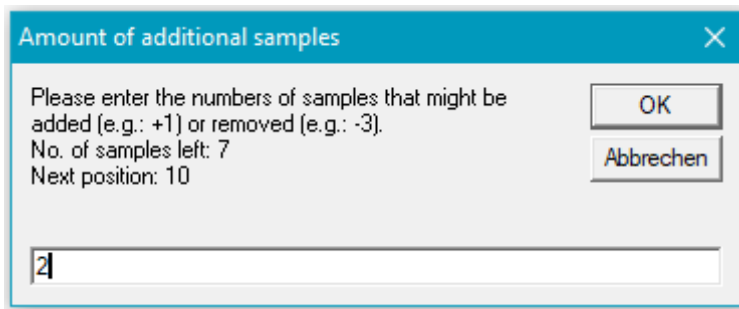
As of version 19/09, samples that can be freely assigned can be prioritized during the pause of a measurement series:

A screenshot of a software dialog box titled "Amount of additional samples". The dialog has a blue header bar with a close button (X) in the top right corner. The main area contains the text: "Please enter the numbers of samples that might be added (e.g.: +1) or removed (e.g.: -3).", "No. of samples left: 7", and "Next position: 10". On the right side, there are two buttons: "OK" and "Abbrechen". At the bottom, there is a text input field containing the number "2".

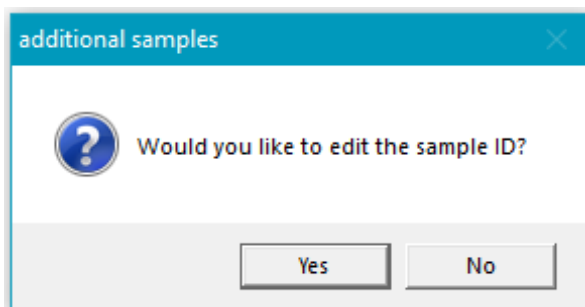A screenshot of a software dialog box titled "additional samples". The dialog has a blue header bar with a close button (X) in the top right corner. The main area contains a question mark icon and the text: "Would you like to edit the sample ID?". At the bottom, there are two buttons: "Yes" and "No".

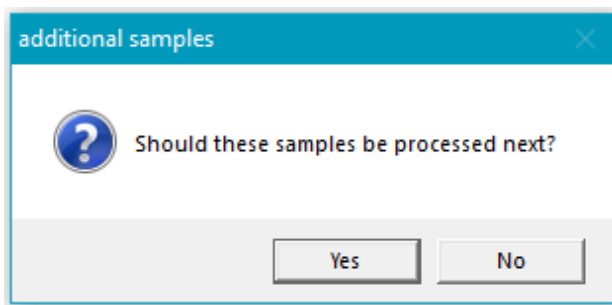
A screenshot of a software dialog box titled "additional samples". The dialog has a blue header bar with a close button (X) in the top right corner. The main area contains a question mark icon and the text: "Should these samples be processed next?". At the bottom, there are two buttons: "Yes" and "No".

With a click on «Yes», these readjusted samples are brought forward in the processing and measured next.


16 Dynamic sample list during measurements

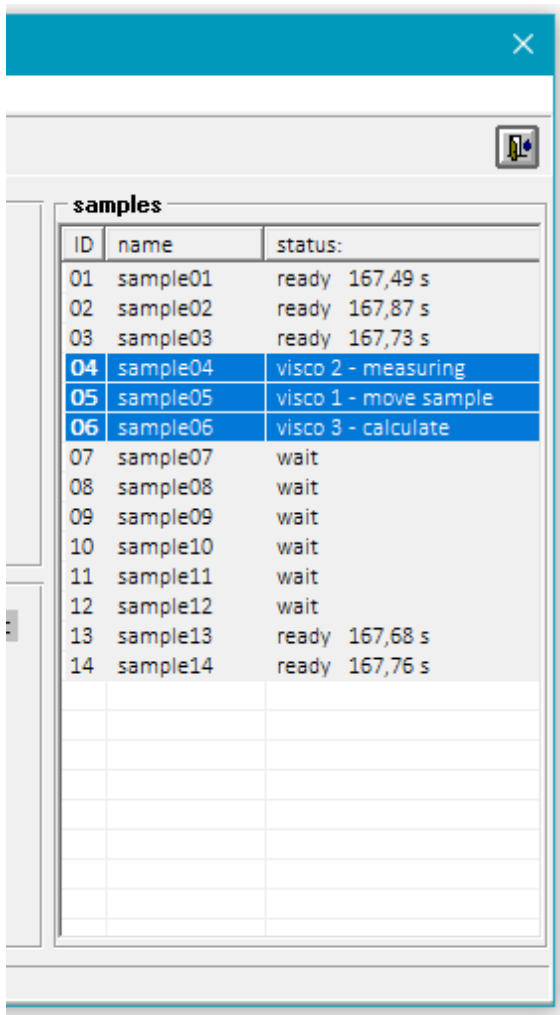From version 19/09 on, a dynamic sample list is displayed. This list shows the current status of each sample involved in the measurement series. For the measured samples, the result value is displayed.

ID	name	status:
01	sample01	ready 167,49 s
02	sample02	ready 167,87 s
03	sample03	ready 167,73 s
04	sample04	visco 2 - measuring
05	sample05	visco 1 - move sample
06	sample06	visco 3 - calculate
07	sample07	wait
08	sample08	wait
09	sample09	wait
10	sample10	wait
11	sample11	wait
12	sample12	wait
13	sample13	ready 167,68 s
14	sample14	ready 167,76 s

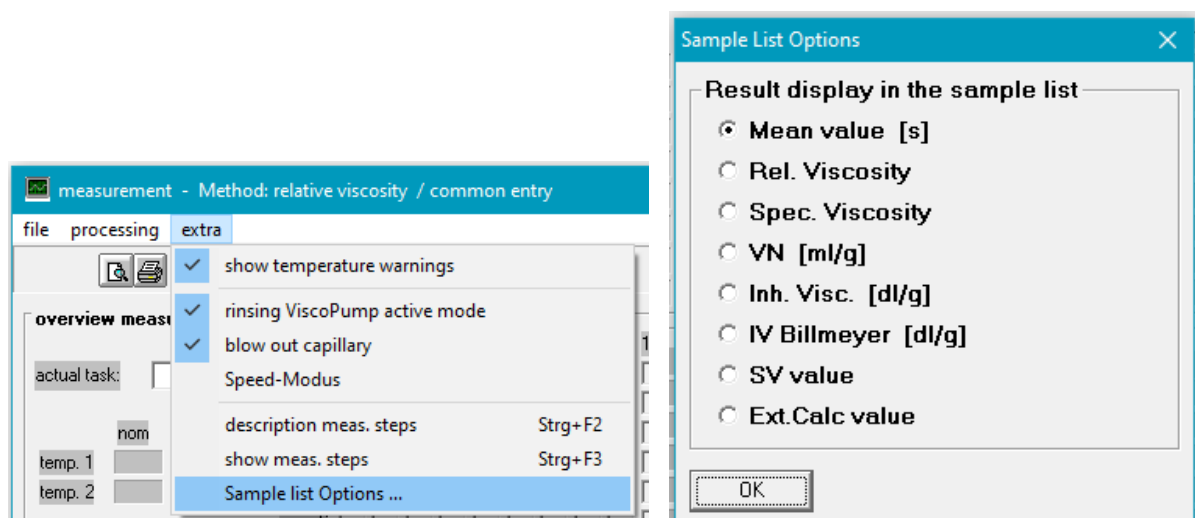The result value to be displayed can be selected with the menu item **«Extras»** → «Sample list options ...»:


17 t_0 editor and multiple sample plates

As of version 22/04, it is optionally possible not only to store one t_0 value for each viscometer, but also to enter up to 10 different values for each solvent used.

Each stored t_0 -flowtime is given a unique name with which it can be assigned to the individual samples in the sample list.

The use of the t_0 -flowtimes list can be activated in the "master data viscosimeter" dialog box with the menu option "processing" => "Use multiple t_0 -flowtimes".

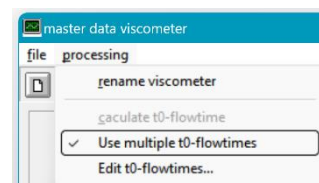By means of the then available menu option "processing" => "Edit t_0 -flowtimes..." the individual t_0 values per connected viscometer can be entered.

t0-flowtime editor

Stored t_0 -flowtime values per connected viscometer

	Sample solved with:	Pos. 1	Pos. 2
1	Cresol neutral	43,23	44,12
2	Cresol acid	41,52	41,65
3			
4			
5			
6			
7			
8			
9			
10			

close

In the sample list, the corresponding t_0 value can be selected for each sample.

sample ID/conc./density (Profile: 'Default')

Sample Tray A [A] [B] [C] [D]

sample 2 of 16

ID: Probe 2

concentration: 1,000 g/dl ☐ valid for whole list

density: 1 g/ml

T0 - Variante: Cresol neutral 1

ID	name/designation	concentration	dimension	density	T0
1	Probe 1	1,000	g/dl	1	2
2	Probe 2	1,000	g/dl	1	1
3	Probe 3	1,000	g/dl	1	0
4	Probe 4	1,000	g/dl	1	0

There are now 4 independent sample lists (A, B, C, and D) available for each profile. This means that another sample list can be prepared for the next series of measurements while one series is being processed. This is particularly interesting if several sample trays are available.

18 Change log

AVS® Pro 22/04 (Build 2.2.0.02)

Fixed:

- If sample 56 was used in individual sample assignment, an error occurred during processing.

AVS® Pro 22/04 (Build 2.2.0.01)

New:

- There are now 4 independent sample lists (A, B, C, and D) to choose from for each profile. This allows another sample list to be prepared for the next series of measurements while one series of prepared for the next series of measurements. This is particularly interesting if several sample trays are available. See video Multi-ProbList.mp4 in the Info folder of the installation CD.
- Optionally, a T₀ editor for the use of different t₀ blank values, depending on the type of sample solvent, can be used. See video t0-editor.mp4 in the Info folder of the installation CD.
- LIMS export files are now kept for 10 days and can be resent during this time.
- (Only AVS® Pro III) Supplement to the rinsing function Vacuum drying: While the dosing unit is being dried in the vacuum port, the visco is also aspirated. The visco suction already starts with the previously executed "Lösem.2 => Waste" rinsing.

Improvement:

- Revised sample tray travel protection concept. (min. MCM firmware 20220308 necessary; only for AVS® Pro II & III)
- (AVS®Pro II & III) MCM firmware vers. 20210717 with monitoring of the Z-axis drive.

Fixed:

- ViscoPump paramer ramp value limited to max. 50%. (previously 100%)
- Sample assignment for additional samples with and without prioritization revised.
- Error If, when using the PreFill option, during the prefill of the dosing device Pause was activated, the Z-axis did not move upwards.
- Error in solvent rinses with pre-rinses.

AVS® Pro 20/09 (Build 2.1.0.31)**New:**

- The LIMS-Output option has been extended:
Data transfer to the LIMS system either via FTP or via SMB to a [network] folder.
There are 3 different data formats available for the file structure.
Other formats can be installed on request
- The LIMS settings are managed separately in each profile.
- The "AVS® Pro Protocol" add-on application can export the result list to an Excel file.

Improvement:

- If, when emptying the dosing unit, the dosing piston has not reached the bottom of the dosing unit, a message will now be issued stating that the withdrawal speed is probably set too high.
- Monitoring of the Z-axis built in. (min. MCM firmware 20200630 necessary)

Change:

- Range of possible COM ports extended to 1 to 16.

Fixed:

- Generation of missing Extcalc.ini files corrected.
- Machine standstill due to unrecognized end of the rinsing process.
- Error at the end of the measurement series if the option "Park in waste bottle" is set.

AVS® Pro 19/09 (Build 2.1.0.16)**New:**

- During a measurement series, a list of all samples involved with the respective processing status is now displayed.
- Reordered samples with free assignment can now be prioritized.
- Automatic check of MCM firmware version for up-to-dateness.
- New calculations "IV Billmeyer" and "SV-Value" integrated.
- The logged-in user is now displayed on the measurement logs.
- Optional "Solvent rinsing after the last sample" for each viscometer.

Improvement:

- No "0" is entered in the measurement protocols for measurements and calculations that were not performed, but nothing.

Change:

- If more than one profile exists, every user can now use the "Profile management" menu option to change the profile. However, only the administrator can still create or delete profiles.
- If the option "Park in waste bottle" is set, the system waits until the journey to the waste bottle is completed before releasing the measurement dialog again.

Fixed:

- When aborting a series of measurements with a 56 sample rack, the safety emptying was carried out at screw speed.
- Unintentional preassignment of the input type with "Determine blank value".
- When changing the measuring method, the heading in the Measuring dialog box was not adjusted.

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Xylem Analytics Germany Sales GmbH & Co.KG