**GEBRAUCHSANLEITUNG**  
Originalversion  
**OPERATING MANUAL**  
**MODE D'EMPLOI**  
**MANUAL DE INSTRUCCIONES**

---

# TitroLine<sup>®</sup> 7500 KF *trace*

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## **Operating Manual ..... Page 83 ... 162**

### **Important notes:**

The operating manual is part of the product. Before initial operation, please carefully read and observe the operating manual and keep it. For safety reasons the product may only be used for the purposes described in these present operating manual. Please also consider the operating manuals for the devices to be connected.

All specifications in this operating manual are guidance values which are valid at the time of printing. However, for technical or commercial reasons or in the necessity to comply with the statutory stipulations of various countries, the manufacturer may perform additions to the product without changing the described properties. A potentially more recent version of this manual is available on our internet website. The German version is the original version and binding in all specifications!

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## TABLE OF CONTENT

<b>1</b>	<b>Technical Specifications of the Titrator TitroLine® 7500 KF trace</b>	<b>85</b>
1.1	Notes to the operating manual	85
1.2	Intended Use	85
1.3	Technical Specifications	86
1.3.1	Titrationstand TM 235 KF	86
1.3.2	Titrationstand TM 235 KF	88
1.4	Warning and safety information	89
1.4.1	General	89
1.4.2	Chemical and biological safety	90
1.4.3	Flammable liquids	90
<b>2</b>	<b>Installation and Commissioning</b>	<b>91</b>
2.1	Unpacking and setting up	91
2.2	Back panel of the titrator TitroLine® 7500 KF trace	92
2.3	Connection and installation of the titrator and the magnetic stirrer TM235/TM235 KF	93
2.4	Setting the language	94
2.5	Installation with magnetic stirrer TM 235 (module 1 and 3)	95
2.6	Installation with magnetic stirrer/pump TM 235 KF (module 2 and 4)	97
2.6.1	Working with the titration stand	98
2.6.2	Troubles	98
2.7	Filling the Titration Vessel	99
2.8	Switch on Device, First Conditioning	99
2.9	Combination with Accessories and Additional Devices	99
2.9.1	Connecting a printer	99
2.9.2	Connecting a USB device	99
2.9.3	Connection of analytical balances	99
<b>3</b>	<b>Working with the Titrator TitroLine® 7500 KF trace</b>	<b>100</b>
3.1	Front Keyboard	100
3.2	Display	100
3.3	External PC Keyboard	101
3.4	Menu Structure	102
3.5	Main Menu	104
3.5.1	Starting a Titration	104
<b>4</b>	<b>Method parameters</b>	<b>107</b>
4.1	Method editing and new method	107
4.2	Default method	108
4.3	Copy Method	108
4.4	Delete Method	109
4.5	Print method	109
4.6	Change Method Parameters	110
4.6.1	Result	110
4.6.2	Formula Editor	117
4.6.3	Titration parameters	123
4.6.4	Sample identification	127
4.6.5	Documentation	128
<b>5</b>	<b>System settings</b>	<b>129</b>
5.1	Global Memory	129
5.2	RS-232 Settings	129
5.3	Date and Time	132
5.4	Password	133
5.4.1	Creation of the first Administrator	133
5.4.2	Creation of additional users	137
5.4.3	Predefined rights and definable rights	138
5.4.4	Delete of users	142
5.5	RESET	143
5.6	Printer	144
5.7	Device Information	144
5.8	System Tones	144
5.9	Data exchange	145
5.10	Software Update	147

<b>6</b>	<b>Network settings</b> .....	<b>149</b>
6.1	General.....	149
6.2	Setup a shared directory .....	150
<b>7</b>	<b>Communication via RS-232 and USB-B interface</b> .....	<b>152</b>
7.1	General Information.....	152
7.2	Chaining multiple devices - "Daisy Chain Concept".....	152
7.3	Instruction Set for RS-Communication.....	152
<b>8</b>	<b>Connection of Analytical Balances and Printers</b> .....	<b>154</b>
8.1	Connection of Analytical Balances.....	154
8.2	Balance data editor .....	155
8.3	Printers .....	156
8.4	Automatic stirrer control .....	157
8.4.1	General.....	157
8.4.2	Basic setting in the system menu.....	157
8.4.3	Set the stirring speed in the method .....	158
8.5	Autosampler .....	159
8.5.1	Connection of sample changer TW 7400.....	159
8.6	Using software TitriSoft .....	159
8.6.1	General.....	159
8.6.2	TitriSoft 3.15 or higher.....	159
<b>9</b>	<b>Maintenance and Care of the Titrator</b> .....	<b>160</b>
<b>10</b>	<b>Guarantee</b> .....	<b>161</b>
<b>11</b>	<b>Storage and transportation</b> .....	<b>161</b>
<b>12</b>	<b>Recycling and Disposal</b> .....	<b>161</b>
<b>13</b>	<b>EC – Declaration of Conformity</b> .....	<b>162</b>

---

## Copyright

© 2022, Xylem Analytics Germany GmbH

Reprinting - even as excerpts - is only allowed with the explicit written authorization.  
Germany. Printed in Germany.

# 1 Technical Specifications of the Titrator TitroLine<sup>®</sup> 7500 KF *trace*

## 1.1 Notes to the operating manual

The provided operating manual will allow you the proper and safe handling of the product. For maximum security, observe the safety and warning instructions in the operating manual!

-  Warning of a general danger:  
Non-compliance results (can result) in injury or material damage.
-  Important information for device use.
-  Refers to another part of the operating manual.

The menu screens shown in this operating manual serve as an example and may differ from what you see!

## 1.2 Intended Use

The TitroLine<sup>®</sup> 7500 KF *trace* is a coulometric titrator and suitable for coulometric KF- titrations or coulometric determination of the bromine number with a maximum of 50 memorisable methods.

The TitroLine<sup>®</sup> 7500 KF *trace* can be used as stand-alone instrument or in combination with a heating oven.

### Solutions to be used:

All common and modern Karl-Fischer reagents can be used for the coulometry process.

-  The device is not intended for use with potentially biohazardous substances.

 **Do not use the device in hazardous locations!**

 **General:**

The safety guidelines that are applicable to the handling of chemicals have to be observed under all circumstances. This applies in particular to inflammable and/or etching liquids.

## 1.3 Technical Specifications

### 1.3.1 Titrator TitroLine® 7500 KF *trace*

Translation of the legally binding German version

(Release: 21. February 2020)

EMC compatibility according to the Council Directive: 2014/30/EU;  
 applied harmonized standards: EN 61326-1  
 Low-voltage directive according to the Council Directive 2014/35/EU;  
 Testing basis EN 61 010-1: for laboratory equipment  
 RoHS Council Directive 2011/65/EU  
 FCC Part 15B and ICES 003

**Country of origin:** Germany, Made in Germany

**The following solvents/titration reagents are allowed to be used:**

- All common and modern Karl-Fischer reagents can be used for the coulometry process.

**Measuring input ( $\mu\text{A}$ ):**

Karl Fischer (Dead-Stop) connector ( $\mu\text{A}$ ) for double platinum electrode.  
 Polarisation voltage variably adjustable from 40 ... 220 mV.  
 Connector: 2 x 4 mm – sockets.

Measurement range I [ $\mu\text{A}$ ]	Display resolution	Measurement accuracy* without sensor probe
0 ... 100	0.1	+/- 0.2 $\mu\text{A}$ $\pm$ 1 Digit

**Display:** 3.5 inches -1/4 VGA TFT display with 320 x 240 pixels.

**Input:** Measurement input  $\mu\text{A}$ :

(Dead-Stop-) connector for double platinum electrode (Connection sockets: 2 x 4 mm)  
 Input for generator electrode (Connection sockets: 2 x 4 mm), colours green (grey) and black

**Power supply:** by external multi-range power supply from 100 – 240 V, 50/60 Hz  
 Input voltage: 12 Volt DC, 2500 mA  
 Power consumption 30 W  
 Corresponds to protection class III:  
 Protection class for dust and humidity IP 50 according to DIN 40 050

**⚠ Only use the power supply TZ 1853 or a power supply approved by the manufacturer!**

**RS-232-C Interface:** separated galvanically through photocoupler, Daisy Chain function available

Data bits: adjustable, 7 or 8 Bit (default: 8 Bit)  
 Stop bit: adjustable, 1 or 2 Bit (default: 1 Bit)  
 Start bit: static 1 Bit  
 Parity: adjustable: even / odd / **none**  
 Baud rate: adjustable: 1200, 2400, **4800**, 9600, 19200 (Default 4800 baud)  
 Address: adjustable, (0 to 15, default: 01)

RS-232-1 for computer, input Daisy Chain

RS-232-2 devices of SI Analytics®:  
 - Titrator TitroLine® 7000 / 7500 / 7500 KF / 7750 / 7800  
 - Sample Changer TW alpha plus, TW 7400  
 - Piston burette TITRONIC® 300 and 500, TITRONIC® 110 *plus*, TITRONIC® *universal*,  
 - Balances of the types Mettler, Sartorius, Kern, Ohaus, (for more, please contact us)  
 - Exit Daisy-Chain

**USB Interface:** 2 x USB-type-A and 1 x USB-type-B

USB-type A for connecting of USB keyboard, - printer, - manual controller, - data media (e.g. USB stick) and USB-Hub

USB-type B for connecting a PC

\* The measurement uncertainty of the sensor probe has to be taken into account as well.

<b>Ethernet Interface:</b>	for connecting a local network (LAN)
<b>Stirrer/pump:</b>	12V DC out, 500 mA power supply for stirrer TM 235 and KF titration stand TM 235 KF
<b>Housing:</b>	
Material:	Polypropylene
Front keyboard:	polyester coated
Dimensions:	15.3 x 45 x 29.6 cm (W x H x D)
Weight:	approx. 2.3 kg for basic unit without stirrer TM 235 or KF titration stand TM 235 KF

**Ambient conditions:**

 **Do not use the device in hazardous locations!**

Climate:	Ambient temperature: + 10 ... + 40 °C for operation and storage Humidity according to EN 61 010, Part 1: Max. relative humidity 80 % for temperatures up to 31 °C, linear decrease down to 50 % relative humidity at a temperature of 40 °C
Altitude:	Device: No restrictions Power supply: up to 5000 m
Pollution degree:	Pollution degree IP 20, indoor use only.

**Software**

Measuring range:	10 µg – 100 mg / 1 ppm – 5 % (recommended)
Measuring speed:	Max. 1.5 mg/min
Number methods:	50
Conditioning:	Automatically with drift determination
End point criteria:	Drift, stop drift tolerance, min. and maximum titration time
Auto start:	Yes, after sample allowance
Statistics:	Mean value, standard deviation and relative standard deviation
Curve printout:	Measuring unit/time
Documentation:	GLP compliant on printer or in PDF format (USB- memory stick)

### 1.3.2 Titrationstand TM 235 KF

Translation of the legally binding German version

(Release: 21. February 2020)

In connection with the titrator TitroLine® 7500 KF *trace*

EMC compatibility according to the Council Directive: 2014/30/EU;  
 applied harmonized standards: EN 61326-1  
 Low-voltage directive according to the Council Directive 2014/35/EU;  
 Testing basis EN 61 010-1: for laboratory equipment  
 RoHS Council Directive 2011/65/EU  
 FCC Part 15B and ICES 003

**Country of origin:** Germany, Made in Germany

**Pump:** Free volume flow- air: flow rate 2.25 l / min  
 Delivery pressure max.: 1,5 bar  
 Flow rate liquid medium: approx. 0,8 l / min

**Stirring speed:** 50 ... 1000 U/min

**Hoses:** PVC- hose (outer diameter 6 x 1 mm)  
 PTFE- hose (outer diameter 4 x 0.5 mm)

#### Connections

Power supply: Low voltage input 12 V / – on the backside of titration stand  
 Plug connection: plug for low voltage connection – phone jack-,  
 Positive pole at pin contact, inside contact  $\varnothing = 2.1$  mm, USA/Japan,  
 Power supply via titrator TitroLine® 7500 KF *trace*

#### Housing:

Material: Polypropylene, polyester coated  
 Dimensions: 80 x 130 x 250 mm (W x H x D), height without stand  
 Weight: 1.0 kg

#### Ambient conditions:

**Do not use the device in hazardous locations!**

Climate: Ambient temperature: + 10 ... + 40 °C for operation and storage  
 Humidity according to EN 61 010, Part 1:  
 Max. relative humidity 80 % for temperatures up to 31 °C,  
 linear decrease down to 50 % relative humidity at a temperature of 40 °C

Altitude: Device: No restrictions  
 Power supply: up to 5000 m

Pollution degree: Pollution degree IP 20, indoor use only.

## 1.4 Warning and safety information

### 1.4.1 General

The device corresponds to protection class III.

It was manufactured and tested according to DIN EN 61 010, Part 1, "**Protective Measures for electronic measurement devices**" and control devices and has left the factory in an impeccable condition as concerns safety technology. In order to maintain this condition and to ensure safe operation, the user should observe the notes and warning information contained in the present operating instructions. Development and production is done within a system which meets the requirements laid down in the DIN EN ISO 9001 standard.

 For reasons of safety, the device must only be used for the range of application described in the present operating manual. Nonobservance of the intended proper use of the device may result in personal injury or damage to property.

 For reasons of safety, the device and the power supply must be opened by authorised persons only; this means, for instance, that work on electrical equipment must only be performed by qualified specialists. **In case of nonobservance of these provisions the titrator and the power supply may constitute a danger: electrical accidents of persons or fire hazard!** Moreover, in the case of unauthorised intervention in the titrator or the power supply, as well as in the case of negligently or deliberately caused damage, the warranty will become void.

 Prior to switching the device on it has to be ensured that the operating voltage matches the mains voltage. The operating voltage is indicated on the specification plate (underside of the device and backside of the power supply). **Nonobservance of this provision may result in damage to the titrator and the power supply, or in personal injury or damage to property!**

 **If it has to be assumed that safe operation is impossible, the device has to be put out of operation and secured against inadvertent putting to operation.** In this case please switch the device off, pull plug of the mains cable out of the power supply, and remove the device from the place of work.

Examples for the assumption that a safe operation is no longer possible,

- if the package is damaged,
- if the device shows visible damages,
- if the power supply shows visible damages,
- if the device does not function properly,
- if liquid has penetrated into the casing.
- if the unit has been altered technologically or if unauthorized personnel tried or succeeded to open the device as attempt to repair it.

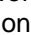
In case that the user operates such a device, all thereof resulting risks are on the user!

 The device must not be stored or operated in humid rooms.

 **The relevant regulations regarding the handling of the substances used have to be observed:** The Decree on Hazardous Matters, the Chemicals Act, and the rules and information of the chemicals trade. On the part of the user it has to be ensured that the persons entrusted with the use of the unit are experts in the handling of substances used in the environment or that they are supervised by specialized persons, respectively.

 For all work with chemicals: **Always wear protective glasses!** Please observe the memorandums of the employer's liability insurance associations and the safety data sheets of the manufacturers.

 The device is equipped with integrated circuits (EPROMs). X rays or other high energy radiation may penetrate through the device's casing and delete the program.

 For working with liquids, not being common titration solvents, especially the chemical resistance of the construction materials of the device have to be considered (see  1.3 Technical Specifications).

 For the use of liquids with high vapour pressure or (mixture of) substances not being mentioned in  1.3 Technical Specifications) as allowed substances, the safe and proper operation of the device has to be guaranteed by the user. When the piston moves upwards within the cylinder, a microfilm of dosing liquid or titration solution will always remain adhered to the inner wall of the cylinder, but this has no influence on the dosing accuracy. This small residue of liquid, however, may evaporate and thus penetrate into the zone underneath the piston, and if non-admitted liquids are being used, the materials of the may be dissolved or corroded (see  9 Maintenance and Care of the Titrator).

### 1.4.2 Chemical and biological safety

 The device is not intended for use with potentially biohazardous substances.

 **The relevant regulations regarding the handling of the substances used have to be observed:** The Decree on Hazardous Matters, the Chemicals Act, and the rules and information of the chemicals trade. On the part of the user it has to be ensured that the persons entrusted with the use of the unit are experts in the handling of substances used in the devices or that they are supervised by specialized persons, respectively.

 When using biohazardous substances, the regulations for handling the substances used must be observed. In such cases, the use is the sole responsibility of the user.

 For all work with chemicals: **Always wear protective glasses!** Please observe the memorandums of the employer's liability insurance associations and the safety data sheets of the manufacturers.

 Dispose of all used solutions in accordance with national regulations and laws. Select the type of protective equipment according to the concentration and quantity of the hazardous substance at the respective workplace.

### 1.4.3 Flammable liquids

When handling flammable liquids, make sure that there is no naked flame in the vicinity of the equipment. Adequate ventilation must be provided. Only small quantities of flammable liquids should be kept in the workplace.

 When working with liquids that do not correspond to common reagents, particular attention must be paid to the chemical resistance of the materials of the devices (vgl.  **Fehler! Verweisquelle konnte nicht gefunden werden.** Technical Specifications).

## 2 Installation and Commissioning

### 2.1 Unpacking and setting up

The device has been put together especially for you (basic unit + corresponding modules and accessories), there may be differences with respect to the delivery and the accessories described in this chapter. The scope of delivery, please refer to the attached packing list. For any questions please contact us directly (see backside of this operating manual).


The device itself as well as all related accessory and peripheral parts have been carefully checked at the factory to ensure their correct function and size. Please ensure that the small accessories are also removed in full from the packaging.

The device may be placed on any flat surface.

Scope of delivery:

- a) Titrator TitroLine® 7500
  - TitroLine® 7500
  - Keyboard TZ 3835
  - Power supply TZ 1853 (100 V ... 240 V) incl. some primary adapter
  - Connection cable for stirrer/pump TZ 1577
  - Stand rod TZ 1748 (10 mm x 280 mm)
  - Piston extraction tool TZ 3813
  
- b) KF accessories
  - KF starter kit TZ 1789 with molecular sieve, glass wool and a set of syringes with needles
  - Electrode KF 1150
  
  - **For module 1 and 3 only:**
    - Magnetic stirrer TM 235 and titration vessel TZ 1751
  - **For module 2 and 4 only:**
    - KF titration stand (pump and stirrer) TM 235 KF with waste (1 L clear glass), solvent (1 L amber glass) and bottle for the desiccant (100 ml) and with all PTFE and PVC hoses
    - Titration vessel TZ 1754
  - **For module 1 and 2:**
    - Generator electrode TZ 1754 without diaphragm
  - **For module 3 and 4:**
    - Generator electrode TZ 1753 with diaphragm
  - **For module 5:**
    - Magnetic stirrer TM 235 and titration vessel TZ 1755

## 2.2 Back panel of the titrator TitroLine® 7500 KF trace

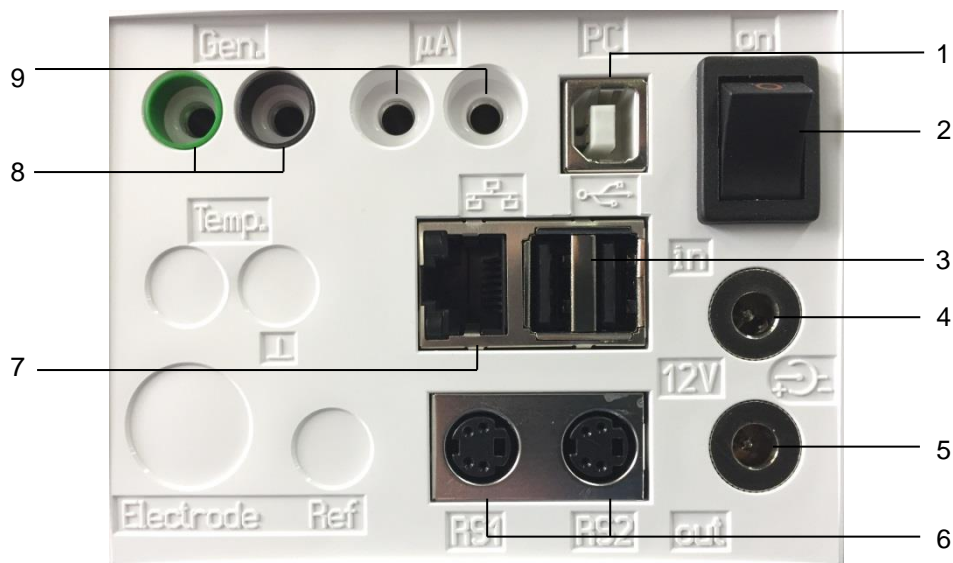**Fig. 1**

The TitroLine® 7500 KF *trace* is equipped with the following connections:

- 1) USB-B interface for connection to a PC
- 2) On/Off switch
- 3) Two USB-A interfaces for connecting USB devices
- 4) Socket "in": Connection of the external power supply TZ 1853
- 5) Socket "out": Connection of the TM 235 / TM 235 KF magnetic stirrer
- 6) Two RS232 ports, 4-channel (Mini-DIN):  
     RS1 for connection to the PC  
     RS2 for connection of a weighing balance and other devices from SI Analytics®
- 7) Ethernet Interface (LAN)
- 8) Input for a generator electrode, coloured green (grey) and black
- 9) µA measurement input for the connection of double platinum electrodes

### 2.3 Connection and installation of the titrator and the magnetic stirrer TM235/TM235 KF

The low voltage cable of the power supply TZ 1853 has to be plugged in to the 12 V socket "in", on the back panel of the titrator. (Fig. 2). Then plug the power supply into the plug socket.

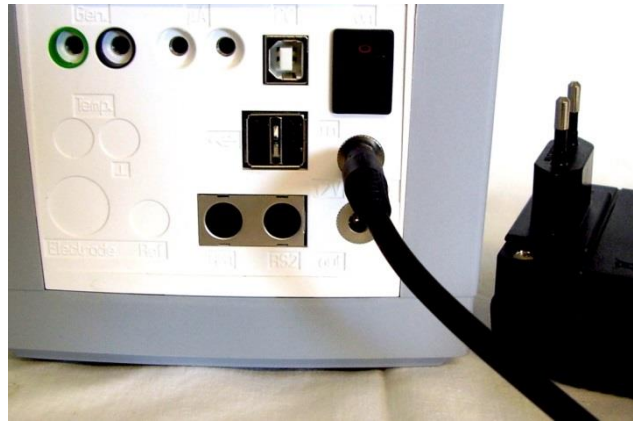
Fig. 2

Fig. 3


**⚠** Place the power supply easily accessible in order to be able to remove the titrator anytime easily from the power circuit.

Place the TM 235/ TM 235 magnetic stirrer to the right of the titrator (Fig. 3) and connect to the 12V out-socket in the rear panel of the piston burette by using the TZ 1577 connection cable.

## 2.4 Setting the language


The ex-factory default language setting is English.

After the device is switched on and the start-up process is complete, the main menu appears (Fig. 4).

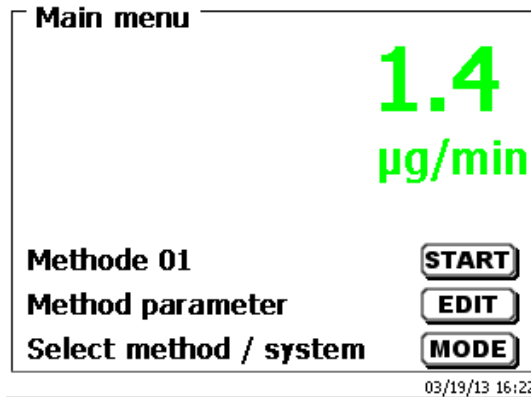
**Fig. 4**

Using <SYS> or <MODE>, you navigate to the system settings («System settings»). The very first menu is to be used for setting the language (Fig. 5).

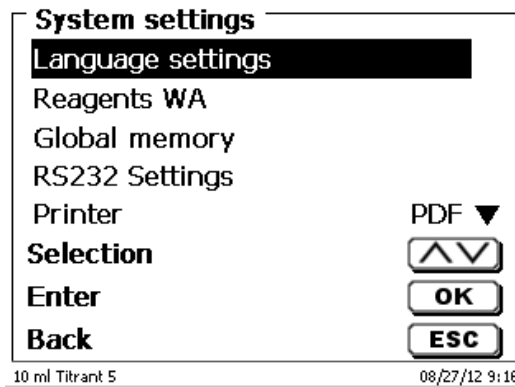**Fig. 5**

Use <ENTER>/<OK> to call the menu.  
Select the language using the arrow keys <↑↓>.  
Confirm with <ENTER>/<OK>.

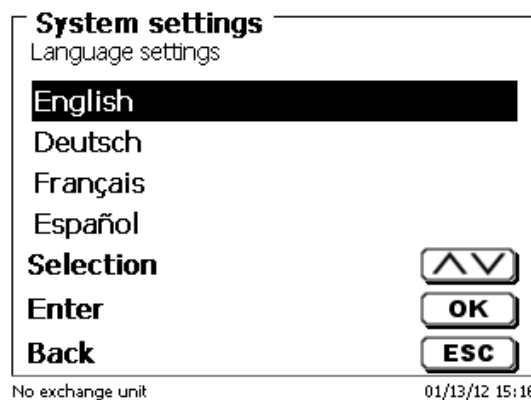**Fig. 6**

The selected language will appear immediately (Fig. 6).  
Pressing <ESC> twice will return the user to the main menu.

## 2.5 Installation with magnetic stirrer TM 235 (module 1 and 3)

Place the TM 235 titration stand to the right of the titrator and connect to the 12V out-socket in the rear panel of the piston burette by using the TZ 1577 connection cable (Fig. 7). Then screw the stand rod into the thread of the titration stand (Fig. 8) and mount the titration clamp (Fig. 9).

Fig. 7

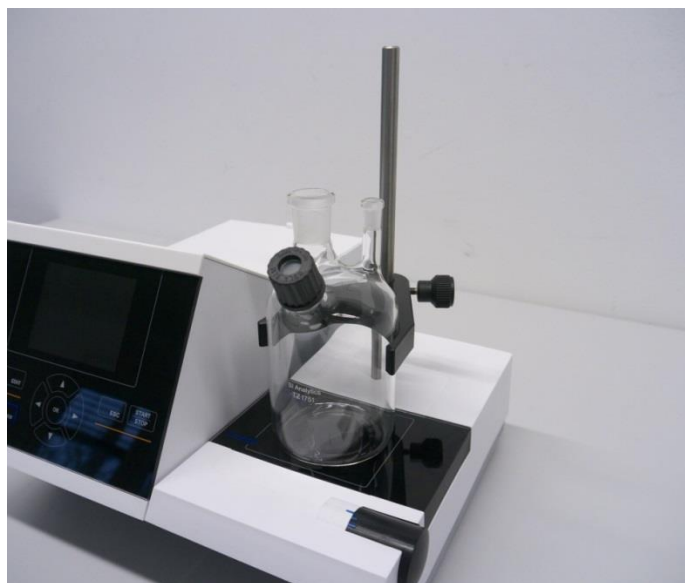
Fig. 8


Fig. 9

Clamp the titration vessel to the retaining clip. Fix the titration vessel in such a way that the bottom of the titration vessel stands directly at the upper surface of the magnetic stirrer (Fig. 10).

**Fig. 10**

Insert the electrode KF 1150 and the generator electrode (TZ 1752 or TZ 1753) into the NS 7 and NS 19 openings of the titration vessel provided for this purpose. Connect the fixed cable of the KF 1150 indicator electrode to the  $\mu\text{A}$  connection. The cable for the generator electrode LB 04 NN has a green (grey) and a black plug and is connected to the sockets of the generator measuring input marked in colour (Fig. 11).

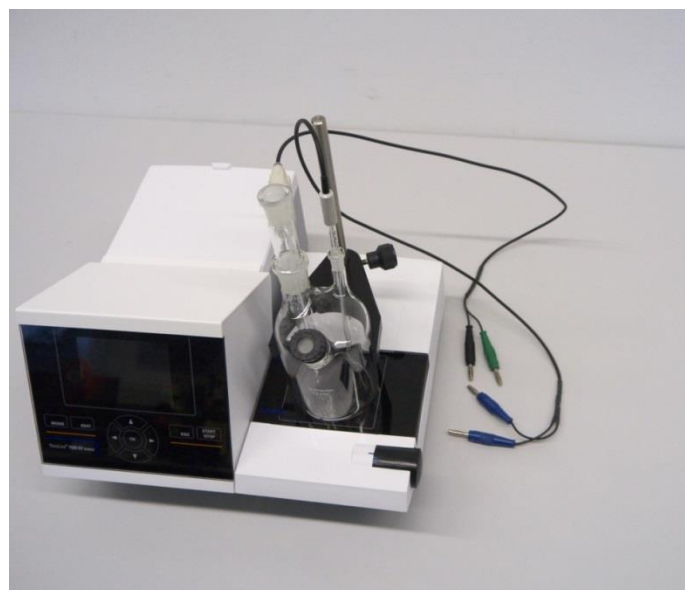**Fig. 11**


## 2.6 Installation with magnetic stirrer/pump TM 235 KF (module 2 and 4)

Place the TM 235 KF titration stand to the right of the titrator and connect to the 12V out-socket in the rear panel of the piston burette by using the TZ 1577 connection cable (see Fig. 7). Then screw the stand rod into the thread of the titration stand (see Fig. 8) and mount the titration clamp (see Fig. 9). Clamp the titration vessel to the retaining clip. Fix the titration vessel in such a way that the bottom of the titration vessel stands directly at the upper surface of the magnetic stirrer (see Fig. 10).

Mount all white plastic adapters and grey screw caps on the dry, waste and solvent bottles. Connect the PVC and PTFE hoses (Fig. 12 - Fig. 14).

The PVC tubes are connected to the connectors at the back side of the TM 235 KF. The long PVC tube is used for the connection of the waste bottle. The two shorter PVC ones are used to connect the moisture bottle and the solvent bottle. The moisture bottle is connected to the right connector (view from above) of the TM 235 KF. The waste (clear) bottle is connected to the left connector (Fig. 12).

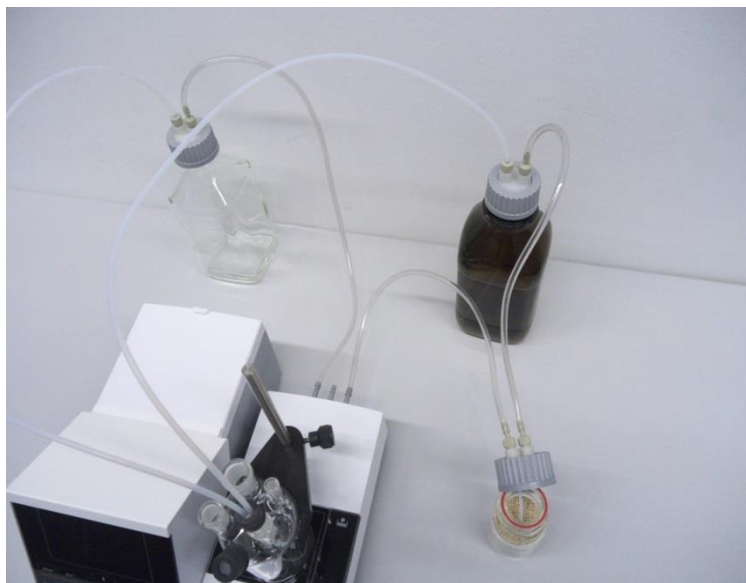
**Fig. 12**

Put the threaded pipe with the NS 14/23 core and the GL 14-thread in one of the NS-14.5-openings. Put both PTFE-tubes through both bore holes of the septum. The PTFE tube from the clear waste bottle is adjusted to the bottom of the titration vessel (tube 1). Please push one tube ("Tube 1") through to the bottom of the titration vessel, the other tube ("Tube 2") only about 1-2 cm (Fig. 13).

**Fig. 13**

Please put the other end of the PTFE tube that touches the bottom of the titration vessel ("tube 1"), through the opening on the cap of the clear square bottle (-> waste bottle). Put the other PTFE tube ("tube 2") through the opening of the cap of the brown reagent bottle. Adjust the dosage and the disposal tubes as depicted in Fig. 14. Then seize the screwing with the tubes on the bottles.

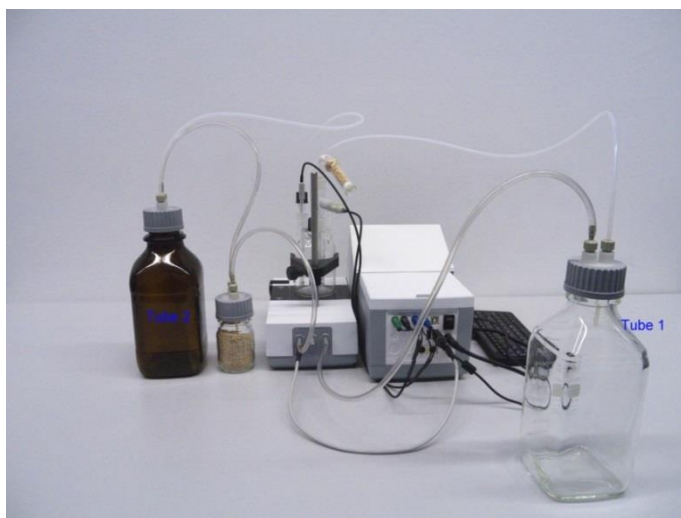**Fig. 14**

Put the indicator electrode KF 1150 and the generator electrode (TZ 1752 or TZ 1753) into the provided openings NS 7,5 and NS 19 and the electrode cables into the colour-coded socket. The indicator electrode has a fixed cable with 2 blue plugs. The cable LB 04 NN has a green and black plug and is connected to the provided port of the generator electrode.

The keyboard is connected to one of the USB-A ports.

**⚠** Place the power supply easily accessible in order to be able to remove the titrator anytime easily from the power circuit.

Fill the KF analyte into the brown 1 l reagent bottle. You can also connect the GL 45 adaptor directly to the analyte bottle (recommended) if it has a GL-45 -thread.

Fill the TZ 1505 glass drying tube with glass wool, then molecular sieve and then again with glass wool. Plug the drying tube onto the generator electrode.

### 2.6.1 Working with the titration stand

- Dosage: Pump the analyte into the titration vessel by pressing the front part of the rocker switch. As long as you press the button the dosing process proceeds.
- Siphon off: Siphon off the solution from the titration vessel by pressing the back part of the rocker switch. As long as you press the button it will be siphoned off.

**⚠ Please always check the fill level of the waste bottle before operating the pump.** Before you siphon off the titrating solution please make sure that the disposal bottle can absorb this amount of solution.

The built-in magnetic stirrer stirs the liquids in the titration vessel. You can adjust the stirring speed at the turning knob at the right upper side of the titration stand or magnetic stirrer.

### 2.6.2 Troubles

Problems might occur if the tubes are not connected properly or the pressure / low pressure system has a leakage. Then, after a few seconds operating time of the pump, no reagents will be conveyed anymore. When checking the tubes please observe that the bottle screwing and all adaptors are leak-proof. The same applies to the tube connections of the drying bottle.

**⚠ Buckled tubes cause incorrect dosages and also the siphon off process does not work properly. Please check the tubes and their connections for leak-tightness on a regular basis. Replace if necessary**

If the reagents continue to run in from the storage bottle after the actual dosage process has ended, position the bottle at lower level than the titration vessel. If the reagents continue to run in from the storage bottle between the dosage process and the siphon off process, please wait a few seconds between the changes.

**i** We recommend to remove the glass adaptor with the dosing and exhaust hose after adding the KF-analyte and to close the ground-in opening with the provided die glass stopper NS 14.5.

## 2.7 Filling the Titration Vessel

Fill the titration vessel approx. up to a half with analyte that are suitable for the respective application using a funnel (Module 1 and 3). You can even use the pump for modules 2 and 4. If you use a generator electrode with diaphragm, it is necessary to fill an ampoule with catholyte into the generator electrode. Here, please use a syringe with needle (Fig. 15).

Fig. 15

## 2.8 Switch on Device, First Conditioning

Set the stirring speed at the TM 235/TM 235 KF. Please do not forget the magnetic stirring bar! Now switch on the device. The mains switch is situated on the left at the back side of the TitroLine<sup>®</sup> 7500 KF *trace*: the device starts conditioning immediately. If the analyte is freshly filled in, the conditioning can last 10 - 20 minutes. In the case of the generator electrode with diaphragm this can also take more than one hour. The drift display is then > 1500 µg/min. When the drift has dropped below 10 µg/min, the TitroLine<sup>®</sup> 7500 KF *trace* is ready for the first measurements.

## 2.9 Combination with Accessories and Additional Devices

### 2.9.1 Connecting a printer

Printers with a USB interface are to be connected to one of the two USB-A interfaces.

**i** These printers **have to** feature HP PCL emulation (3, 3 enhanced, 5, 5e). So-called GDI printers cannot be used!

Alternatively the thermo-compact printer Seiko S445 can be connected.

### 2.9.2 Connecting a USB device

The following USB devices can be connected to the USB-A interfaces:

- PC-keyboard
- TZ 3880 manual controller
- Printer
- USB storage devices, e.g. USB sticks
- USB-Hub
- USB barcode scanners

### 2.9.3 Connection of analytical balances

Analytical balances are to be connected to the RS232-2 using an appropriate cable.

### 3 Working with the Titrator TitroLine® 7500 KF trace

#### 3.1 Front Keyboard

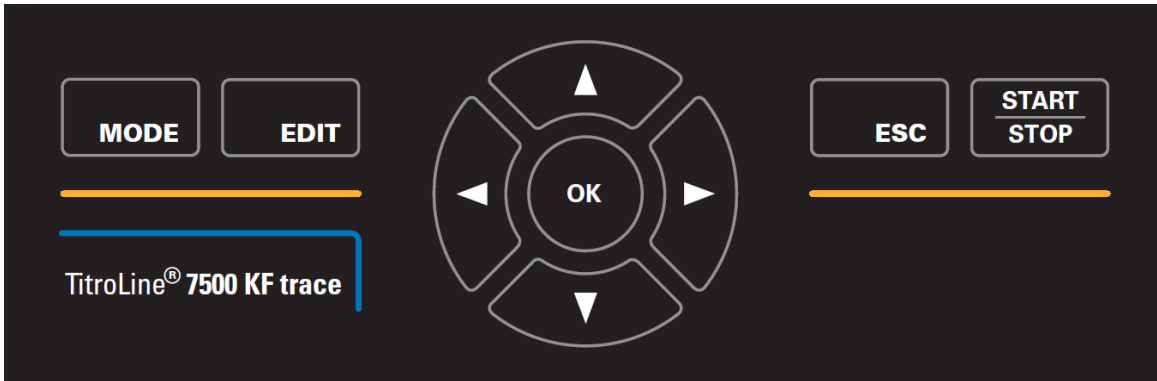
Fig. 16

**i** Apart from alphanumeric input (a-z, A-Z, 0-9) and a few other functions, almost all functions can be performed using the front keyboard (Fig. 16).

- <Mode>: Methods selection, rinsing, system settings
- <EDIT>: Changing the current method, new method, copy and delete method
- <ESC>: <ESC> will take you back to the previous menu level
- <START/STOP>: Start and Stop of a current method

The individual functions are described in detail in [3.3 External PC Keyboard](#).

#### 3.2 Display

The display (Fig. 17) consists of a graphical LCD display with a resolution of 320 x 240 pixels. It also offers the possibility to display graphics, e.g. the measuring curve while or after the titration is/was running.

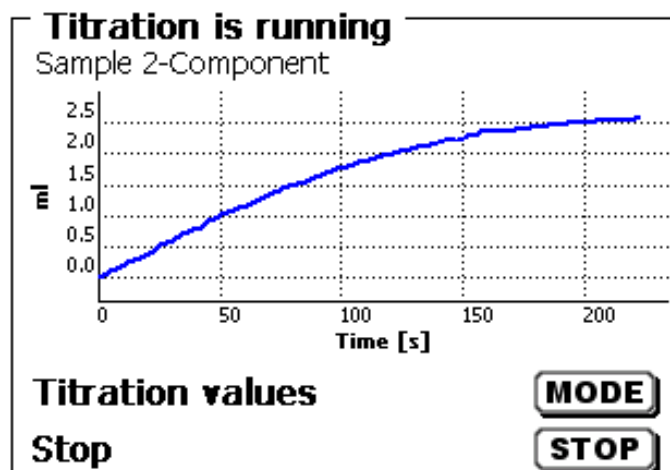
Fig. 17

### 3.3 External PC Keyboard

Tasten	Funktion
<ESC>	<ESC> will take the user to the previous level on the menu
<F1>/<START>	Start of a selected method
<F2>/<STOP>	Stop of the current method
<F3>/<EDIT>	Change of the current method, new method, copy method
<F4>/<FILL>	Without function
<F5>/ 	Display and modification of the balance data. With <Shift + F5> display and modification of the global memories
<F6>/<MODE>	Selection of method, rinsing, system settings
<F7>/<SYS>	System settings (language selection, time/date ...)
<F8>/<CAL>	Without function
<F9>/+ / -	Change of sign
<F10>/<DOS>	Start dosing menu
Num/ Scroll Lock/ Lock	Without function
Prt Sc Sys Rq	Without function
< ↑ > < ↓ > < ← > < → >	Selection of individual menus and numeric values
0...9	Input of numeric values
<ENTER>	Confirmation of input parameters
< ←Backspace >	Deletion of one input digit / an input character to the left of the flashing cursor
Letters, ASCII-symbols	Alphanumeric input possible, Uppercase and lowercase possible
All other keys	Do not have any function


### 3.4 Menu Structure

**i** The menu screens shown in this manual serve as an example and may differ from what you see!

There are 4 selection menus:


- Start or main menu
- Method parameters
- Method selection
- System settings.

After power-up, the main menu is always the first menu to appear.  
The method displayed will always be the last method that was used (Fig. 18).

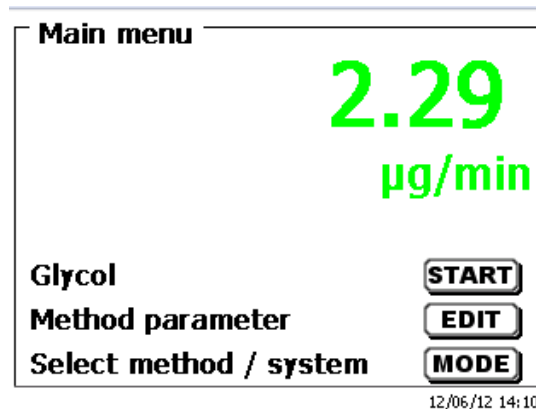**Fig. 18**

Pressing <START> will result in the immediate execution of the method shown.  
<EDIT> will take you to the method parameters (Fig. 19).

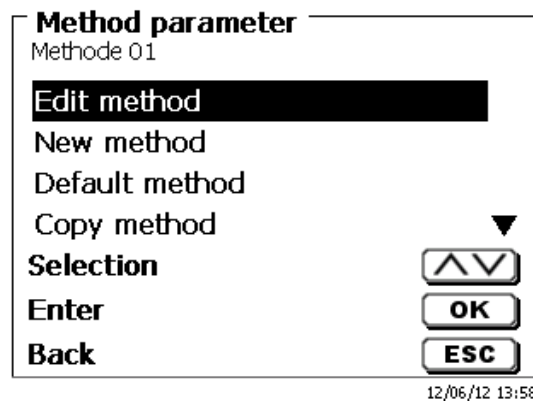
**Fig. 19**

At this point you can

- modify the current method
- create a new method
- call and memorise standard methods
- copy or delete an existing method

Use <↓> and <↑> to select the submenus.  
Confirm your selection with <ENTER>/<OK>.  
<ESC> will take you back to the main menu.


<MODE>/F6 leads you to the select method menu (Fig. 20).

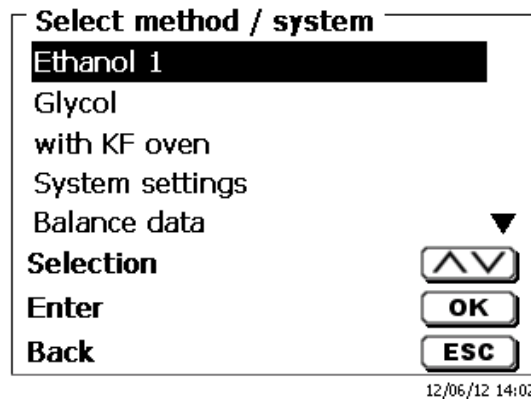
**Fig. 20**

Existing methods can be selected by pressing <↓> and <↑> and confirming the selection with <ENTER>/<OK>. Once the selection made, you will return to the main menu with the newly selected method. If no method is selected <ESC> will also take you back to the main menu.

To navigate directly to the system settings (Fig. 21 and Fig. 22) you can use the <SYS>/F7 key; you can also navigate there through the method selection menu.

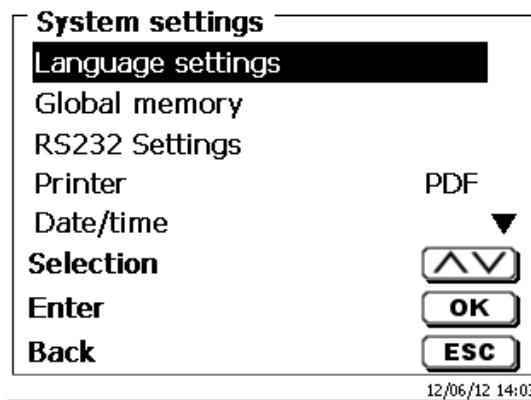**Fig. 21**

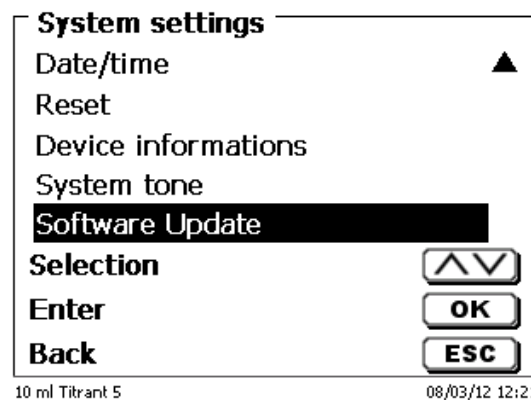**Fig. 22**

### 3.5 Main Menu

After power-up, the main menu is always the first menu to appear. The method displayed will always be the last method that was used (Fig. 23).

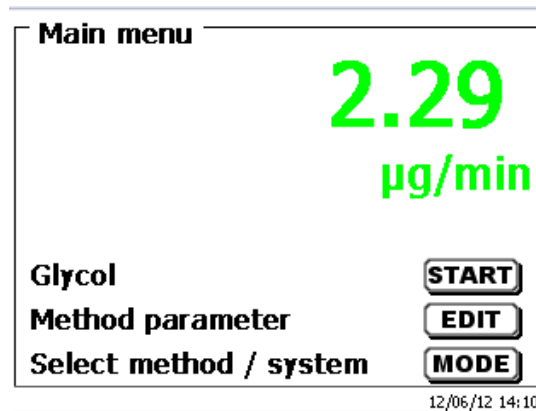
Fig. 23

#### 3.5.1 Starting a Titration

Once all preparations have been finished, you can start to titrate samples. The titrator starts automatically with the conditioning process when it switched on and analyte is present in the titration vessel. Here the titration cell is titrated until it is dry. This includes the entire water of the analyte and even adherent liquids in the titration vessel.

**i** If the start drift value is not reached (normally 10 µg/min) the method can't be started.

Start the selected method with <START> in the main menu. If the drift value is below a set value (normally 10 µg/min) you will be prompted to dose the sample (Fig. 24), for the sample identification (Fig. 25) and the weighed-in quantity (Fig. 26).

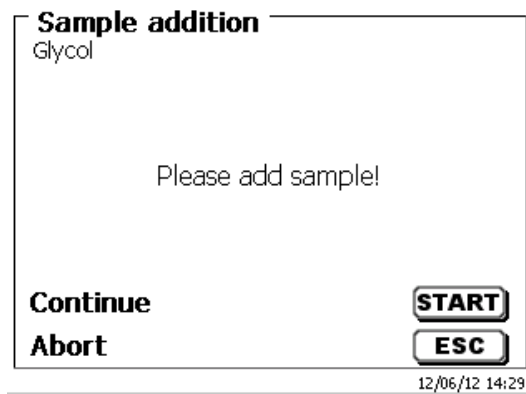
Fig. 24


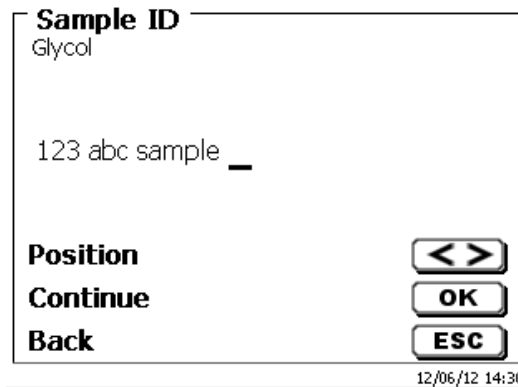
Fig. 25

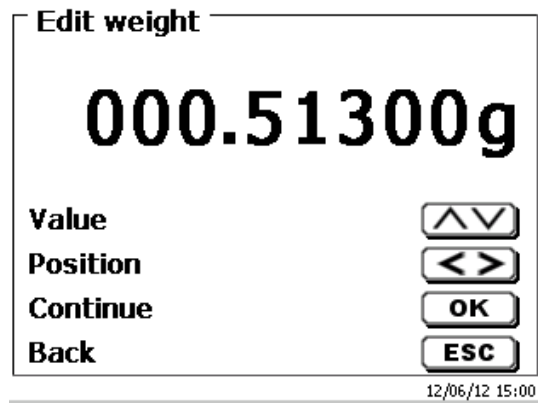
**Fig. 26**


The balance data can be entered using the front keyboard or an external keyboard. The input is to be confirmed with **<ENTER>/<OK>**.

In the case of an automatic acceptance of the balance data, the weighed-in quantities will be read in from a memory. If the memory does not contain any balance data, a message will appear to indicate that no balance data are present (Fig. 27).

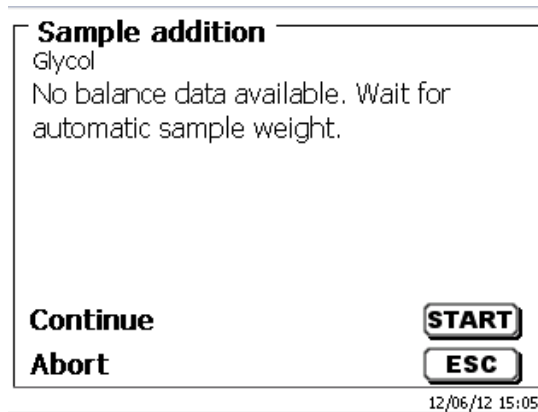
**Fig. 27**

Pressing the Print key will transfer the balance data, too. Titration will then begin directly after the transfer of the balance data without any further confirmation being necessary. The display will show the  $\mu\text{A}$ -value, the drift value and the current consumption. The top of the display will show the "Titration is running" status indication and the method being used (Fig. 28).

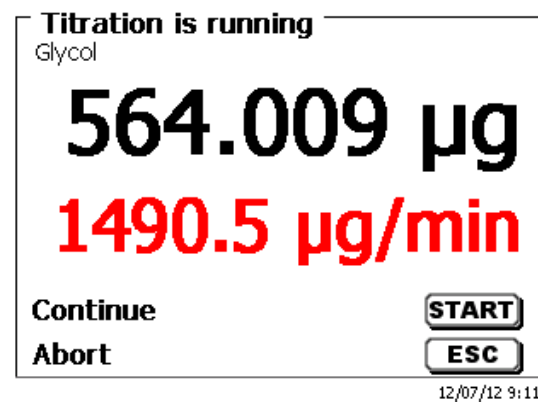
**Fig. 28**

Pressing **<F6>** will cause the titration curve to be displayed (Fig. 29).

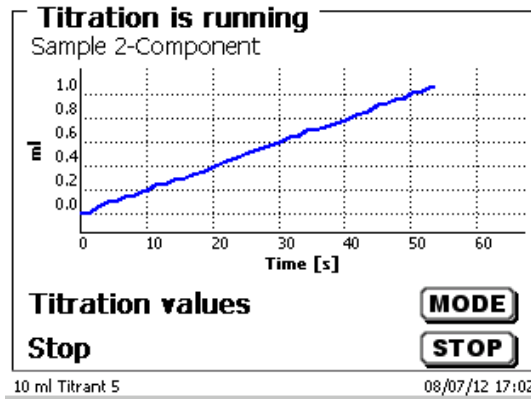
**Fig. 29**

Scaling of the chart will be done automatically. The result will be displayed at the end of the titration (Fig. 30).

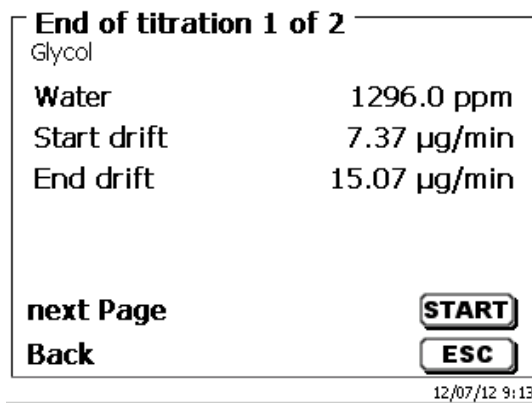**Fig. 30**

<MODE> can be used to view the titration curve or further results (Fig. 31).

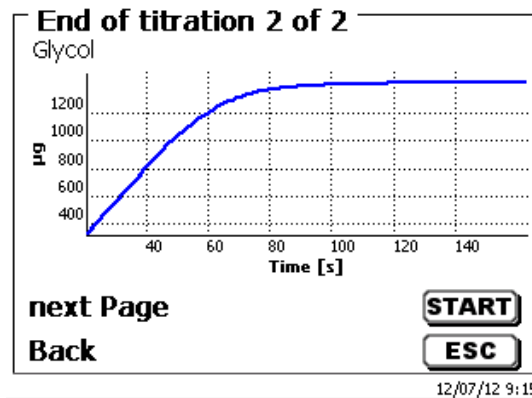**Fig. 31**

If a printer is connected, the results will either be printed according to the settings made for the method, or else they will be memorised in the form of a PDF- and CSV-file file on a connected USB stick. If no printer or USB stick is connected, you get a message on display.

<ESC> will take you back to the main menu where you can start the next titration immediately.

## 4 Method parameters

From the main menu, <EDIT> will take you to the method parameters (Fig. 32).

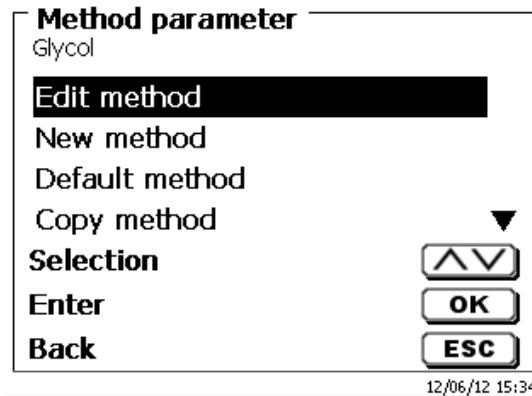
Fig. 32

### 4.1 Method editing and new method

If you select «**edit method**» or «**new method**» you will be taken to the modification or new creation of a method.

Selecting «**new method**» will always lead to the prompt for the input of a method name (Fig. 33). This prompt will not appear in the case of the modification of an already created method.

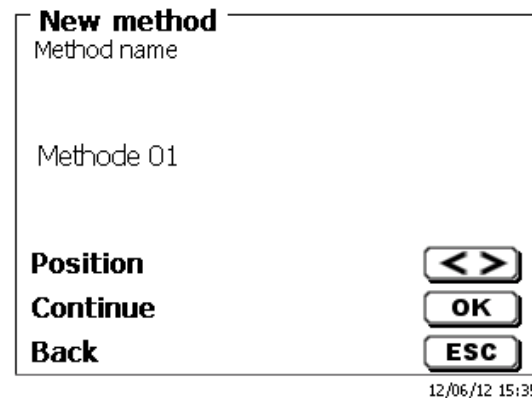
Fig. 33

The method name can contain up to 21 characters. Special characters are also possible.


**i** If no keyboard is connected, the method name being displayed **has to be** adopted.

Numbering of methods will occur automatically. Press <ENTER>/<OK> to confirm the input. The method name can be changed at any time.

Please continue at this point with  4.6 Change Method Parameters.


## 4.2 Default method

The «**Default methods**» item of the device contains a series of ready-made standard methods which can be conveniently selected (Fig. 34).

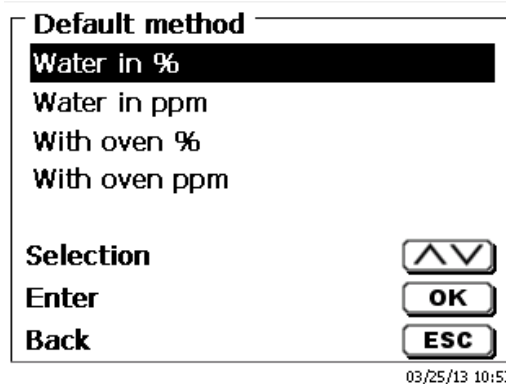**Fig. 34**

Once the selection made, you are directly prompted for the input of the method name (Fig. 35).

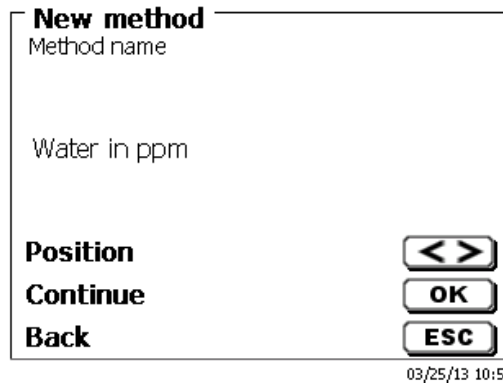
**Fig. 35**

The standard name may be adopted or modified. Subsequently, you will be taken to «**Change method parameters**».

Please continue at this point with  4.6 Change Method Parameters.

## 4.3 Copy Method

Methods can be copied or stored with a new name (Fig. 36). If you select this function, the current method will be copied and you can include a new name.

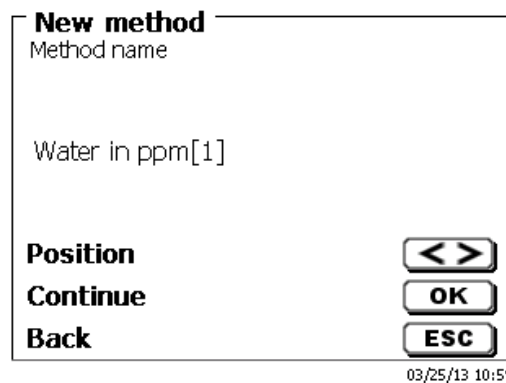**Fig. 36**

 A new name with the suffix [1] is assigned automatically in order to avoid the existence of two methods having the same name. Subsequently, you will be taken to <**Change method parameters**>.

Please continue at this point with  4.6 Change Method Parameters.

#### 4.4 Delete Method

In this function you will be prompted to know whether the current method is actually to be deleted (Fig. 37). You have to reply «Yes» in explicit terms and also confirm this reply with <ENTER>/<OK>.

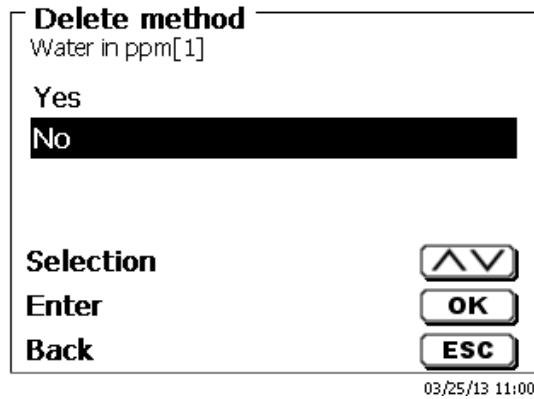
Fig. 37

#### 4.5 Print method

The currently selected method can be printed on a connected printer or stored on an USB drive as PDF file (Fig. 38).

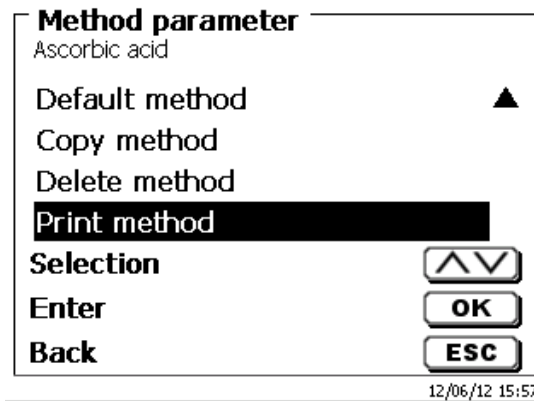
Fig. 38

## 4.6 Change Method Parameters

The input or modification of the method name was already described in  4.1 and 4.3

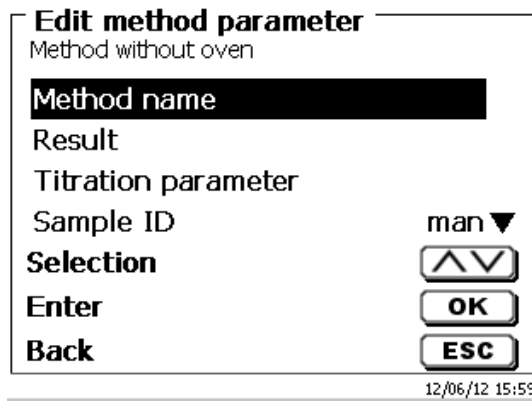
Fig. 39

### 4.6.1 Result

Result offers the following settings (Fig. 40)

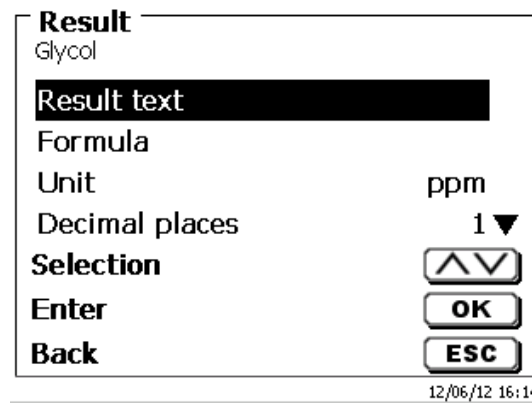
Fig. 40

The «**Result text**» may contain up to 21 alphanumeric characters including special characters (Fig. 41).


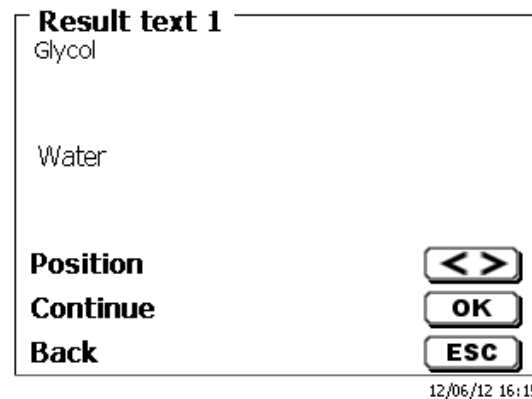
Fig. 41

Please confirm your input with <ENTER>/<OK>.

**4.6.1.1 Calculation Formula**

The appropriate calculation formula is selected on the Formula selection submenu (Fig. 42).

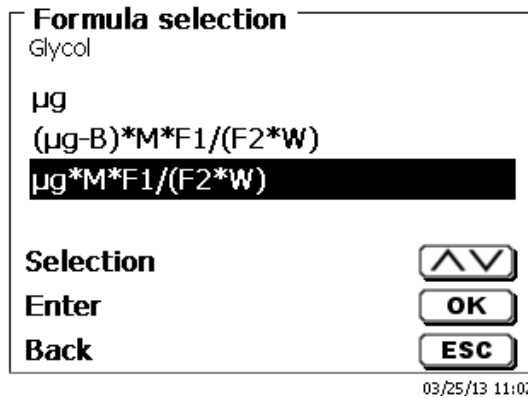**Fig. 42**

The following calculation formulae are available for automatic titration mode:


Formula	Additional information
µg	Formula for calculating only the absolute water content in µg
$(\mu\text{g}-B)*M*F1/(W*F2)$	Formula for calculating the concentration of a sample taking into account a blank value in terms of µg.
$\mu\text{g}*M*F1/(W*F2)$	Formula for calculating the concentration of a sample

The abbreviations used here have the following meaning

- µg: absolute water content
- B: Blank value in µg. Mostly determined by way of titration
- M: Mol; mol- or equivalence weight of the sample
- F1 – F2 Factor 1 - 2 conversion factors
- W "Weight", weighed-in quantity in g or volume in ml

Confirm your selection with <ENTER>/<OK>.

The values for the blank value, the titers and factors F1 – F2 can be entered or read from a global memory (Fig. 43).

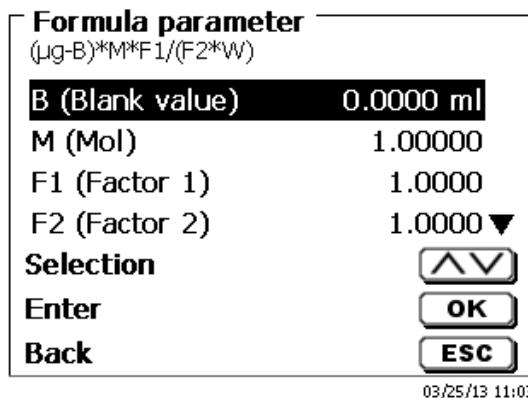**Fig. 43**

The values from the global memory were defined in advance by a titration or were manually entered (Fig. 44 and Fig. 45).

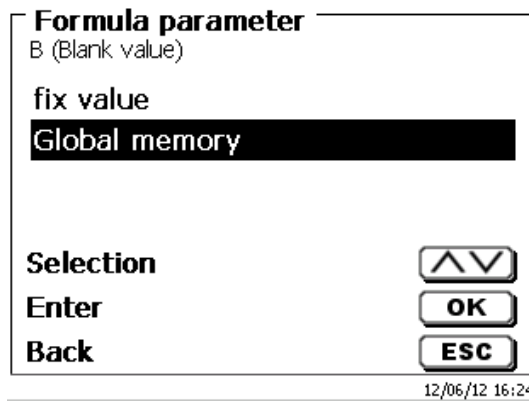
Fig. 44

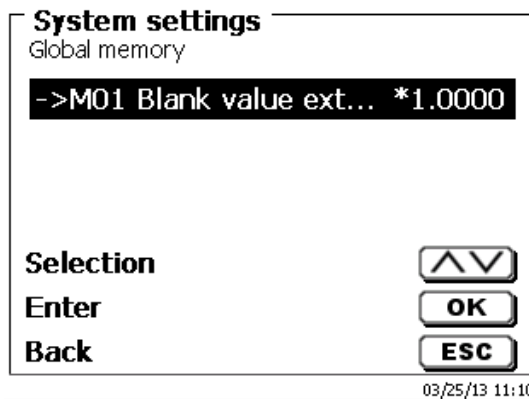
Fig. 45

The global memory used is displayed (Fig. 46).

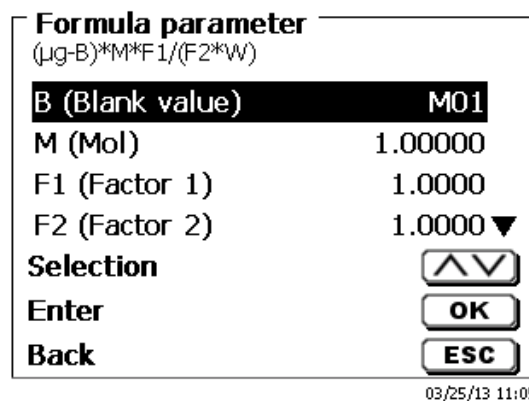
Fig. 46

Storing results in global memories is described in [4.6.1.6](#).

The values of the individual parameters of the selected calculation formula can be input one by one (Fig. 47).

**Formula parameter**  
F2 (Factor 2)

**+01000.0000**

**Value**

**Position**

**Continue**

**Back**

12/06/12 16:27

Fig. 47

#### 4.6.1.2 Sample weight and volume (sample quantity)

The Sample Quantity (W) item (Fig. 48) is used to select whether one is wishing to use a sample weight or a sample volume for titration or solution preparation (Fig. 49).

**Formula parameter**  
(µg-B)\*M\*F1/(F2\*W)

**M (Mol)** 1.00000 ▲

**F1 (Factor 1)** 1.0000

**F2 (Factor 2)** 1.0000

**W (Amount)** man

**Selection**

**Enter**

**Back**

03/25/13 11:06

Fig. 48

**Formula parameter**  
Amount

**Weight manual**

**Weight automatic**

**Fixed weight**

**Manuel Volume** ▼

**Selection**

**Enter**

**Back**

12/06/12 16:28

Fig. 49

You have the following options:

- **«Manual sample weight»:** The sample weight is enquired by a prompt at the start of the method and manually input.
- **«Automatic sample weight»:** The sample weight is automatically transferred by a connected balance.
- **«Fixed sample weight»:** A fixed sample weight is input in g. This weight will then automatically be used for each start of the method.
- **«Manual sample volume»:** The sample volume in ml is prompted at the start of the method and manually input.
- **«Fixed sample volume»:** A fixed sample volume is input in ml. This volume will then automatically be used for each test of the method.

### 4.6.1.3 Formula unit

The formula unit can be selected in the «Unit» submenu (Fig. 50).

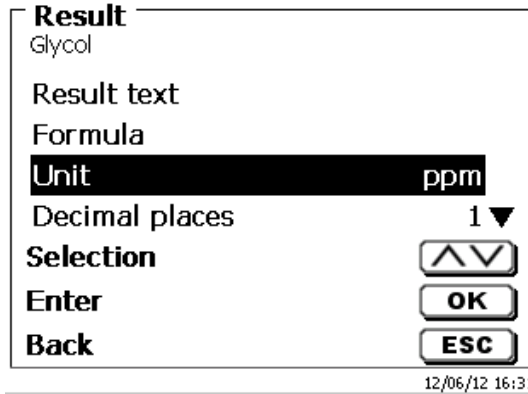
Fig. 50

Once the selection made (e.g. «%»), the unit will also be displayed as piece of information (Fig. 51).

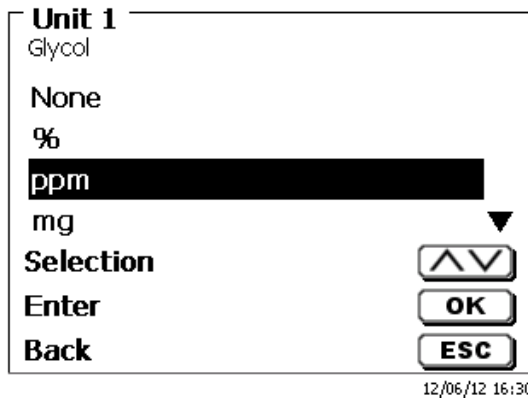
Fig. 51

By pressing the «INS» (Insert) key on the external keyboard, you can also add new units.

### 4.6.1.4 Decimal digits

To conclude, it is possible to determine the number of decimal digits from 0 - 6. The standard setting is 1 (Fig. 52).

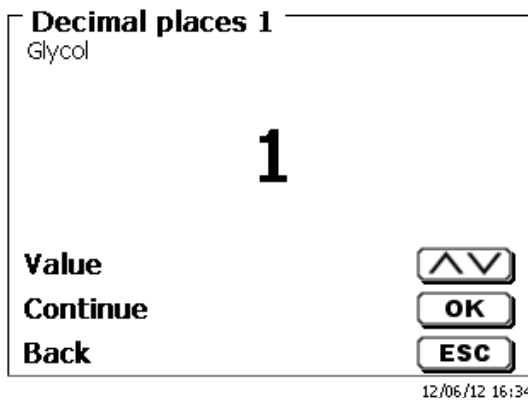
Fig. 52

### 4.6.1.5 Statistics

The mean value and relative standard deviation can be automatically calculated and documented by using the statistics (Fig. 53).

**Result**  
Glycol

Formula ▲

Unit ppm

Decimal places 1

**Statistics** None ▼

Selection ▲▼

Enter OK

Back ESC

12/06/12 16:33

**Fig. 53**

The calculation of the mean value is already possible from 2 individual values, the calculation of the relative standard deviation is only possible from 3 single values (Fig. 54). The maximum quantity is 10.

**Statistics Formula 1**  
Glycol

None

2

**3**

4 ▼

Selection ▲▼

Enter OK

Back ESC

12/06/12 16:33

**Fig. 54**

#### 4.6.1.6 Global Memories

Results of titrations can be written into one of the 50 global memories (M01 - M50) for additional calculations (Fig. 55).

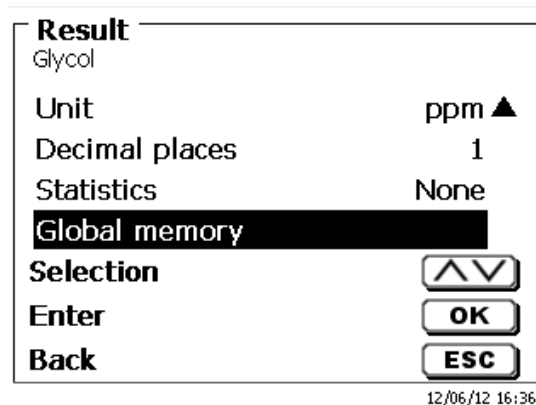
Fig. 55

The mean value is written into the global memory when the statistic is switched on. You enter the submenu with <ENTER>/<OK>. If a global memory has not been created, a memory can be created by using the insert key <INS>. The titrator proposes a memory name, such as **M01** (M01 - M50). The name of the memory can be changed in reference to the application (Fig. 56). This simplifies later the allocation of the global memory in another method.

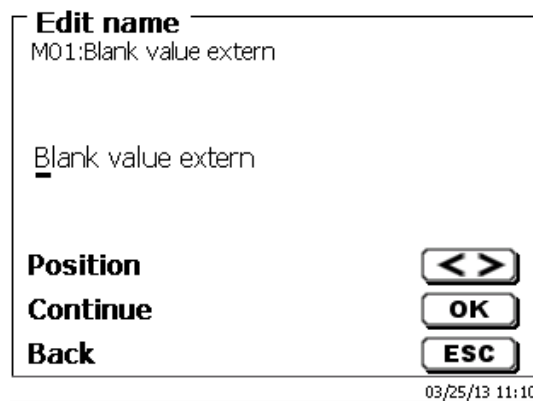
Fig. 56

**Example:** The blank value of a chloride titration is defined with the support of an extra method. The result in ml is thereby automatically written into global memory M01 by using the name "Blanc value" (Fig. 57). The blank value is then automatically deducted from the titrant consumption within the chloride method.

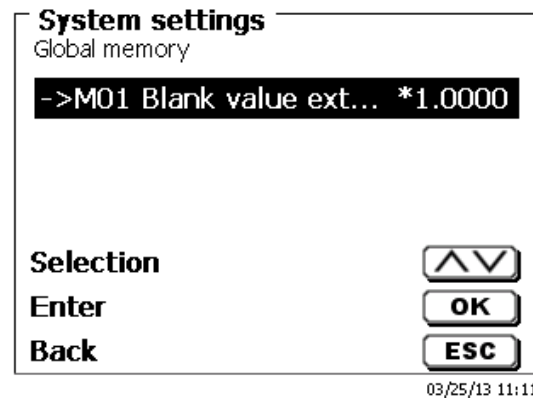
Fig. 57

The menu for the global memory can always be accessed by pressing <Shift> or via system settings. The name or values can be changed by using <EDIT> and have the methods shown that are used in the global memories (Fig. 58).

**Global memory**  
M01: blank external extr.

**Edit name**

**Edit value**

writing method

reading method ▼

**Selection** ▲▼

**Enter** OK

**Back** ESC

12/06/12 16:41

Fig. 58

## 4.6.2 Formula Editor

The formula editor is intended to complement the existing standard formulas. The standard formulas are loaded and can then be modified. The original standard formula itself is never changed.

### 4.6.2.1 Start and Work with the Formula Editor

With <EDIT> you are going to «Edit method», «New method» or «Default methods» and select then «Result» (Fig. 59).

**Edit method parameter**  
Methode 01

**Method name**

**Method type** auto

**Mode** Dynamic

**Result** ▼

**Selection** ▲▼

**Enter** OK

**Back** ESC

20 ml NaOH 02/13/17 7:58

Fig. 59

Confirm the selection «Result» with <ENTER>/<OK> (Fig. 60):

**Result**  
Methode 01

**Calculation options** 1 EQ

**Formula**

**Selection** ▲▼

**Enter** OK

**Back** ESC

20 ml NaOH 02/13/17 7:59

Fig. 60

Select "Formula" with <ENTER>/<OK> (Fig. 61).

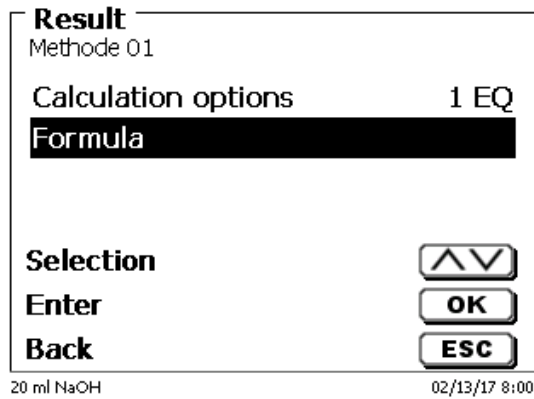
Fig. 61

The following selection appears (Fig. 62).

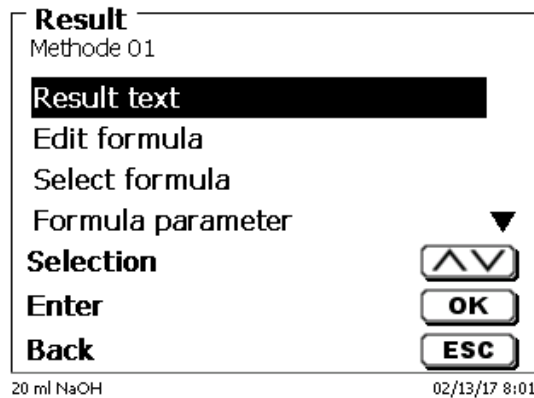
Fig. 62

You can select the existing methods with <↓> and <↑> and confirm the selection with <ENTER>/<OK> (Fig. 63).


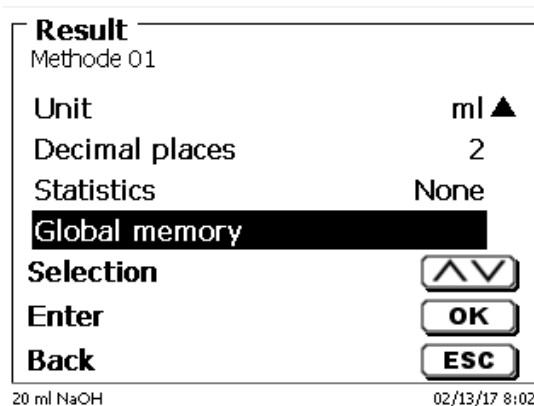
Fig. 63

«Result text», «Select fomula», «Formula parameter», «Unit», «Decimal places», «Statistics» and «Global memory» don't differ from previous versions.

**i** New is the menu item «Edit formula»!


If you select «**Edit formula**» and confirm it with <ENTER>/<OK> the currently selected formula is displayed (Fig. 64).

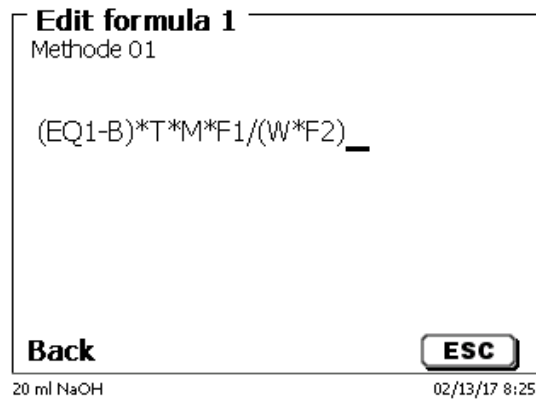
**Fig. 64**

The formula **(EQ1-B)\*T\*M\*F1/W\*F2)** can now be changed and confirmed with <ENTER>/<OK> after the change. If you leave the editor with <ESC> the formula remains unchanged.


You can use the backspace key ← to delete the formula characters from the end (Fig. 65) or use the cursor left and right keys to select the locations and then use the <DELETE> key to delete the selected formula character or a value (Fig. 66 and Fig. 67).

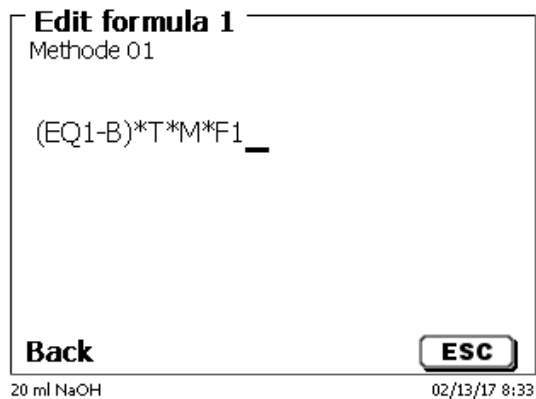**Fig. 65**

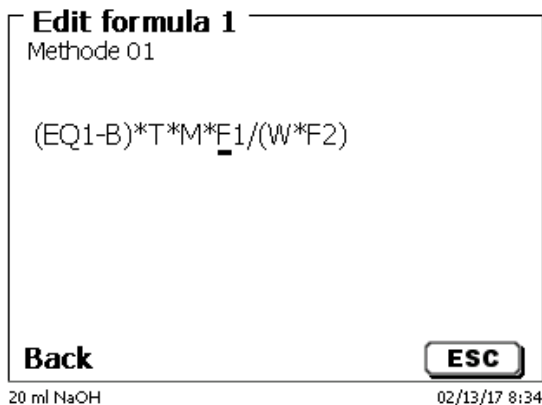
**Fig. 66**

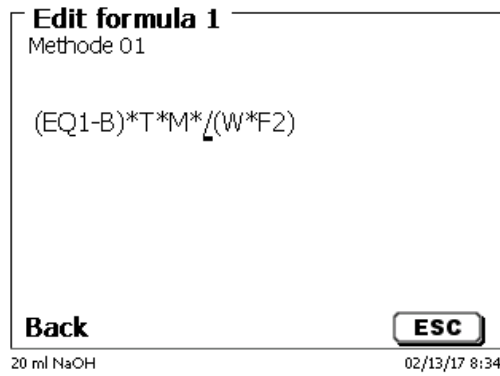
**Fig. 67**

Instead of the formula character **F1** you can now use e.g. directly enter a numeric value (Fig. 68).

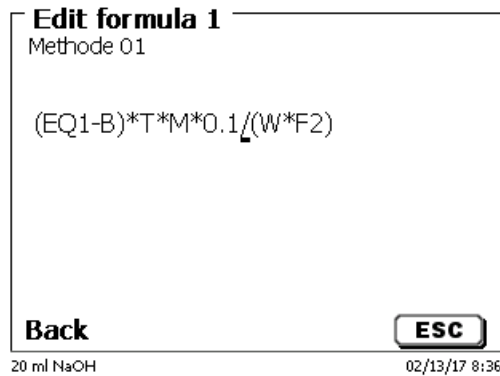**Fig. 68**

The decimal point of the numeric value can be entered as a point or a comma. Press <ENTER>/<OK> to leave the editor. The formula is automatically saved. Under «**Formula parameter**», the values can be entered as before (Fig. 69).

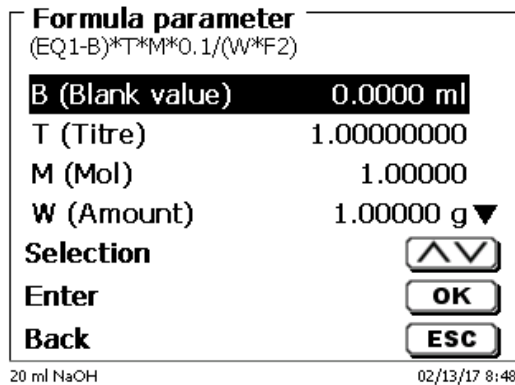**Fig. 69**

#### 4.6.2.2 Applicable Formula Characters, Arithmetic Operations and Values

The following arithmetic operations can be used:

Arithmetic operations	Formula character
• Addition	+
• Subtraction	-
• Multiplication	*
• Division	/
• Calculations with brackets to 25 levels	()
• Logarithm to base 10	L
• Exponential function	^

The following formula characters are available:

Formula characters	Meaning
EP1, EP2, EQ1, EQ2	Results from a titration like e.g. EQ1, EQ2 etc
F1 –F10	Values which can contain fixed, manual, global memory or results of other formulas.
T	Titer of the titration burettes
W	Weight sample
B	Blanc value
D	Density
S	Slope in ml/s of a pH stat-application
EV	End- or total volume of a titration. Is needed if you want to calculate the difference between an equivalence point EQ or end point EP to the total (end) volume
M	Molecular mass or equivalent weight
M01-Mxx	Global memories
R1-2	Result of a formula calculated previously in the application.

**i** If a global memory Mxx is used, which is not created, this is created automatically and assigned the default value 1.

**i** Only results of the preceding formulas can be used. This is checked in the syntax check.

#### 4.6.2.3 Syntax check

The syntax check is performed each time the formula is saved by the formula editor.

It is checked,

- whether the number of opening brackets is equal to that of the closing ones.
- whether the entered variables and calculation operations are allowed.

If an error occurs in the syntax, an error message is displayed (Fig. 70 and Fig. 71).

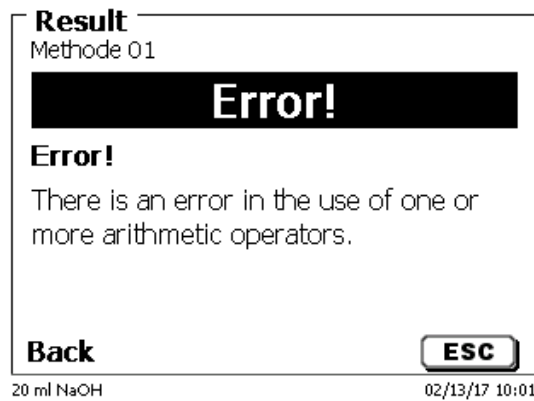
Fig. 70

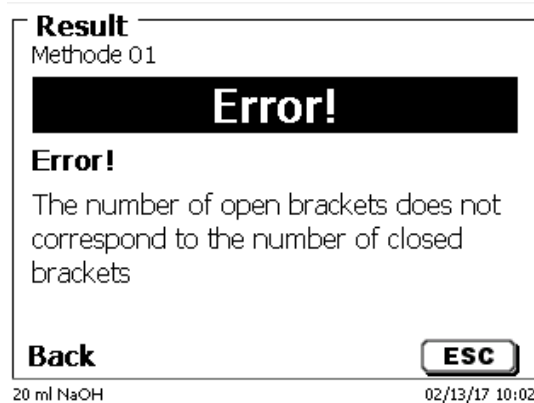
Fig. 71

### 4.6.3 Titration parameters

The «**Titration parameter**» submenu is used to determine the actual parameters of the method (Fig. 72).

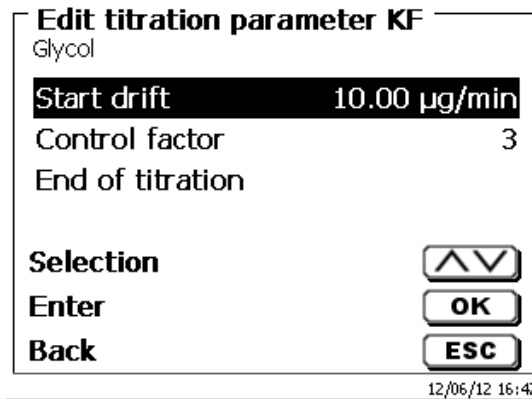
Fig. 72

#### Generally applicable titration parameters

The following parameters are valid:

- Start drift
- Control factor
- End criteria:
  - Maximum titration time
  - Minimum titration time
  - Nachlaufzeit Stop delay time
  - Working point
  - Stop drift (delta)
  - Stop drift tolerance

#### 4.6.3.1 Start Drift

The value of the «**start drift**» in µg/min must be equal or even lower in order to start a titration. If the value is exceeded, the TitroLine® 7500 KF *trace* is in the conditioning mode. The standard value is 10.00 µg/min. This value can be entered from 0.01 to 99 µg/min (Fig. 73).

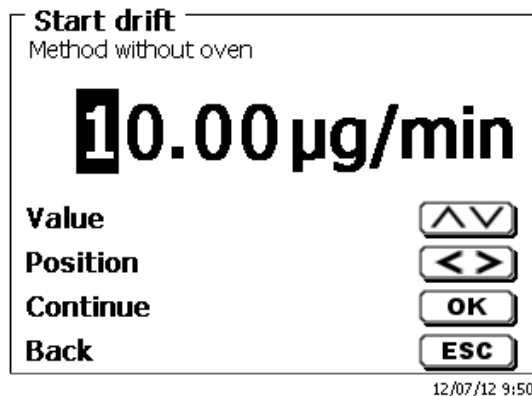
Fig. 73

#### 4.6.3.5 Stop delay time

Is the time in seconds in which the drift-stop criteria are being checked. The «stop delay time» can be set between 0 and 60 seconds. The pre-set standard value is 5 s (Fig. 77).

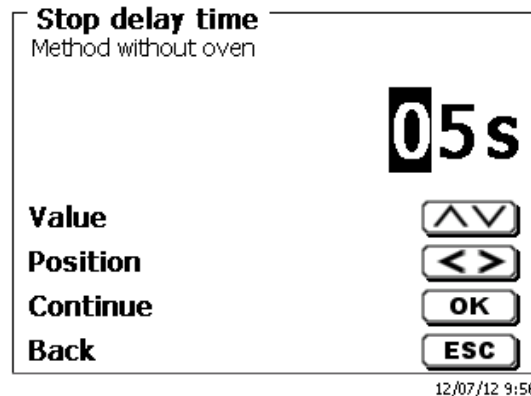
Fig. 77

#### 4.6.3.6 Working point

The «working point» in mV is the base value for the indicator electrode. It can be set from 1 to 1000 mV. 300 mV is pre-set as standard value and works for many reagent/reagent combinations. A higher value is maybe sometimes necessary (Fig. 78).

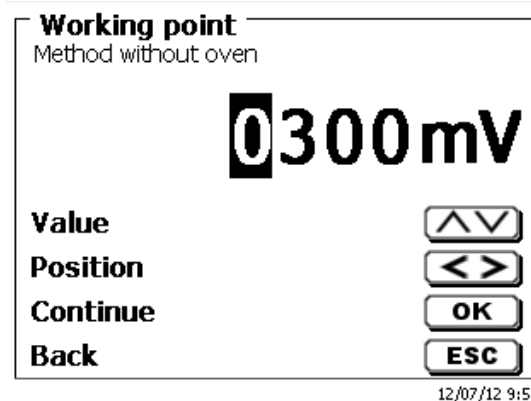
Fig. 78

#### 4.6.3.7 Stop Drift (delta)

The «stop drift» in  $\mu\text{g}/\text{min}$  is not an absolute end value. The end value of a titration is always the currently measured drift + stop drift (Fig. 79).


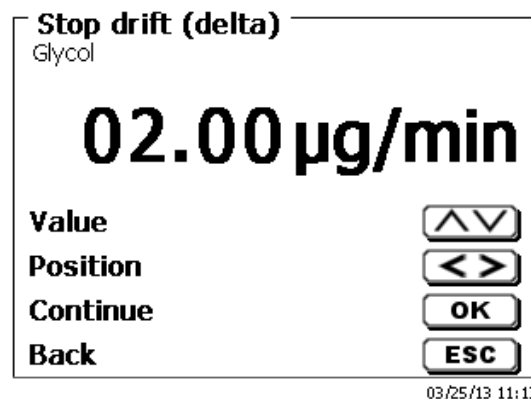
Fig. 79

**An example:**

If the current drift shows a value of 1.5 µg/min and the stop drift is set to 2.0 µg/min, then the end drift actually to be reached is 3,5 µg/min.


This means: The lower the entered stop drift, the longer takes the measurement. If a high stop drift (e.g. 20 µg/min) is entered, the measurement ends significantly faster. Low value = exact measurement, high value = inexact measurement. As stop drift 2 µg/min are pre-set as standard value. The value can be entered from 0.01 to 25 µg/min.

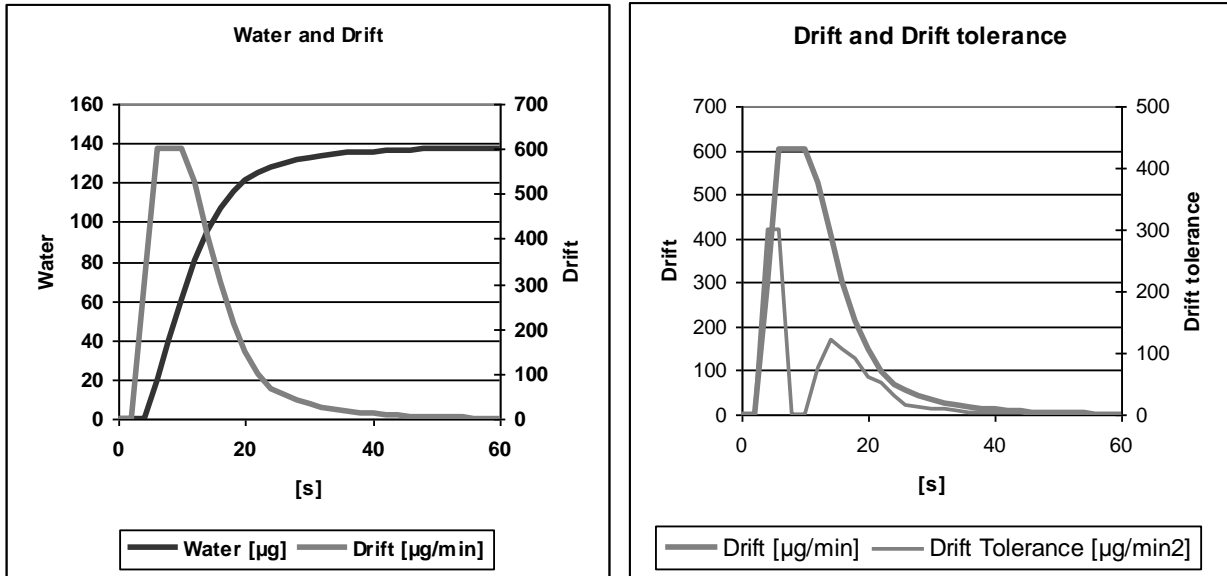**Diagram 1** shows on the left the titration curve with the measured variable water per time and the derived variable drift per time.


**4.6.3.8 Stop Drift Tolerance**

As stop drift tolerance 0.02 µg/min is pre-set as standard value. The value can be entered from 0.01 to 25.00 µg/min (Fig. 80).

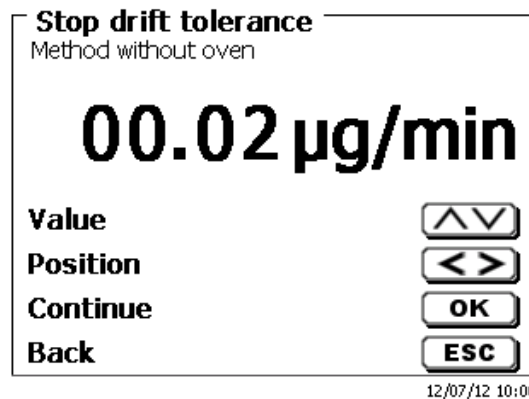**Fig. 80**


In case the drift stop has not been reached as criterion, as new stop criterion is the stop drift tolerance will be used. It is the derivative of the drift with respect to time. Diagram 2 shows the typical progression of the stop drift tolerance. Thus, the automatic stop of titrations with side reactions is possible

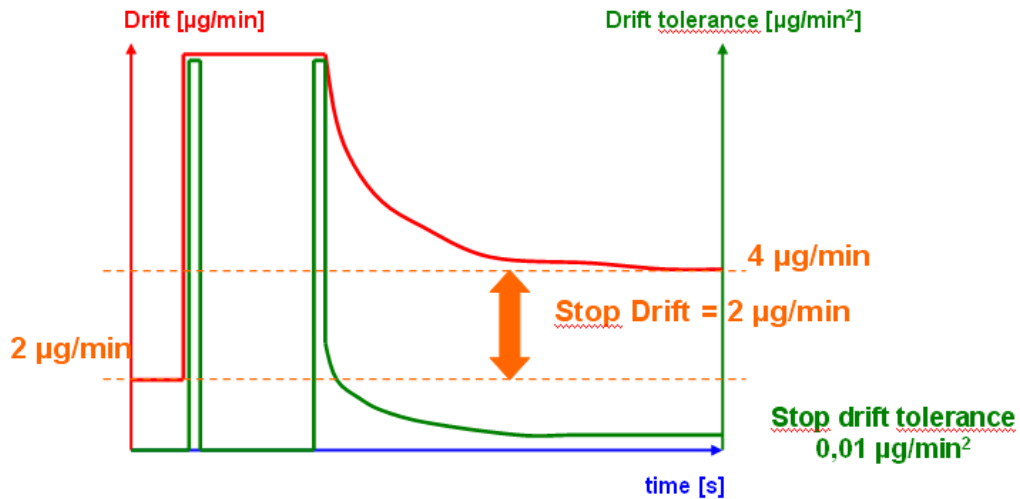**Diagram 2:** The criterion stop drift as difference to the start drift, the stop drift tolerance as stability criterion for the drift. Only one of the two end criteria stop drift and stop drift tolerance must be met to end the measurement.


#### 4.6.4 Sample identification

Here you determine whether a «sample identification» is used. You can set a «manual», «automatic» and «no sample name» (Fig. 81).

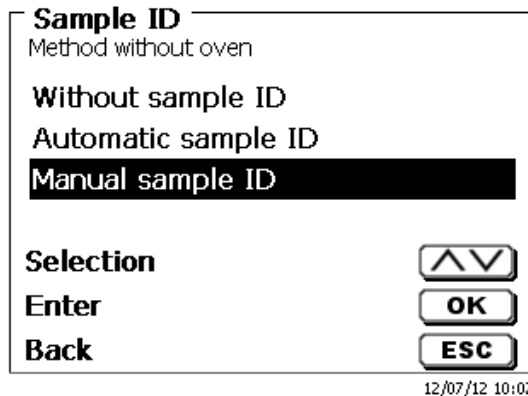**Fig. 81**

With manual sample identification, the sample identification is always asked for after the start of the method (see also 3.5 Main Menu). With the automatic sample identification a root identification is defined (here sample, see Fig. 82), which is then automatically numbered starting with 01.

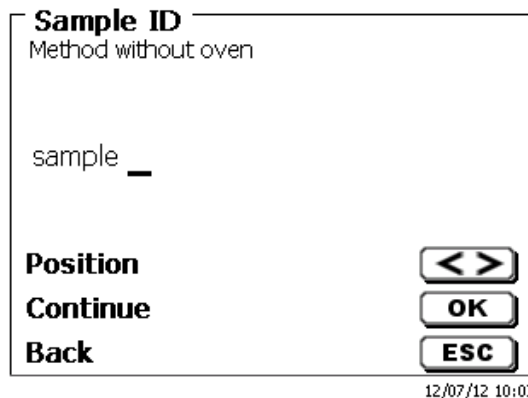**Fig. 82**

After a new power-up, numbering will resume with 01.

### 4.6.5 Documentation

Three different format settings are available for documentation (Fig. 83) on a printer or USB device: «**short**», «**standard (with curve)**» and «**GLP**» (Fig. 84).

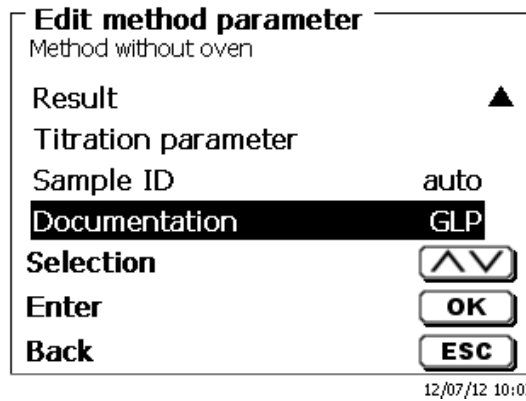
Fig. 83

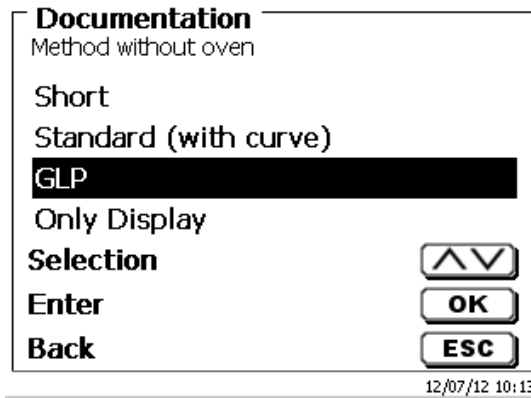
Fig. 84

Short documentation	Standard documentation	GLP-documentation
Method name, date, time, duration of titration, sample description, weight/volume, starting and end measurement values, results and calculation formula	Same as 'Short documentation' + titration curve	Same as 'Standard documentation' + method contents

## 5 System settings

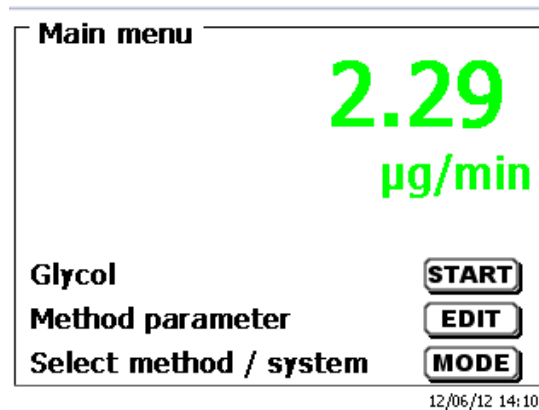
Fig. 85

From the main menu (Fig. 85) you can access the system settings with <SYS> (Fig. 86).

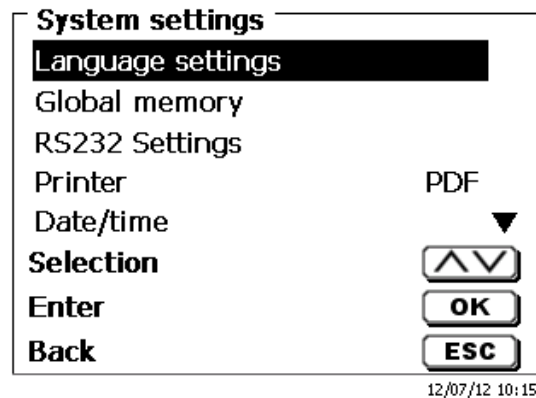
Fig. 86

Setting the national language was already described in [2.4](#).

### 5.1 Global Memory

Working with global memories was already described in [4.6.1.6 Global Memories](#).

### 5.2 RS-232 Settings

The «RS232 settings» item can be used to determine the device address of the TitroLine® 7500 KF *trace* and set the parameters of the two RS-232 interfaces independent from each other (Fig. 87).

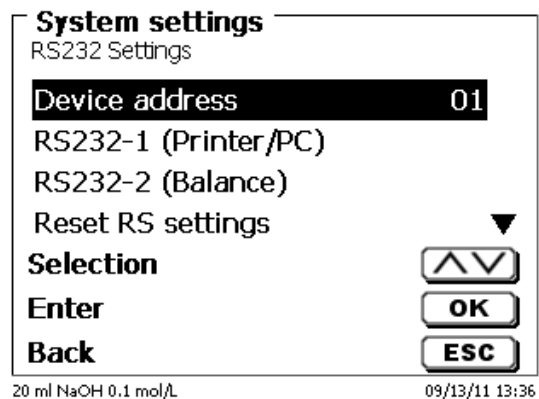
Fig. 87

The device address can be set from 0 - 15. Address 1 is the default setting (Fig. 88).

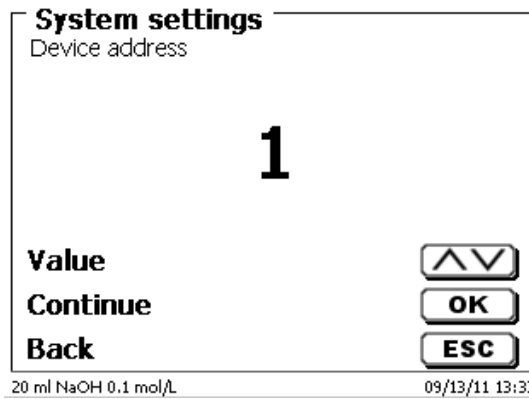
Fig. 88

The baud rate is preset to 4800 (Fig. 89).

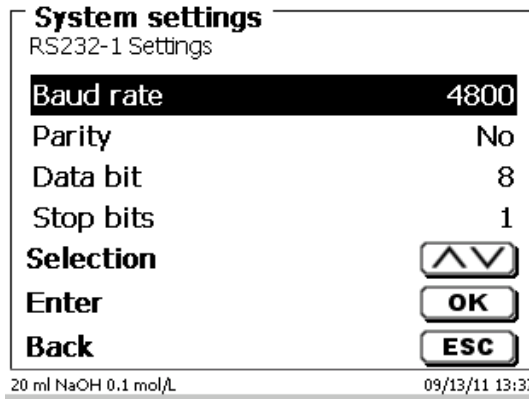
Fig. 89

It may be set to 1200 - 19200 (Fig. 90).



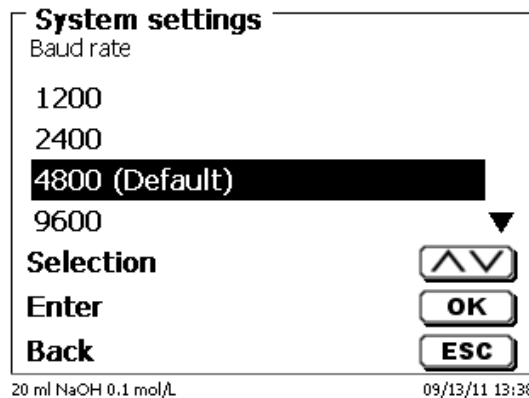
Fig. 90

The parity can be selected amongst «No», «Even» and «Odd». «No» is the default setting (Fig. 91).

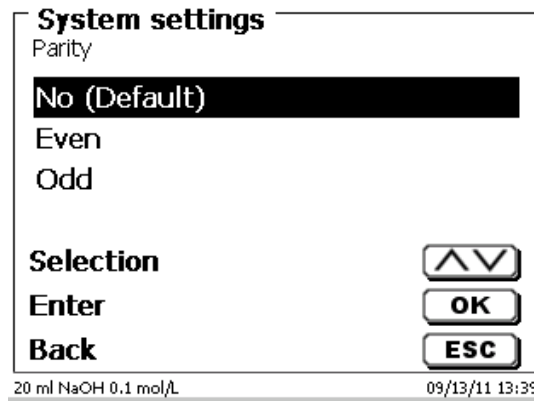**Fig. 91**


You may select between 7 and 8 data bits. 8 bits is the default setting (Fig. 92).

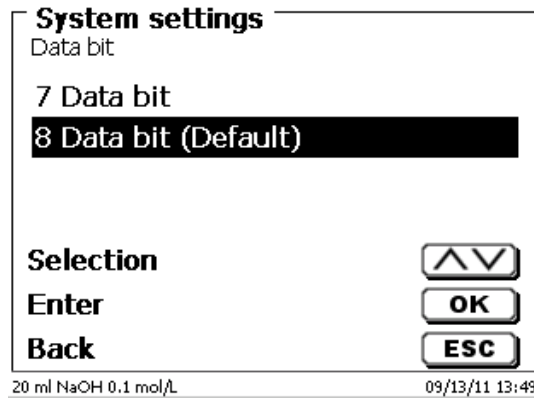**Fig. 92**

**i** The RS232 parameters can be set to the factory settings.

The RS232-1 can be converted from RS on USB (Fig. 93 and Fig. 94).  
In this case, the titrator via the USB PC connection to the PC is connected.

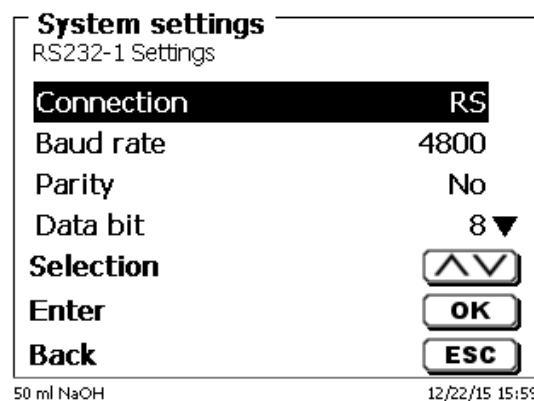**Fig. 93**

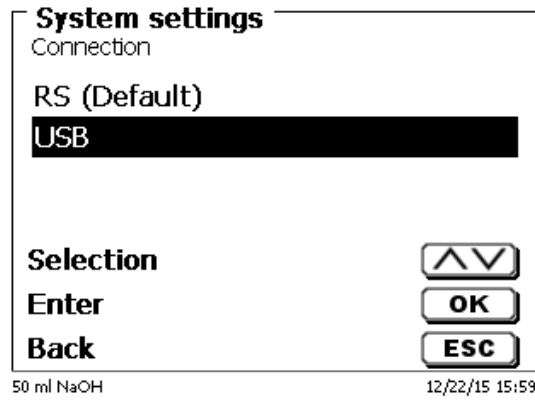
Fig. 94

For the USB connection, a driver must be installed on the PC side.

**i** The driver can be downloaded from the manufacturer website.

### 5.3 Date and Time

The factory time setting is Central European Time. This setting may be changed, where necessary (Fig. 95).

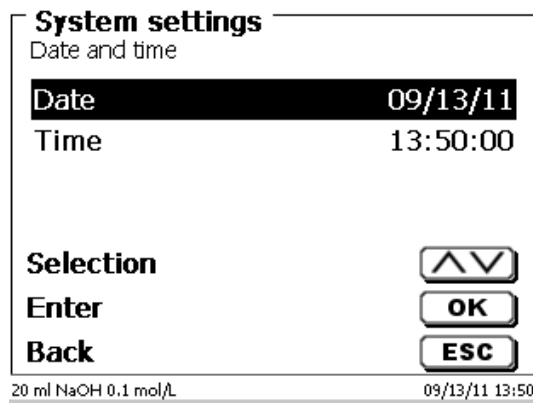
Fig. 95

## 5.4 Password

**i** Please read the instructions before you activate the password!

**i** When you activate the user management the first time, a user with administrator rights are created automatically. **Important for this first Administrator: Please note your password and user name. If you forget it, you do not have access to the device anymore!** In this case, please contact the service (see backside of this document).

The administrator can create new users with different access levels to the instrument software.

**i** The TITRONIC® 500 and TitroLine® 6000 allow maximum 5 users and all 7XXX titrators up to 10 users.

### 5.4.1 Creation of the first Administrator

Go to «System settings» and select «User management» (Fig. 96).  
Confirm the selection with <ENTER>/<OK>.

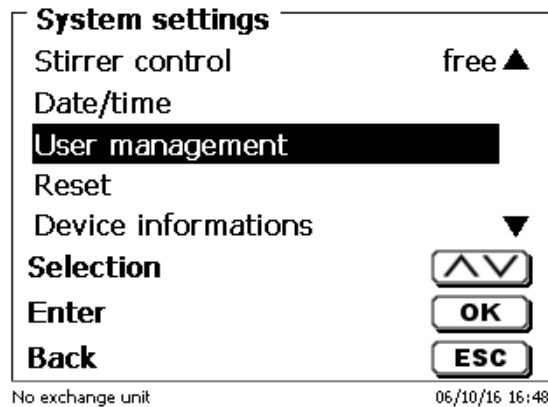
Fig. 96

«Activate» the User management with <ENTER>/<OK> (Fig. 97).

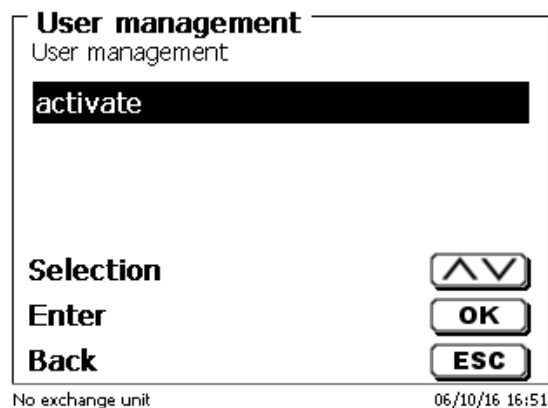
Fig. 97

Enter a user name (Fig. 98).

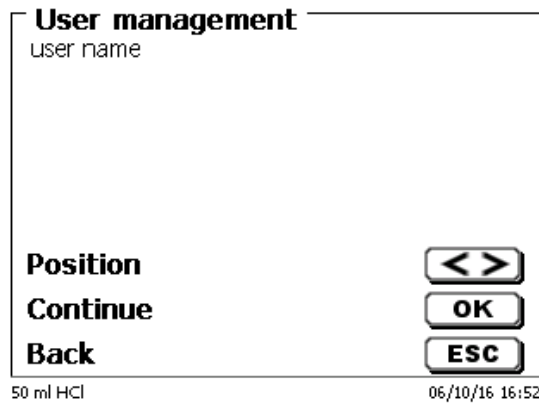
Fig. 98

It could be your first name, also the function like «admin» or more simple like «ad» (Fig. 99).

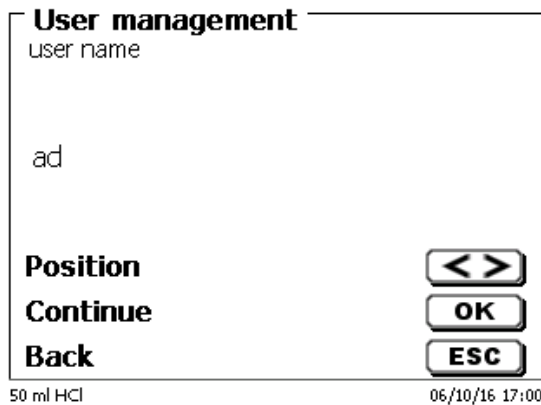
Fig. 99

Confirm with <ENTER>/<OK>.

You have to enter now your full user name (full name) and then your password (Fig. 100).

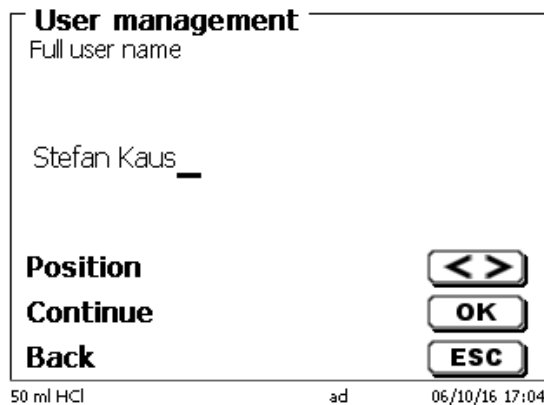
Fig. 100

The password must have at least **5 characters**.  
Allowed are all alphanumeric signs in **lower** and also **capital** letters.  
A simple example is:

**Abc12**


**i** When you activate the user management the first time, a user with administrator rights are created automatically. **Important for this first Administrator: Please note your password and user name. If you forget it, you do not have access to the device anymore!** In this case, please contact the service (see backside of this document). We need only the serial number of the device. Then we can create a master password for the device which is valid for one week

If you do not enter the password an error message appears (Fig. 101).


**Fig. 101**

Go back with <ESC> and enter then a password (Fig. 102).

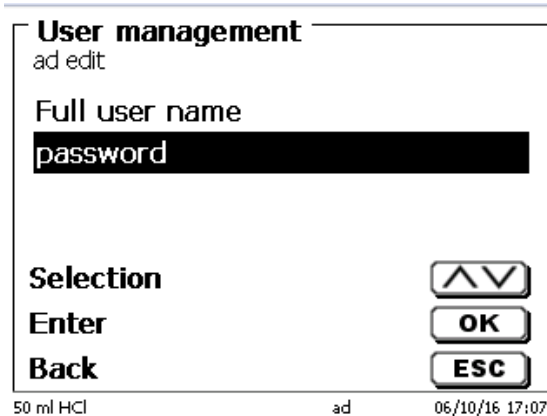**Fig. 102**

Confirm the selection with <ENTER>/<OK> (Fig. 103).

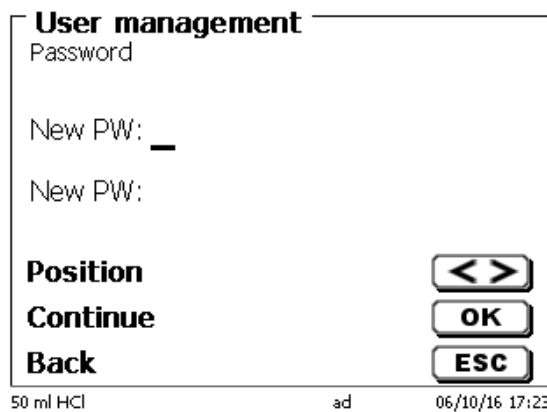**Fig. 103**

Enter the password two times and confirm with <ENTER>/<OK> (Fig. 104).

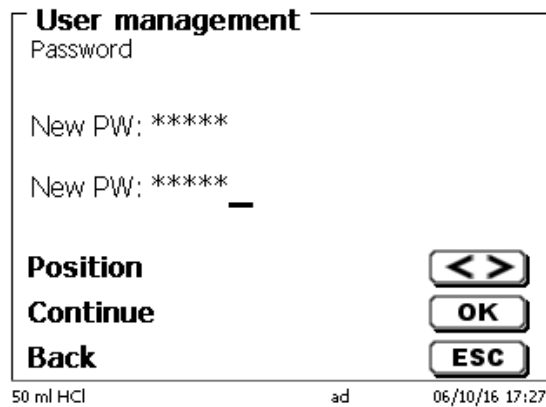
Fig. 104

Go back to the main menu with <ESC>.

You are logged in as administrator and have full access to all levels and menus.

You can see the user name at the bottom line of the display. Here in the example (Fig.10) it is «ad» (Fig. 105).

Fig. 105

As administrator you have the rights to create new users with different levels. If you start the titrator now you have to activate the user with **ctrl+L**.

**i** Without one active user it is not possible to work properly with the device!

Possible are only

- the change of the exchange heads
- the FILL function works
- and the F10 DOS function works

When you have entered the user name and password you have full access to all menus.


### 5.4.2 Creation of additional users

The administrator has the rights to create additional new users (Fig. 106).


**Fig. 106**

Confirm with **<ENTER>/<OK>**. Enter the user name of the new user. The minimum characters are here two. In the example (Fig. 107) it is "Michael".

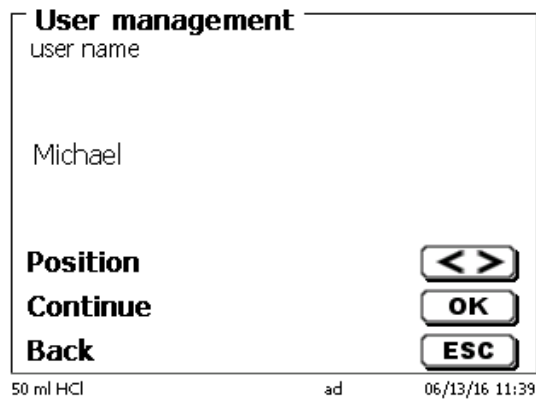**Fig. 107**

You have to enter the full user name. Possible are between 2 and 20 characters (Fig. 108 and Fig. 109). Confirm with **<ENTER>/<OK>**.

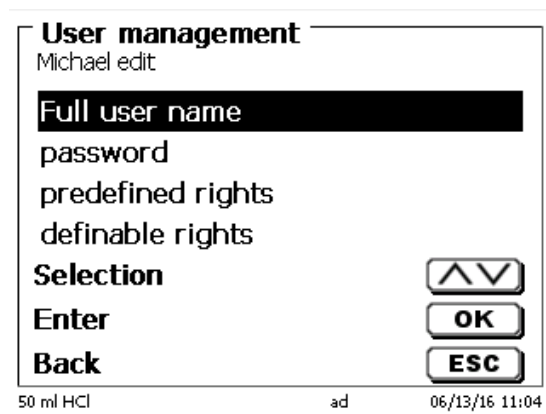**Fig. 108**

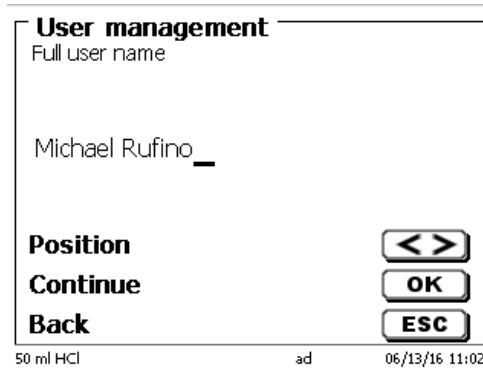
Fig. 109

You have to enter the password (Fig. 110 and Fig. 111).  
Confirm with <ENTER>/<OK>.

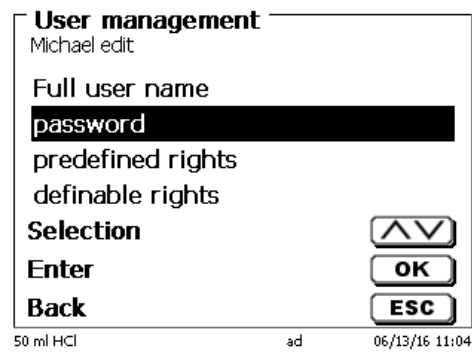
Fig. 110

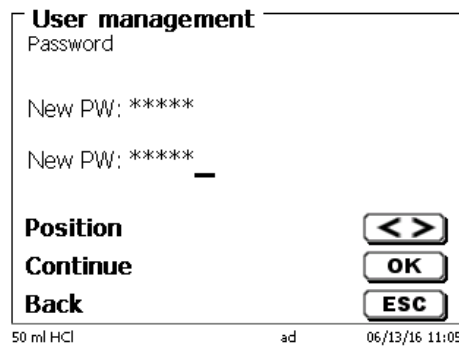
Fig. 111

### 5.4.3 Predefined rights and definable rights

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**Fig. 113**

The «**extended user**» has similar rights as the «**administrator**» but do not have access to the user management and not able to delete existing methods but can be edit methods.

The <**user**> has limited rights and no access to systems settings.  
The edit of existing methods is not possible with the user rights.

It is possible to change the access rights for all three levels of user (see [5.4.3.2 Definable rights](#)).

**i** Not possible is to change the rights from the first Administrator!

The table below shows the access rights for the three predefined users:

Menu access/functions	User	Extended user	Administrator
System settings	No	Yes	Yes
User management	No	No	Yes
RS settings	No	Yes	Yes
In / export	No	Yes	Yes
Exchange unit	No	Yes	Yes
Electrode menu	No	Yes	Yes
Global memory	No	Yes	Yes
Method selection	Yes	Yes	Yes
Edit, new, default, copy methods	No	Yes	Yes
Print methods	Yes	Yes	Yes
Delete methods	No	No	Yes
Start method	Yes	Yes	Yes
Start CAL	Yes	Yes	Yes
FILL	Yes	Yes	Yes
Update	No	Yes	Yes
Dose with F10	Yes	Yes	Yes
Output/print	Yes	Yes	Yes
Rinsing	Yes	Yes	Yes
New calculation	Yes	Yes	Yes
Edit balance data	Yes	Yes	Yes
Printer	No	Yes	Yes
Communication via RS	Yes	Yes	Yes
Network setting	No	No	Yes

Yes = access  
No = no access

### 5.4.3.2 Definable rights

If you have created a new user, you can define all rights in the menu «definable rights» (Fig. 114).

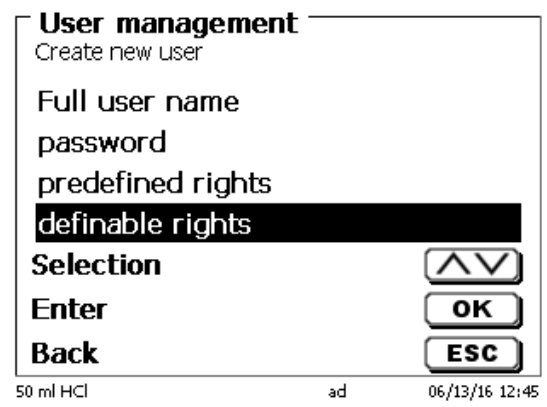
Fig. 114

Confirm the selection with <ENTER>/<OK>.

The default settings are always from a **user** if you do not have selected the extended user before.

**X** means **no access**, **W** means **access**. You can change with <ENTER>/<OK> from **X** to **W**. Below you can see all possible definable rights (Fig. 115 - Fig. 120).

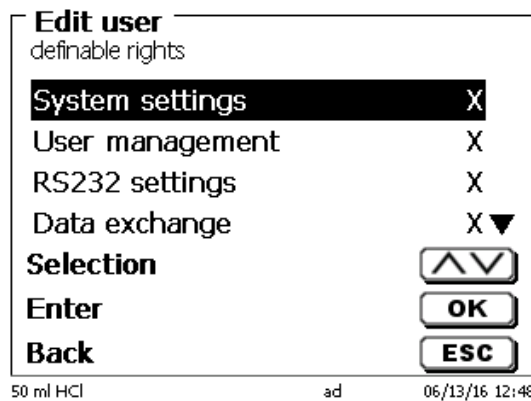
Fig. 115

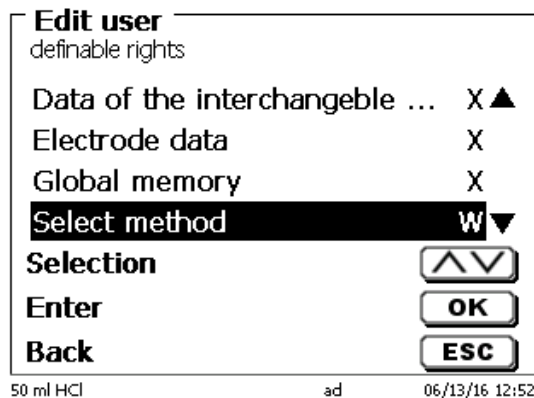
Fig. 116

### 5.4.4 Delete of users

It is possible to delete a single user with the <DEL> key on the external keyboard. You select the user with the up and down keys and then press on <DEL> (Fig. 121).

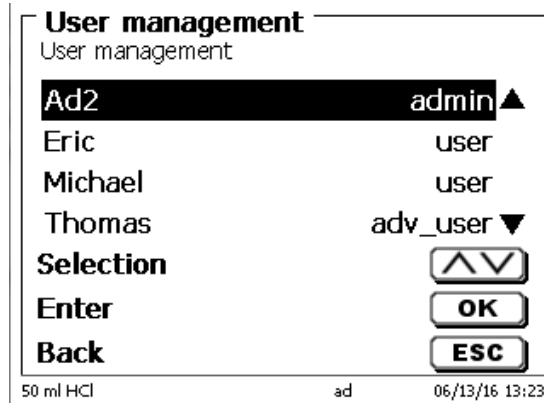
Fig. 121

After <DEL> the user is immediately deleted without any additional request (Fig. 122).

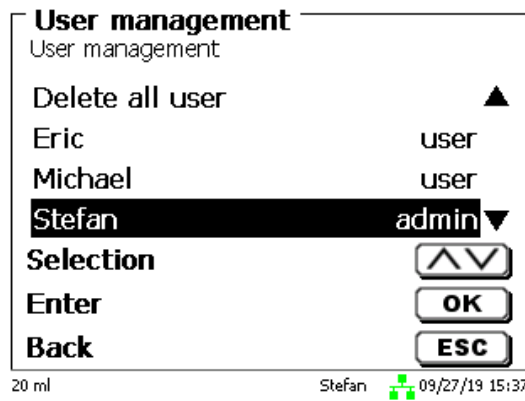
Fig. 122

You can delete all users with «delete all users» (Fig. 123).


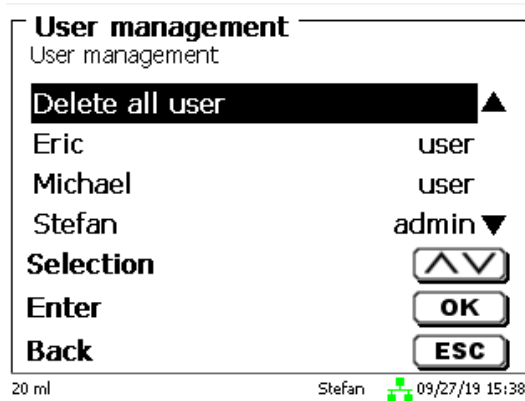
Fig. 123

Confirm with <ENTER>/<OK>.

You have to confirm the delete of all users with «Yes» (Fig. 124).

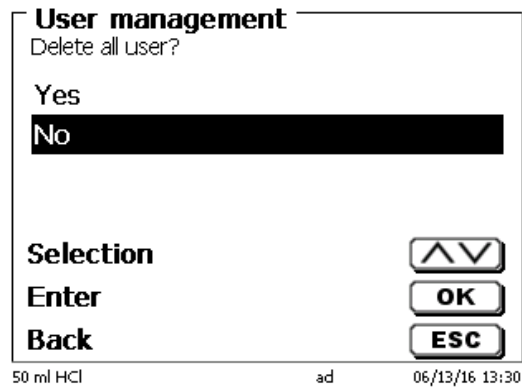**Fig. 124**

At the end only the first Administrator is active (Fig. 125).

**Fig. 125**

You can deactivate and activate the user management if you want easily.  
The first administrator is still there.


**i** Only a RESET will delete the first administrator!

## 5.5 RESET

RESET will reset all settings to the factory setting.

**i** All methods will also be deleted! So please print the methods or export/copy them to a connected USB storage medium (this will be possible with a higher update!).

The RESET has to be confirmed separately once again (Fig. 126).

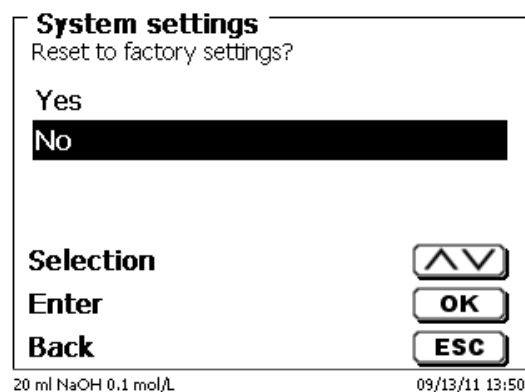**Fig. 126**

## 5.6 Printer

For connecting printers (Fig. 127) please refer to  8.3 Printers.

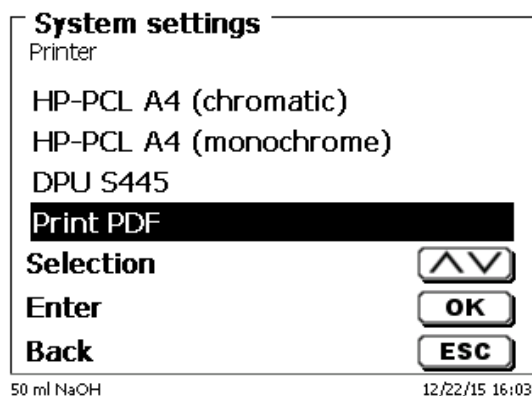
Fig. 127

## 5.7 Device Information

This point contains information about the device (Fig. 128).

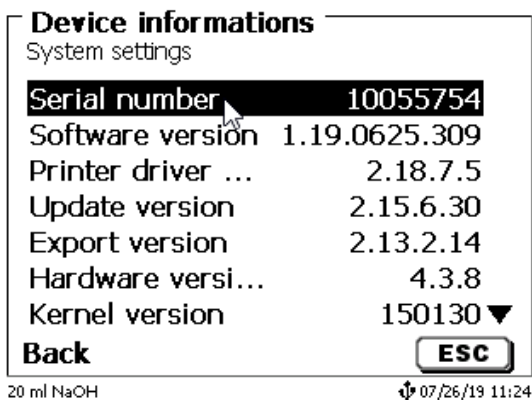
Fig. 128

## 5.8 System Tones

This is the point to set the volume of the system sounds and the front keyboard of the device (Fig. 129). The system sounds become audible e.g. at the end of the titration or in case of an erroneous operation. The keys of the front keyboard produce a clicking sound if the key was used successfully.

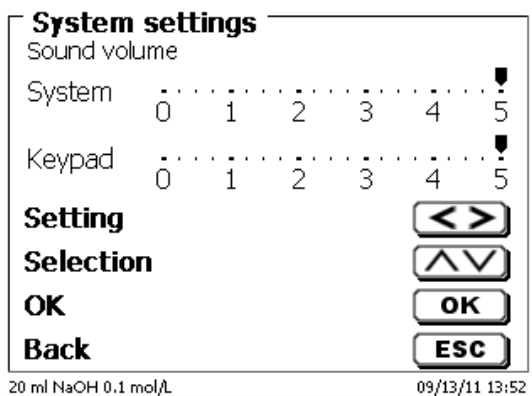
Fig. 129

 No sounds will occur when the external keyboard is used.

## 5.9 Data exchange

All methods with all parameter settings and global memories can be stored and restored on a connected USB-memory. It is also possible to transfer the settings from one titrator to another one. The backup will be started with «**Settings backup**» (Fig. 130).

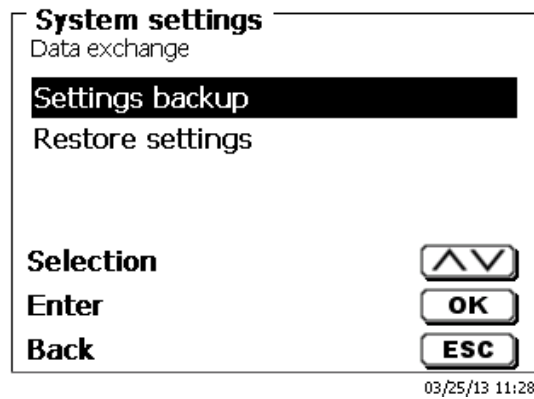
Fig. 130

“Backup settings” is displayed during the backup in blue (Fig. 131).

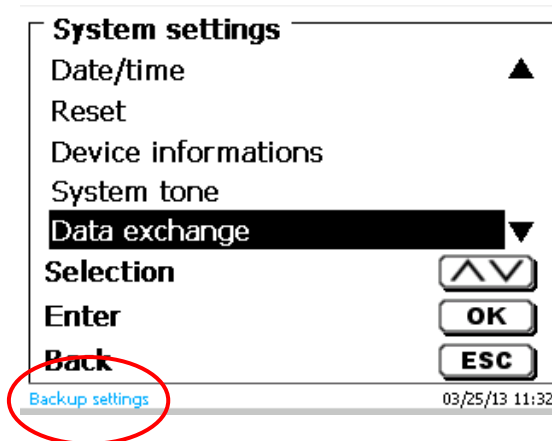
Fig. 131

After a Reset or a maintenance case it is possible to restore the backup with «**Restores settings**» (Fig. 132)

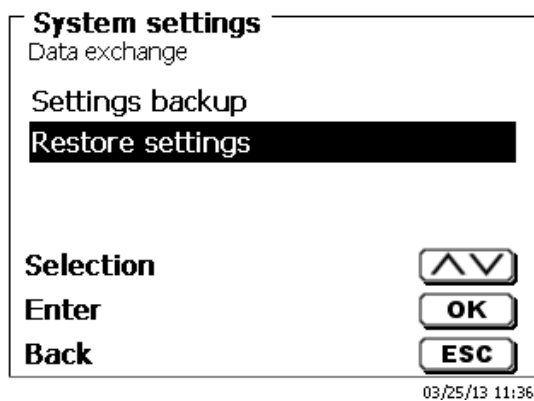
Fig. 132

The backup folder on the USB-memory Stick starts with the backup date (Fig. 133).

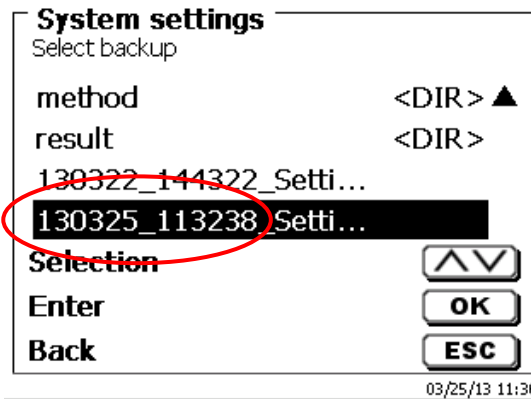
Fig. 133

Confirm the selection with <ENTER>/<OK>.

"Settings are being restored" is displayed during the restoring process of the backup in blue (Fig. 134).



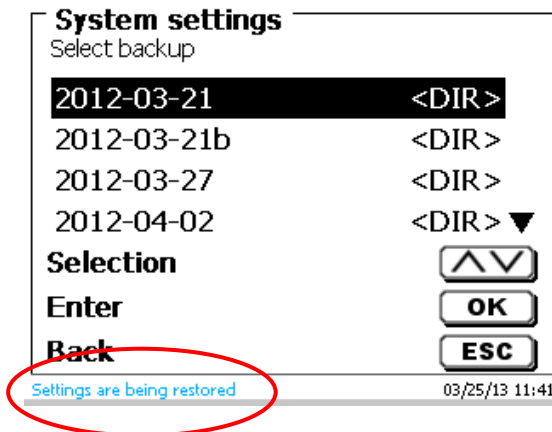
Fig. 134

## 5.10 Software Update

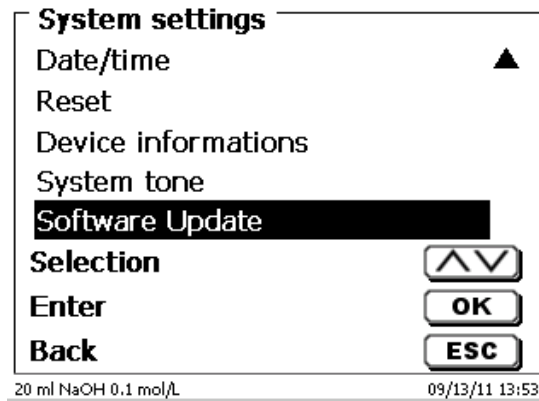**Fig. 135**


An update of the device software (Fig. 135) requires a USB stick containing a new version. For this operation, the two files that are needed have to be located in the root directory of the USB device (Fig. 136).

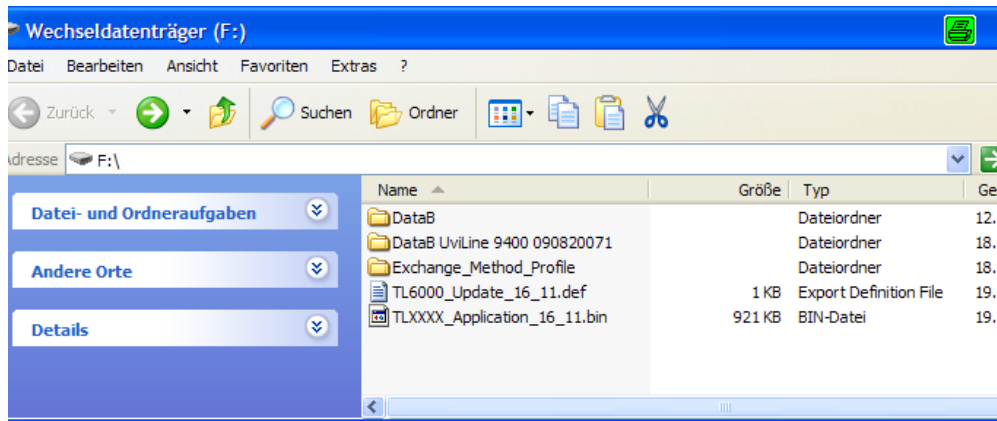**Fig. 136**

Plug the USB device into a free USB-A port, wait for some seconds, and then select the Software Update function. The valid software updates will be shown on the display.

In the present case (Fig. 137) this is Version "15\_50" from week 50 and year 2015.

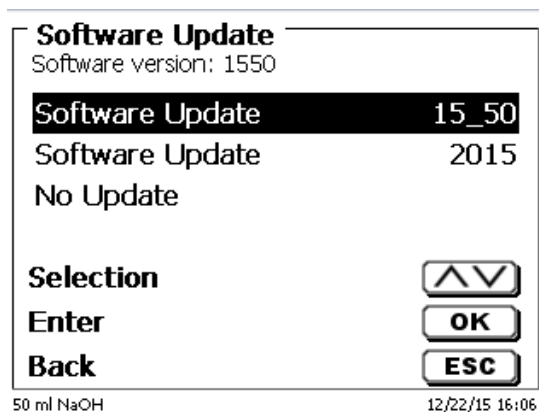**Fig. 137**

After starting the update using <ENTER>/<OK>, next thing to appear is the following graphic (Fig. 138),

## TitroLine® 7500 KF trace

Waiting for system readiness...

Vers.2.15.6.30.20

**Fig. 138**

which will change after a few seconds to the following display (Fig. 139).

## TitroLine® 7500 KF trace

System is updating. Please wait...


Vers.2.15.6.30.20

**Fig. 139**

Upon completion of the update (approx. 4 - 5 minutes), the device will shut down the software completely and proceed to a new start.

**i** In the course of an update, the methods will not be deleted! You can continue to use them.

If no valid update file is stored on the USB stick, a message will appear (Fig. 140)

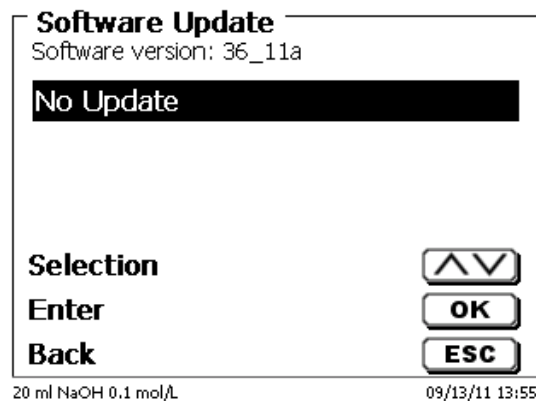**Fig. 140**

## 6 Network settings

### 6.1 General

Via the network/Ethernet interface it is possible to save the results in PDF and CSV -format on shared directories of a network. Instead of saving results to a network directory, you can also set the output on a network printer.

Connect the titrator to your network with a suitable network cable.

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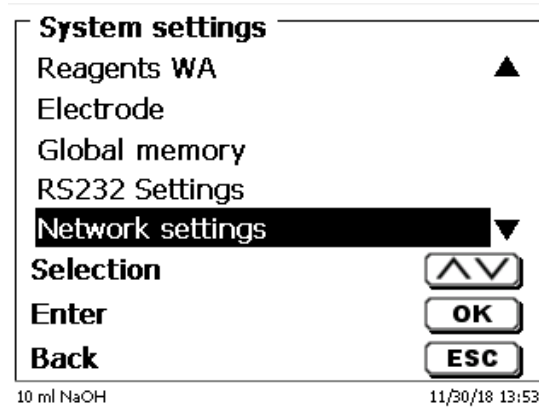
Fig. 141

confirm the selection with <ENTER>/<OK>.

As a rule, the titrator automatically obtains an IP address from the network when DHCP is activated (Fig. 142).

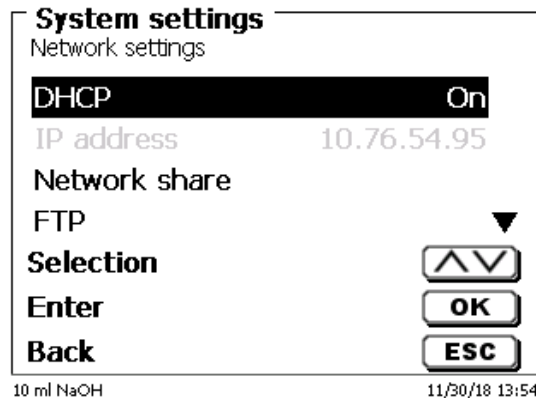
Fig. 142

If DHCP is disabled, you can also enter the relevant network data manually (Fig. 143).

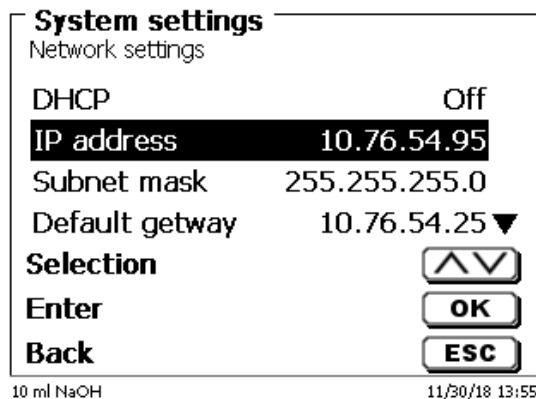
Fig. 143

## 6.2 Setup a shared directory

Select «**Network share**» and confirm your selection with <ENTER>/<OK> (Fig. 144).

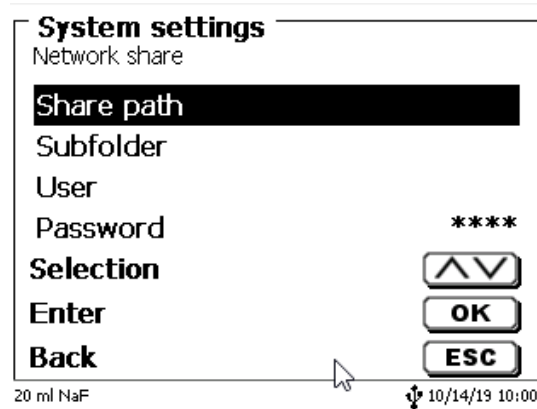
Fig. 144

Enter the «**share path**» (Fig. 145).  
Please ask your IT specialist what exactly this path is.

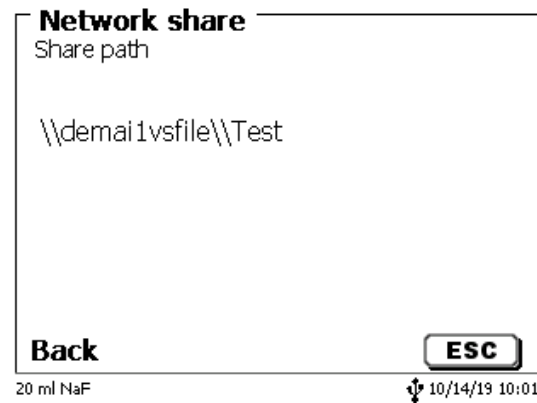
Fig. 145

Complete the entry with <ENTER>/<OK>.  
Now enter your «**Username**» and «**Password**» for your corporate network (Fig. 146).


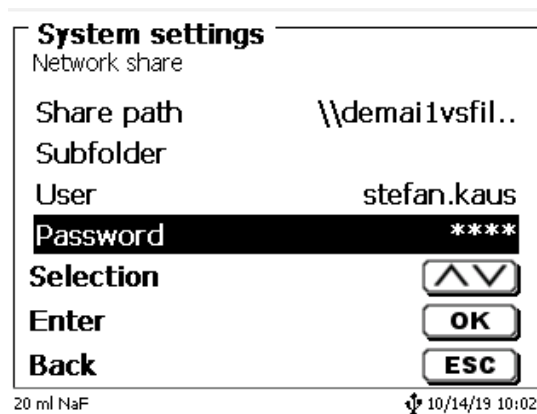
Fig. 146

After leaving the network menu short a window appears with information about the connection to the network.


Under «**User**» and «**Password**» a combination authorized for the folder must be entered. If access is denied or the share can not be reached then this will be displayed when exiting the menu.

Now go back one step with <ESC> to the system settings.  
Go to «**Printer selection**» (Fig. 147).

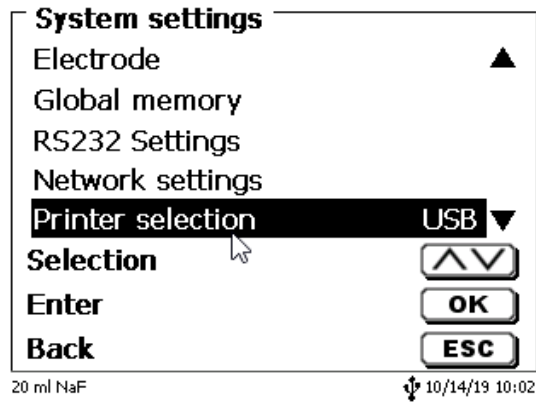**Fig. 147**

And select «**Network share**» (Fig. 148).

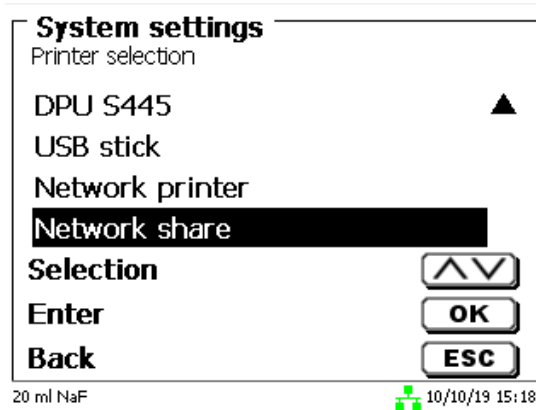**Fig. 148**

PDF and CSV files are now automatically saved on the shared network drive.

**i** You can also select a network printer instead of the network share. The network printer must understand the HP-PCI 3, 4, 5, or 5e printer language.

## 7 Communication via RS-232 and USB-B interface

### 7.1 General Information

The TitroLine® 7500 KF *trace* has two serial RS-232-C interfaces to communicate data with other devices. By means of these two interfaces it is possible to operate several devices on one computer (PC) interface. In addition to that, the TitroLine® 7500 KF *trace* also has an alternatively USB-B interface, which can only be used to connect a PC. RS-232-C-1 establishes the connection to a connected computer or to the previous device of the “Daisy Chain”. At the RS-232-C-2 it is possible to connect additional devices (Daisy Chain Concept).

PIN assignment of the RS-232-C interfaces:

PIN No.	Meaning / Description
1	T x D Data output
2	R x D Data input
3	Digital mass

### 7.2 Chaining multiple devices - “Daisy Chain Concept”

In order to activate several devices in a chain individually, each device must have an own device address. For this it is at first necessary to establish a connection from the computer to the RS-232-C interface 1 of the first device in the chain by means of a RS-232-C data cable, e.g. Type No. TZ 3097. With the additional RS-232-C data cable, Type No. TZ 3094, the RS-232-C- interface 2 of the first device is connected with the RS-232-C-interface 1 of the second device. At interface 2 of the second device it is possible to connect an additional device.

The TitroLine® 7500 KF *trace* can also be connected via USB cable TZ 3840 (type A (M) - type B (M), 1.8m) to a USB interface of a PC. To accomplish this connection, a driver has to be installed on the PC. Then the USB-B interface takes over the function of the RS232-1 interface.

The address always consists of two characters: e.g. address 1 of the two ASCII- characters <0> and <1>. The addresses can be set from **00** to **15**, i.e. 16 possibilities. It must be ensured that the devices in a chain have different addresses. If a device is addressed with its address, this device will process this command without sending it to another device. The reply to the computer has also an own address. The addresses are allocated as described in  **5.2 RS-232 Settings**.

The TitroLine® 7500 KF *trace* receives commands from a PC at the interface **1** (USB- B) if the computer knows the address. It also sends the answer via this interface. If the address of the incoming command does not match the device address, the complete command will be forwarded to interface **2**. Interface 2 is connected to interface 1 of another device. This device checks the address as well and reacts to the command as the first TitroLine® 7500 KF *trace* did before.

All information (data strings) which arrive at interface 2 of the TitroLine® 7500 KF *trace* will immediately be send to the computer via interface 1 (or USB-B interface). Thus, the computer receives the data of all devices. In practice it is possible to connect up to 16 devices to one computer- (PC-) interface.

### 7.3 Instruction Set for RS-Communication

The commands consist of three parts:

Adresse, two-digit aa	e.g. <b>01</b>
Command	e.g. <b>DA</b>
Variable, if necessary	e.g. <b>14</b>
an end of command	<b>&lt;CR&gt; &lt;LF&gt;</b>

 **Every command must be completed with the ASCII - sign <CR> and <LF>** (Carriage Return and Line Feed). Only if the respective action has ended the answers will be returned to the computer.

Example:

The command to dose 12.5 ml shall be sent to the TitroLine® 7500 KF *trace* with the address 2.

The command consists of the characters:

<b>02DA12.5&lt;CR LF&gt;</b>	in detail:
02	= Device address
DA	= Dosage command with filling and zero points of the display
12.5	= Volume in ml to be dosed
<CR LF>	= Control character as command end

Command	Description	Reply
aaAA	automatic allocation of device address	aaY
aaMC1...XX	choosing a method	aaY
aaBF	"filling burette"	aaY
aaBV	output of dosed volume in ml	aa0.200
aaDA	dose volume without filling, with adding the volume	aaY
aaDB	dose volume without filling, reset of the volume	aaY
aaDO	dose volume with filling, without adding the volume	aaY
aaGDM	dosing speed in ml/min	aaY
aaGF	filling time in seconds (min is 20, default 30)	aaY
aaEX	"exit" function.back to main menu	aaY
aafd	$\mu$ A "dead stop" measurement function	aaY
aafp	pH measurement function	aaY
aaft	temperature measurement function	aaY
aafv	mV measurement function	aaY
aaagdm	dosing speed in ml/min (0.01 – 100 ml/min)	aaY
aaGF	filling time in sec (adjustable 20 – 999 seconds)	aaY
aaGS	output serial no. Of device	aaGS08154711
aaLC	output of the CAL parameters	
aaLD	output of the measurement data	aaY
aaLR	output report (short report)	aaY
aaM	output of the preset measurement value (pH/mV/ $\mu$ A)	aaM7.000
aaRH	request of identification	aaIdent: TitroLine® 7500 KF <i>trace</i>
aaRC	send last command	aa"last command"
aaRS	report status possible answers are: "STATUS:READY" for ready "STATUS:dosing" dosing "STATUS:filling" filling "ERROR:busy" if no interchangeable unit has been attached	aaStatus:"text"
aaSM	start selected method	aaY
aaSEEPROM	EEPROM reset to factory defaults	aaY
aaSR	stop the actual function	aaY
aaSS	titration start with the transfer of the pH end value	aaY
aaVE	Version number of the software	aaVersion

## 8 Connection of Analytical Balances and Printers

### 8.1 Connection of Analytical Balances

As it often happens that the sample is weighed in on an analytical balance, it makes sense to connect this balance to the TitroLine® 7500 KF *trace*. The balance must have a RS-232-C-interface and the connection cable must be configured accordingly. For the following types of balances there are already assembled connection cables:

Balance	TZ-Number
Sartorius (all type with 25-pole RS-232), partially Kern	TZ 3092
Mettler, AB-S, AG, PG, Sartorius with USB-Port	TZ 3099
Precisa XT-Serie	TZ 3183
Kern with 9-pole RS-232	TZ 3180

For all other types of balances it is possible to obtain an already assembled connection cable (on demand). For this we need detailed information about the RS-232-C-interface of the balance used.

The connection cable is to be connected to the RS-232-C-interface 2 of the TitroLine® 7500 KF *trace*. This side of the connection cables always consists of a 4-pole mini-plug. The other side of the cable can, depending on the type of balance, be a 25-pole plug (Sartorius), a 9-pole plug (Mettler AB-S) or a 15-pole specialised plug (Mettler AT) etc.

In order to allow the balance data to be sent to the TitroLine® 7500 KF *trace*, the data transmission parameters of the titrator and the balance must correspond to each other. Additionally, it is necessary to carry out some more standard settings on the side of the balances:

- The balance is to send the balance data via RS-232-C only by means of a print command
- The balance is to send the balance data only after the display standstill
- The balance should never be set to “automatic sending” and/or “send continuously”
- “Handshake” on the balance must be set to “off”, or even “Software Handshake” or “Pause”

No special characters such as **S** or **St** are allowed to be used as prefix in the balance data of the balance data string. In such a case it might be possible that the TitroLine® 7500 KF *trace* cannot process the balance data correctly.

After you have connected the balance with the appropriate cable and have adjusted all settings in the balance software, and possibly in the TitroLine® 7500 KF *trace*, you can now test the data transfer of the balance very easily. Start the one method. Confirm the sample designation. Then, the display asks you:

- a) to press the print-button at the balance  
→ Parameters to “weighted sample automatically”
- b) to enter the weighted sample → then the parameters are still set to “weighted sample manually”

Put an object onto the balance and press the print button.

After the standstill of the balance display there will be beep and the transmitted balance data appear:

- a) the display changes automatically into the measuring display.
- b) the weighted sample must again be confirmed with **<ENTER>/<OK>**.

## 8.2 Balance data editor

Pressing «F5/balance symbol» will invoke the so-called balance data editor. A list with the existing balance data will appear (Fig. 149).

List of balance data				
3 Weights				
002	M	10.42980	g	13:59:57
003	M	0.87360	g	14:00:10
004	M	4.37650	g	14:00:21
Selection				<input type="button" value="^v"/>
Enter				<input type="button" value="OK"/>
Back				<input type="button" value="ESC"/>
20 ml NaOH 0.1 mol/L			09/13/11 14:00	

Fig. 149

The balance data can be edited one by one. Following a change, a star will appear opposite the weighed-in quantity (Fig. 150).

List of balance data				
3 Weights				
002	M	10.42980	g	13:59:57
003	*M	0.86360	g	14:00:10
004	M	4.37650	g	14:00:21
Selection				<input type="button" value="^v"/>
Enter				<input type="button" value="OK"/>
Back				<input type="button" value="ESC"/>
20 ml NaOH 0.1 mol/L			09/13/11 14:00	

Fig. 150

Weights may be deleted or added individually. It is also possible to delete all weights at one stroke (Fig. 151).

Balance data	
003 *M 0.86360 g	
Edit weight	
Delete weight	
Add weight	
Delete all?	
Selection	<input type="button" value="^v"/>
Enter	<input type="button" value="OK"/>
Back	<input type="button" value="ESC"/>
20 ml NaOH 0.1 mol/L	
09/13/11 14:01	

Fig. 151

If no balance data is available, the «No balance data found» message will appear (Fig. 152).

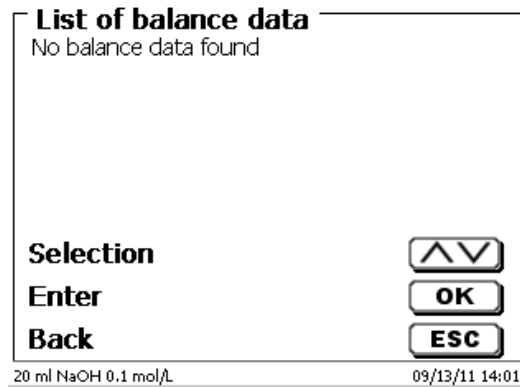
Fig. 152

### 8.3 Printers

The results, calibration data and methods can be printed on the following media

- HP PCL compatible printer (A4)
- Seiko DPU S445 (Thermo paper 112 mm width)
- On the USB stick in PDF- and CSV-format

To connect the printers to the burette please use the USB socket.

When printing, please check whether the correct printer is connected. It is not possible to print “HP” printer layouts on another thermal printer or vice versa. The printer settings should always be checked and adjusted after changing the printer (Fig. 153).

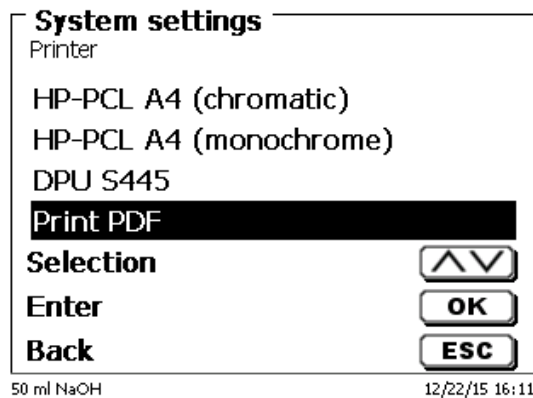
Fig. 153

**i** Only one printer should be connected for one titrator because automatic printer recognition is not activated. «Print PDF» is the default setting.

## 8.4 Automatic stirrer control

### 8.4.1 General

If the magnetic stirrer TM 235 or TM 235 KF is connected via USB, the stirrer can be controlled via the titrator. A suitable connection cable is included with the TM 235 / TM 235 KF.

### 8.4.2 Basic setting in the system menu

Connect the magnetic stirrer with the USB cable to one of the two USB A sockets. Under «**System Settings**», select «**Stirrer Control**» (Fig. 154).

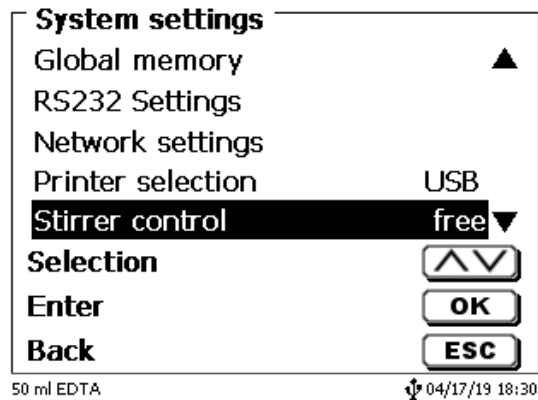
Fig. 154

Confirm the selection with <ENTER>/<OK>. The default setting is set to «free». The stirrer control thus only works with the thumb wheel on the magnetic stirrer (Fig. 155).

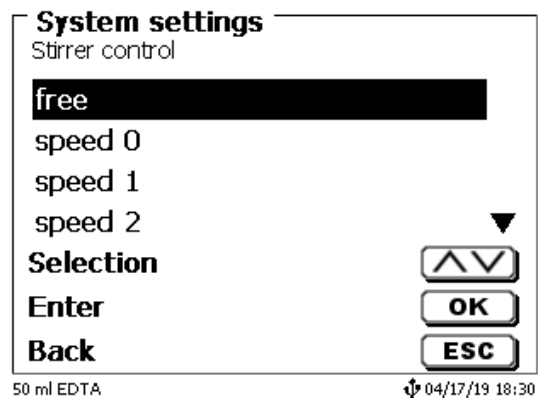
Fig. 155

If you want to deactivate the stirring speed when switching on, you must select the speed «0» level (Fig. 156).

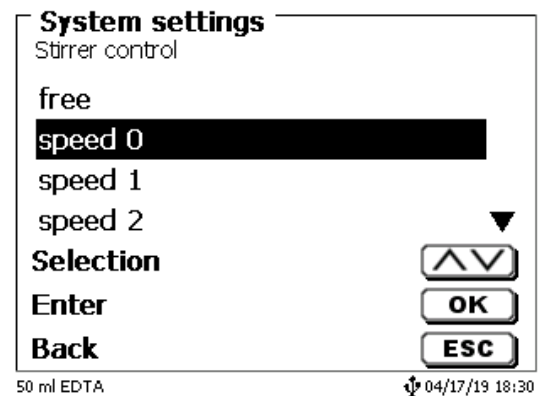
Fig. 156

### 8.4.3 Set the stirring speed in the method

Thereafter, an individual stirring speed in the titration parameters can be set for each method (Fig. 157 and Fig. 158).

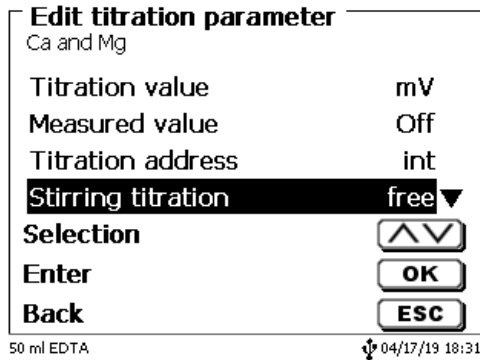
Fig. 157

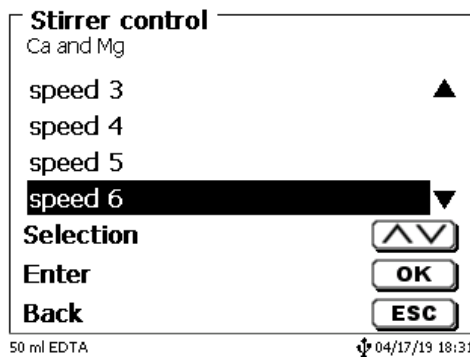
Fig. 158

The stirring speed can also be set individually for the individual pre-dosing steps, the pre-titration step and the following waiting times (Fig. 159 and Fig. 160)

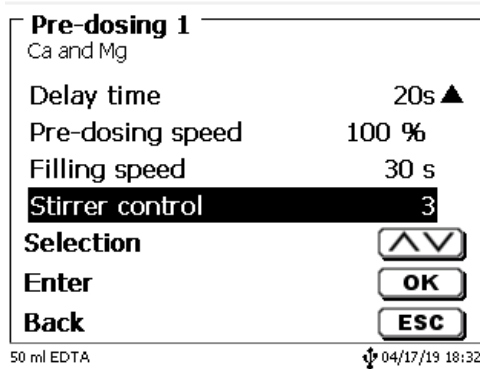
Fig. 159

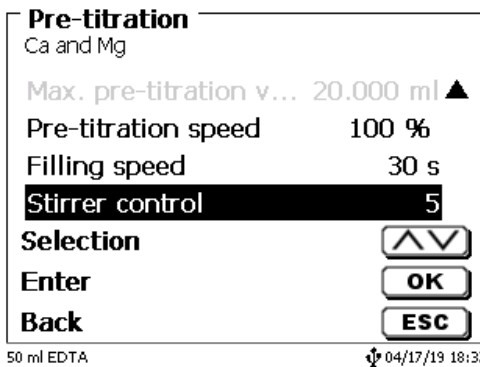
Fig. 160

## 8.5 Autosampler

### 8.5.1 Connection of sample changer TW 7400

The sample changer TW 7400 plus is connected to RS232-2 (RS2) of the titrator by cable **TZ 3987**.

**i** The settings of the RS232-2 interface do not have to be changed. They can remain at 4800, No. 8.1.

## 8.6 Using software TitriSoft

### 8.6.1 General

The titrator is connected to the PC via the RS232 or USB-1-B interface. Cables TZ 3097 and TZ 3091 can be used via RS232-1 for the connection.

### 8.6.2 TitriSoft 3.15 or higher

When using the new software 3.15 or higher TitriSoft, the factory settings of the RS232-1 can be maintained.

Reading and writing the intelligent exchange units and ID electrodes is possible with TitriSoft 3.15. For more information, please refer to the operating manual of TitriSoft.

## 9 Maintenance and Care of the Titrator

 The preservation of the proper functioning of the device requires testing and maintenance work to be performed on a regular basis. Regular inspections are essential prerequisites for the correctness of the volume and the proper functioning.

The accuracy of the volume is determined by all chemicals-carrying components (piston, cylinder, valve, titration tip and hoses). These parts are subject to wear and tear. The piston and cylinder are subject to particular strain, hence they require special attention.

### Heavy strain:

Use of e.g. concentrated solutions, reagents and chemicals (> 0.5 mol/L); chemicals attacking glass, such as fluorides, phosphates, alkali solutions; solutions with a tendency to crystallising out; Fe (III) chloride solutions; oxidising and corroding solutions such as iodine, potassium permanganate, Cer (III), Karl-Fischer titration agent, HCl; solutions with a viscosity of > 5 mm<sup>2</sup>/s; frequent, or even daily use.

### Normal strain:

Use of solutions, reagents and chemicals (up to 0.5 mol/l) which do not attack glass, crystalize out or corrode.

### Interrupted use:

If the dosing system is not in use for more than two weeks, we recommend emptying and cleaning the dosing unit [6]. This applies in particular under the operating conditions referred to in the „Heavy strain” section. If this recommendation is not adhered to, the piston of the valve may become leaking, this may result in damage to the piston burette.

 If the liquid is left within the system, you will also have to reckon with corrosion and an alteration of the solutions used over time, which includes e.g. crystallisation. Considering that as of the state of the art there are no plastic hoses available for the use in titration equipment which would be perfectly free of diffusion phenomena, particular attention is to be paid to the range of the hose lines.

### We recommend the following inspection and maintenance work:

	Heavy strain	Normal strain
Simple cleaning: <ul style="list-style-type: none"> <li>Wiping off splashed chemicals from the outer surface [1]</li> </ul>	Whenever required in operation	Whenever required in operation
Sight check: <ul style="list-style-type: none"> <li>Check for leakage in the area of the dosing system. [2]</li> <li>Is the piston tight? [3]</li> <li>Is the valve tight? [4]</li> <li>Titration to clear? [5]</li> </ul>	Weekly, when putting back into operation	Monthly, when putting back into operation
Basic cleaning of the dosing system: <ul style="list-style-type: none"> <li>All parts of the dosing system to be cleaned separately. [6]</li> </ul>	Every three months	Whenever necessary
Technical inspection: <ul style="list-style-type: none"> <li>Check for air bubbles in the dosing system. [7]</li> <li>Visual inspection</li> <li>Check of the electrical connections. [8]</li> </ul>	Semi-annually, when putting back into operation	Semi-annually, when putting back into operation
Verification of the volume according to ISO 8655 <ul style="list-style-type: none"> <li>Perform basic cleaning</li> <li>Inspection according to ISO 8655 Part 6 or Part 7. [9]</li> </ul>	Semi-annually	Annually

 Depending on the respective application, there may be different specifications for the entirety of the inspection and maintenance work to be performed. The individual intervals may be extended if no complaints occur, but they will have to be shortened again as soon as any problem has arisen

The inspection of the metrological reliability including maintenance work is offered as a service (including a manufacturer's certificate, if so ordered). In this case the titration device is to be sent in. Please contact the service (see backside of this manual).

## Detailed description of the inspection and maintenance work

- [1] Wipe off using a soft cloth (and some water with a normal household detergent).
- [2] Leaking connections can be identified by moisture or crystals at the threaded connections of the hoses, at the sealing lips of the piston inside the dosing cylinder or at the valve.
- [3] If any liquid becomes visible below the first sealing lip, it has to be checked at short timely intervals whether any liquid will build up under the second sealing lip, too. In this case both the piston and the glass cylinder have to be replaced immediately. It is easily possible that in operation small liquid droplets build up under the first sealing lip, but they may also disappear again. This phenomenon alone is no reason for replacement.
- [4] The valve has to be removed from its housing for inspection. In this process, the hoses remain connected to the valve. Please check for moisture underneath the valve. When reinserting the valve, please make sure that the small cam at the rotating axis is fitted into the corresponding groove again.
- [5] The titration tip must be free of sedimentation or crystals which might obstruct the dosing process or falsify the results.
- [6] Remove the cylinder, take the valve out of the valve housing, unscrew the hoses and then rinse all parts carefully with distilled water. For the assembly of the cylinder, hoses and other parts of the interchangeable unit, please refer to the operating instructions.
- [7] Dose one burette volume, then refill. Air bubbles will gather at the tip of the cylinder and in the titration hose where they can be detected easily. If bubbles become visible, please re-tighten all connections finger tight, and then repeat dosing. If air bubbles still remain within the system, [6] please check the valve and replace the hose connections. The air bubbles may also occur at the interface between the sealing lip of the piston and the cylinder. If a reduction of the filling speed will not do, the dosing unit has to be replaced.
- [8] Check the electrical plug contacts for corrosion and mechanical damage. Defective parts have to be repaired or replaced by new parts.
- [9] Please refer to the application "Burette inspection according to ISO 8655 Part 6".

## 10 Guarantee

We provide guarantee for the device described for two years from the date of purchase. This guarantee covers manufacturing faults being discovered within the mentioned period of two years. Claim under guarantee covers only the restoration of functionality, not any further claim for damages or financial loss. Improper handling/use or illegitimate opening of the device results in loss of the guarantee rights. The guarantee does not cover wear parts, as lobes, cylinders, valves and pipes including the thread connections and the titration tips. The breach of glass parts is also excluded. To ascertain the guarantee liability, please return the instrument and proof of purchase together with the date of purchase freight paid or prepaid.

## 11 Storage and transportation

If the TitroLine® 7500 KF *trace* or the interchangeable units have to be stored over some time, or to be dislocated, the use of the original packing will be the best protection of the devices. However, in many cases this packing will not be available anymore, so that one will have to compose an equivalent packaging system. Sealing the lower section in a foil is hereby recommended. The devices should be stored in a room with a temperature between +10 and +40°C, and the (relative) humidity of the air should not exceed 70 %.

 If the interchangeable have to be stored over some time, or to be dislocated, the fluids inside the system, especially aggressive solution have to be removed.

## 12 Recycling and Disposal

Please observe the applicable local or national regulations concerning the disposal of "waste electrical and electronic equipment".

The TitroLine® 7500 KF *trace* and his packaging are manufactured as far as possible from materials which can be disposed of environmental-friendly and recycled in a technically appropriate manner. If you have any question regarding disposal, please contact the service (see backside of this manual).

 The main printed board carries a lithium battery (type CR 2430). Batteries should not to be disposed of with the normal domestic waste. They will be taken back and recycled or disposed of properly by the manufacturer at no cost.

### **13 EC – Declaration of Conformity**

The corresponding declaration of conformity of the device can be found on our homepage. It will also be made available to you on request.

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SI Analytics is a trademark of Xylem Inc. or one of its subsidiaries.
**Xylem Analytics Germany Sales GmbH & Co.KG**
© 2022 Xylem, Inc. Version 220407 M 827 231 5