**BEDIENUNGSSANLEITUNG**  
Originalversion  
**OPERATING MANUAL**  
**MODE D'EMPLOI**  
**MANUAL DE INSTRUCCIONES**

---

# ScienceLine IDS<sup>®</sup>

---

**SI Analytics**  
a xylem brand

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## **Operating Manual ..... Page 11 ... 18**

### **Important notes:**

The operating manual is part of the product. Before initial operation, please carefully read and observe the operating manual and keep it. For safety reasons the product may only be used for the purposes described in these present operating manual. Please also consider the operating manuals for the devices to be connected.

All specifications in this operating manual are guidance values which are valid at the time of printing. However, for technical or commercial reasons or in the necessity to comply with the statutory stipulations of various countries, the manufacturer may perform additions to the product without changing the described properties. A potentially more recent version of this manual is available on our internet website. The German version is the original version and binding in all specifications!

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## TABLE OF CONTENT

<b>1</b>	<b>General information, Technical Data</b>	<b>13</b>
1.1	General information	13
1.2	Technical Data	13
1.2.1	General data	13
1.2.2	Measurement and application characteristics	13
1.2.3	Shaft dimensions, material, electrical connection	13
1.2.4	Connection cable	14
1.2.5	Accuracy of the IDS <sup>®</sup> measuring technique	14
<b>2</b>	<b>Commissioning, calibration and measuring</b>	<b>14</b>
2.1	Commissioning	14
2.2	Calibration and measurement	15
2.2.1	General rules	15
2.2.2	Subsequent calibrations	15
<b>3</b>	<b>Storage</b>	<b>16</b>
3.1	During short measuring breaks	16
3.2	Overnight or longer	16
<b>4</b>	<b>Aging</b>	<b>16</b>
<b>5</b>	<b>Maintenance and cleaning</b>	<b>17</b>
5.1	Cleaning	17
5.2	After cleaning	17
<b>6</b>	<b>Wear parts and accessories</b>	<b>18</b>

### Copyright

© 2021, Xylem Analytics Germany GmbH

Reprinting - even as excerpts - is only allowed with the explicit written authorization.

Germany. Printed in Germany.

# 1 General information, Technical Data

## 1.1 General information

The sensor electronics with the stored sensor data is in the connecting head of the electrode. The data include, among other things, the sensor type and series number. With each calibration, the calibration data is written in the sensor and the calibration history is recorded. The data is recalled by the meter when the sensor is connected and is used for measurement and for measured value documentation. Storing the calibration data in the sensor ensures that the correct slope and asymmetry are automatically used if the sensor is operated with several meters.

The digital transmission technique guarantees the failure-free communication with the meter even with long connection cables. If the sensor firmware is enhanced by SI Analytics<sup>®</sup>, it can be updated via the meter.

## 1.2 Technical Data

### 1.2.1 General data

SI Analytics <sup>®</sup> Model	Reference electrolyte	Junction	NTC	Special features
ScienceLine A6880 IDS <sup>®</sup>	KCl 3 mol/l, Ag <sup>+</sup> -free	Ceramic	Yes	for cut-in measurement
ScienceLine A157 IDS <sup>®</sup>	KCl 3 mol/l, Ag <sup>+</sup> -free	Platinum	Yes	Micro
ScienceLine A161/A162 IDS <sup>®</sup>	KCl 3 mol/l, Ag <sup>+</sup> -free	Platinum	Yes	Glass shaft
ScienceLine A7780 IDS <sup>®</sup>	Gel	Ceramic	Yes	Triple ceramic

### 1.2.2 Measurement and application characteristics

SI Analytics <sup>®</sup> Model	pH measuring range	Allowed temperature range	Membrane resistance at 25 °C	Typical application
ScienceLine A6880 IDS <sup>®</sup>	0 ... 14	- 5 ... 80 °C	< 400 MOhm	Laboratory/Food
ScienceLine A157 IDS <sup>®</sup>	0 ... 14	- 5 ... 100 °C	< 500 MOhm	Laboratory
ScienceLine A161/A162 IDS <sup>®</sup>	0 ... 14	- 5 ... 100 °C	< 400 GOhm	Laboratory
ScienceLine A7780 IDS <sup>®</sup>	0 ... 14	- 5 ... 100 °C	< 400 GOhm	Laboratory/Field

### 1.2.3 Shaft dimensions, material, electrical connection

SI Analytics <sup>®</sup> Model	Shaft			Electrical connection		
	Length [mm]	Ø [mm]	Material	Combination electrode connection	Meter connection	Cable length
ScienceLine A6880 IDS <sup>®</sup>	70/50	12/8	Glass	Fixed cable	Digital plug	1.5 m
ScienceLine A157 IDS <sup>®</sup>	70/130	12/5	Glass	Fixed cable	Digital plug	1.5 m
ScienceLine A161/A162 IDS <sup>®</sup>	120 (A162) 170 (A161)	12	Glass	Fixed cable	Digital plug	1.5 m
ScienceLine A7780 IDS <sup>®</sup>	120	12	Glass	Fixed cable	Digital plug	1.5 m

### 1.2.4 Connection cable

<b>Diameter</b>	4,3 mm
<b>Smallest allowed bend radius</b>	Fixed installation: 20 mm Flexible use: 60 mm
<b>Plug type</b>	Socket, 4 pins

### 1.2.5 Accuracy of the IDS<sup>®</sup> measuring technique

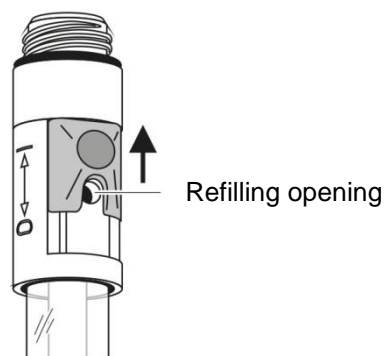
Measured parameter	Accuracy ( $\pm 1$ Digit)
pH	$\pm 0.004$
U [mV]	$\pm 0.2$
T [°C]	$\pm 0.1$

## 2 Commissioning, calibration and measuring

### 2.1 Commissioning

Prepare the electrode for measuring as follows:

- Remove the watering cap from the electrode tip. Possible salt deposits in the area of the watering cap do not affect the measuring characteristics and can easily be removed with deionized water.
- Open the refilling opening for the reference electrolyte solution (except ScienceLine A7780 IDS<sup>®</sup>)

**i** The refilling opening must always be open during calibration and measurement!

**i** Please keep the watering cap. It is required for the electrode to be stored. Always keep the watering cap clean.

## 2.2 Calibration and measurement

### 2.2.1 General rules

**i** Please keep the watering cap. It is required for the electrode to be stored. Always keep the watering cap clean.

- Remove any gas bubbles behind the pH membrane by shaking.
- Connect the electrode to the meter.
- Calibrate the electrode according to the operating manual of the meter and observe the following rules while doing so:
  - Avoid the carryover of any solution (sample or buffer solution) from one measurement to the next by taking the following measures:
    - Shortly rinse the calibration and sample beakers with the solution the beakers are to be filled with next.
    - Between measurements, rinse the electrode with the solution that follows. Alternatively, you can also rinse the electrode with deionized water and then carefully dab it dry.
  - To measure in aqueous solutions, it is recommended to immerse the electrode in a vertical or slightly tilted position.
  - Observe the correct depth of immersion and make sure the contact between the junction and test sample is thorough. The junction is in the area of the bottom end of the shaft (see arrow):

**Caution: Only the shaft part of the combination electrode may be immersed!**

- For measurement in aqueous solutions, provide approximately the same stirring conditions for measuring as for calibrating.

### 2.2.2 Subsequent calibrations

The frequency of subsequent calibrations depends on the application. Many meters provide an option where you can enter a calibration interval. After the calibration interval has expired, the meter will automatically remind you of the due calibration.

## 3 Storage

### 3.1 During short measuring breaks

Immerse the electrode (except ScienceLine A7780 IDS®) in reference electrolyte (KCl 3 mol/l, Ag<sup>+</sup>-free). Prior to the next measurement, shortly rinse the electrode with the test sample or deionized water.

**i** Do not scratch the pH membrane.

### 3.2 Overnight or longer

Put the clean electrode in the watering cap that is filled with reference electrolyte (KCl 3 mol/l, Ag<sup>+</sup>-free).

**i** pH electrodes must not be stored dry or in deionized water. The electrode could be permanently damaged by this. If the liquid in the watering cap has dried up, condition the electrode in reference electrolyte (KCl 3 mol/l, Ag<sup>+</sup>-free) for at least 24 hours.

**i** During longer storing periods, salt sediments may develop on the watering cap. They do not affect the measuring characteristics and can easily be removed with deionized water when the electrode is put into operation again.

## 4 Aging

Every pH electrode undergoes a natural aging process. With aging, the responding behavior becomes slower and the electrode slope and asymmetry change. Moreover, extreme operating conditions can considerably shorten the lifetime of the electrode. These are:

- Strong acids or lyes, hydrofluoric acid, organic solvents, oils, fats, bromides, sulfides, iodides, proteins
- High temperatures
- High changes in pH and temperature.

The warranty does not cover failure caused by measuring conditions and mechanical damage.

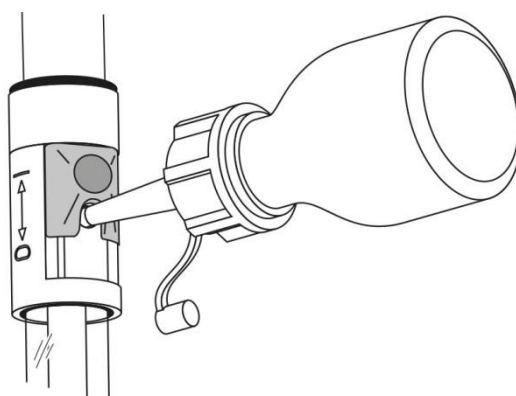
## 5 Maintenance and cleaning

### 5.1 Cleaning

During operation, a small amount of reference electrolyte leaks through the junction from the combination electrode into the test sample. If the level of reference electrolyte becomes too low with time, refill it (except ScienceLine A7780 IDS<sup>®</sup>) through the refilling opening. Refilling is very easy with the enclosed dropping bottle. Proceed as follows:

- Cut off the tip of the dropping bottle at a right angle until the opening in the tip can be seen
- Open the refilling opening of the combination electrode
- Press the tip of the dropping bottle into the refilling opening while turning it slightly
- Pump the reference electrolyte in the shaft using the dropping bottle
- Pull the dropping bottle out of the refilling opening while turning it slightly if necessary.

- Remove water-soluble contamination by rinsing with deionized water. Other types of contamination have to be removed as follows while the contact time with the detergents should be kept as short as possible:

Contamination	Cleaning procedure
Fat and oil	Rinse with water containing household washing-up liquid
Lime and hydroxide deposits	Rinse with citric acid (10 % by weight)
Proteins	Immerse in pepsin cleaning solution PEP/pH for approx. 1 hour. <b>Note:</b> Make sure the level of the reference electrolyte is above that of the cleaning solution.

**i** Hydrofluoric acid, hot phosphoric acid and strong alkaline solutions destroy the glass membrane.

### 5.2 After cleaning

Rinse the electrode with deionized water and condition it in reference electrolyte solution for at least 1 hour. Then recalibrate the electrode.

## 6 Wear parts and accessories

Description	Model	Order no.
Reference electrolyte solution 50 ml (KCl 3 mol/l, Ag <sup>+</sup> - free)	L3008	285138505
Pepsin cleaning solution	L5104	285138295

**i** Information on our wide range of buffer solutions and more accessories is given in the Meters and Electrodes catalog.

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SI Analytics is a trademark of Xylem Inc. or one of its subsidiaries.
### **Xylem Analytics Germany Sales GmbH & Co.KG**
© 2021 Xylem, Inc. Version 211008 **M T814254**