## QUICK START GUIDE

ba77037y02 06/2019

# HandyLab 100

pH METER

**SI Analytics**  
a xylem brand

**Copyright**

© 2019, SI Analytics GmbH

Reprinting - even as excerpts - is only allowed with the explicit written authorization of SI Analytics GmbH, Mainz.

Printed in Germany.

## HandyLab 100 - Contents

EN English .....	2
DE Deutsch.....	4
CS Český .....	6
DA Dansk .....	8
EL Ελληνικά .....	10
ES Español.....	12
FI Suomi .....	14
FR Français .....	16
HU Magyar .....	18
IT Italiano .....	20
JA 日本語 .....	22
KO 한국어.....	24
NL Nederlands .....	26
NO Norsk .....	28
PL Polski.....	30
PT Português .....	32
RO Română .....	34
RU Русский .....	36
SK Slovenský .....	38
SV Svenska .....	40
TR Türkçe .....	42
ZH 中文.....	44

## HandyLab 100

### Overview


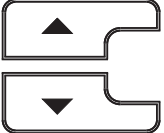
The HandyLab 100 meter enables you to perform measurements (pH) quickly and reliably.

The HandyLab 100 meter provides a maximum degree of operating comfort, reliability and measuring certainty.

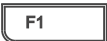
### Keypad

Keys are indicated by brackets <..>.


The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec). A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER\_\_>).

	<F1>: <F1__>:	Situation dependent function (is shown on the display) e.g.: <F1>/[Menu]: Opens the menu for measurement settings <F1__>/[Menu]: Opens the menu for system settings
	<On/Off>:	Switches the meter on or off
	<M>:	Selects the measured parameter
	<CAL>: <CAL__>:	Calls up the calibration procedure Displays the calibration data
	<STO>:	Saves a measured value manually
	<RCL>:	Displays the manually stored measured values
	<▲><▼>:	Menu control, navigation
	<ENTER>: <ENTER__>:	Opens the menu for measurement settings / confirms entries Opens the menu for system settings
	<AR>:	Freezes the measured value (HOLD function) Switches the AutoRead measurement on or off

## Display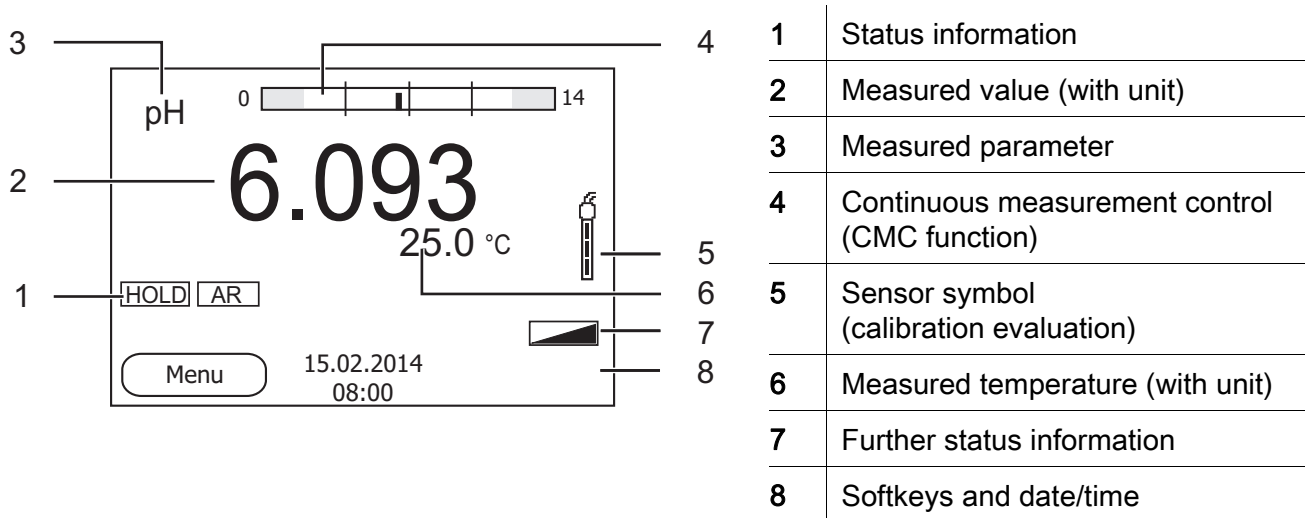

## Jack field /connections

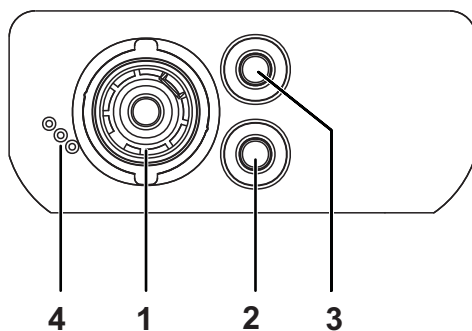1	pH electrode
2	Reference electrode
3	Temperature sensor
4	Service interface

## Initial commissioning

- 1 Insert the supplied batteries. Make sure that the poles are in the correct position.
- 2 Switch on the meter with the <On/Off> key.

### Caution

Follow the safety instructions of the used sensors. The operating manuals of the sensors are available on the supplied CD.

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# Xylem |'zīləm|

- 1) The tissue in plants that brings water upward from the roots;
- 2) a leading global water technology company.

We're a global team unified in a common purpose: creating advanced technology solutions to the world's water challenges. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. Our products and services move, treat, analyze, monitor and return water to the environment, in public utility, industrial, residential and commercial building services settings. Xylem also provides a leading portfolio of smart metering, network technologies and advanced analytics solutions for water, electric and gas utilities. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise with a strong focus on developing comprehensive, sustainable solutions.

**For more information on how Xylem can help you, go to [www.xylem.com](http://www.xylem.com).**

## SI Analytics

a xylem brand

### Hersteller

(Manufacturer)

**Xylem Analytics Germany GmbH**

Am Achalaich 11  
82362 Weilheim  
Germany

SI Analytics

Tel. +49.(0)6131.66.5111

Fax. +49.(0)6131.66.5001

E-Mail: [si-analytics@xylem.com](mailto:si-analytics@xylem.com)

[www.xylemanalytics.com](http://www.xylemanalytics.com)

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(Service and Returns)

**Xylem Analytics Germany Sales GmbH & Co. KG**

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Fax. +49.(0)881.183.414

E-Mail: [wtw.rma@xylem.com](mailto:wtw.rma@xylem.com)

SI Analytics is a trademark of Xylem Inc. or one of its subsidiaries.  
© 2021 Xylem, Inc.