QUICK START GUIDE

ba75783y03 05/2017

Cond 3110

- EN** The latest version of the detailed operating manual is available on the Internet under www.wtw.com.
- DE** Die aktuelle Version der ausführlichen Bedienungsanleitung finden Sie im Internet unter www.wtw.com.
- FR** La version actuelle du mode d'emploi détaillé est disponible sur Internet à l'adresse www.wtw.com.
- ES** La versión actual de la manual de instrucciones detallado lo encuentra Ud. en el internet (www.wtw.com).

Copyright © 2017, Xylem Analytics Germany GmbH
Printed in Germany.

Contents

EN English	4
DE Deutsch.....	8
CS Český	12
DA Dansk	14
EL Ελληνικά	16
ES Español.....	18
FI Suomi	22
FR Français	24
HU Magyar	28
IT Italiano	30
JA 日本語	34
KO 한국어.....	36
NL Nederlands	38
NO Norsk	40
PL Polski.....	42
PT Português	44
RO Română	48
RU Русский	50
SK Slovensky	52
SV Svenska	54
TR Türkçe	56
ZH 中文.....	58


Conductivity meter, Cond 3110

Overview

The Cond 3110 compact precision conductivity meter enables you to perform conductivity measurements quickly and reliably. The Cond 3110 handheld meter provides the maximum degree of operating comfort, reliability and measuring certainty for all applications.


Keypad

Keys are indicated by brackets <.>. The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec) in this operating manual. A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER_>).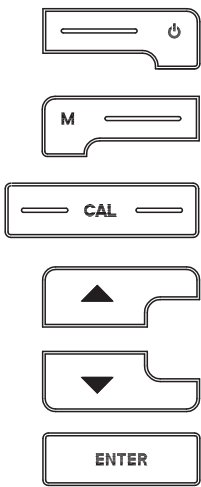

<On/Off>:	Switches the meter on/off
<On/Off_>:	Resets calibration data
<M>:	Selects the measured parameter
<M_>:	Opens the setting menu for calibration and measurements
<CAL>:	Calls up the calibration procedure
<CAL_>:	Displays the calibration data
<>:	Increments values, scrolls
<>:	Decrements values, scrolls
<ENTER>:	Confirms entries
<ENTER_>:	Opens the setting menu for system settings


Display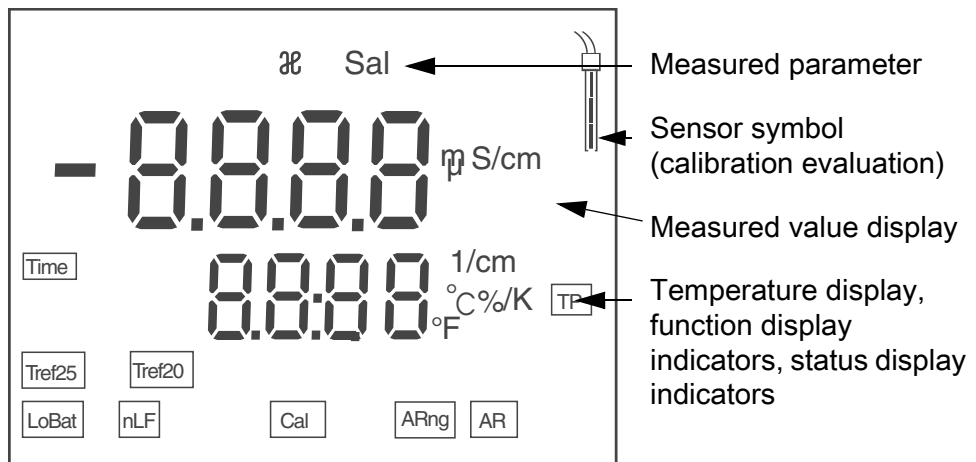

Status display indicators

AR	Stability control (AutoRead) is active
ARng	Automatic range switchover; meter measures with highest possible resolution
Cal	Calibration
LoBat	With battery operation: Batteries almost empty
nLF	Nonlinear temperature compensation
TP	Temperature measurement active
Tref20	Reference temperature of 20 °C
TRef25	Reference temperature of 25 °C
TIME	Setting of calibration interval

Socket field connections

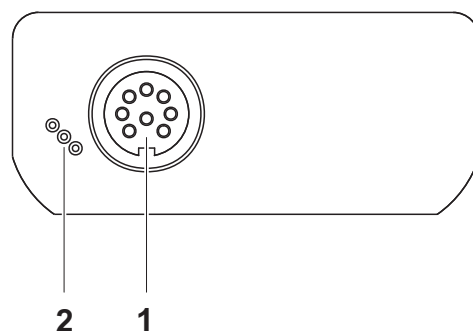1	Conductivity measuring cell
2	Service interface

Initial commissioning

1	Insert the supplied batteries. When doing so make sure that the poles of the batteries are positioned correctly.
2	Switch on the meter with the <On/Off> key.

**Caution**

Follow the safety instructions of the used sensors.
The operating manuals of the sensors are available on the supplied CD and on the Internet under www.WTW.com.

Determining/ setting up the cell constant [C]

Due to aging, the cell constant slightly changes. As a result, an inexact measured value is displayed. Calibration determines the current value of the cell constant and stores this value in the meter.
Thus, you should calibrate at regular intervals.

You can either determine the cell constant of the conductivity measuring cell in the range $0.800 \dots 0.880 \text{ cm}^{-1}$ by calibration in the control standard 0.01 mol/l KCl or adjust it manually in the range $0.800 \dots 1.200 \text{ cm}^{-1}$.

Determining the cell constant (calibration in control standard)

1	Press <CAL> repeatedly until <i>CAL CELL</i> is displayed.
2	Press <ENTER> or <CAL__> to confirm the selection of <i>CAL CELL</i> . The cell constant of the last calibration is displayed.
3	Immerse the measuring cell in the control standard solution, 0.01 mol/l KCl .
4	Start the calibration with <ENTER> . The determination of the cell constant with stability control starts. The [Auto] display flashes until a stable signal is available. The cell constant determined is displayed. The meter automatically stores the cell constant.
5	Switch to the measuring mode with <ENTER> . The determined cell constant is used.

Using the last calibrated cell constant

Prerequisite: No valid calibration exists.

1	Press <CAL> repeatedly until <i>USE CELL</i> is displayed.
2	Press <ENTER> or <CAL__> to confirm the selection of <i>USE CELL</i> .
3	If necessary, press <CAL> repeatedly until <i>CAL</i> and the last calibrated cell constant is displayed.
4	Confirm the selection with <ENTER> . The displayed cell constant is used. The meter switches to the measured value display.

Setting the cell constant manually

The cell constant to be set must either be taken from the operating manual of the measuring cell or is printed on the measuring cell.

1	Press <CAL> repeatedly until <i>USE CELL</i> is displayed.
2	Confirm the selection with <ENTER> or <CAL__> . The cell constant that was set last is displayed.
3	If necessary, press <CAL> repeatedly until a cell constant in the range 0.800 ... 1.200 cm ⁻¹ is displayed.
4	Set the cell constant to be used with <>< > , e.g. 0.846 cm ⁻¹ .
5	Confirm the selection with <ENTER> . The new cell constant is used from now on. The meter switches to the measured value display.

Measuring the conductivity

6	Immerse the conductivity measuring cell in the test sample.
7	If necessary, scroll with <M> until the measured parameter \mathcal{X} with the unit mS/cm or μ S/cm is displayed.
8	Wait for a stable measured value. The [Auto] display indicator flashes as long as the measured value is not yet stable.

Calibration interval (*Int.C*)

The calibration interval reminds you to calibrate regularly. After the specified calibration interval (*Int.C*) has expired, the symbol for the calibration evaluation flashes. It is still possible to measure.

1	Open the menu for measurement settings with <M__> .
2	Confirm all settings with <ENTER> until <i>Int.C</i> is displayed.
3	Set the calibration interval with <>< > .
4	Confirm the setting with <ENTER> .

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Xylem | 'zīləm|

- 1) The tissue in plants that brings water upward from the roots;
- 2) a leading global water technology company.

We're a global team unified in a common purpose: creating advanced technology solutions to the world's water challenges. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. Our products and services move, treat, analyze, monitor and return water to the environment, in public utility, industrial, residential and commercial building services settings. Xylem also provides a leading portfolio of smart metering, network technologies and advanced analytics solutions for water, electric and gas utilities. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise with a strong focus on developing comprehensive, sustainable solutions.

For more information on how Xylem can help you, go to www.xylem.com.

Service and Returns:

Xylem Analytics Germany
Sales GmbH & Co. KG
WTW
Am Achalaich 11
82362 Weilheim
Germany

Tel.: +49 881 183-325
Fax: +49 881 183-414
E-Mail wtw.rma@xylem.com
Internet: www.xylemanalytics.com

Xylem Analytics Germany GmbH
Am Achalaich 11
82362 Weilheim
Germany