## QUICK START GUIDE

ba75787y03 05/2017

# pH 3310

a xylem brand

- EN** The latest version of the detailed operating manual is available on the Internet under [www.wtw.com](http://www.wtw.com).
- DE** Die aktuelle Version der ausführlichen Bedienungsanleitung finden Sie im Internet unter [www.wtw.com](http://www.wtw.com).
- FR** La version actuelle du mode d'emploi détaillé est disponible sur Internet à l'adresse [www.wtw.com](http://www.wtw.com).
- ES** La versión actual de la manual de instrucciones detallado lo encuentra Ud. en el internet ([www.wtw.com](http://www.wtw.com)).

**Copyright** © 2017, Xylem Analytics Germany GmbH  
Printed in Germany.

## Contents

EN English .....	4
DE Deutsch .....	6
CS Český .....	8
DA Dansk .....	10
EL Ελληνικά .....	12
ES Español .....	14
FI Suomi .....	16
FR Français .....	18
HU Magyar .....	20
IT Italiano .....	22
JA 日本語 .....	24
KO 한국어 .....	26
NL Nederlands .....	28
NO Norsk .....	30
PL Polski .....	32
PT Português .....	34
RO Română .....	36
RU Русский .....	38
SK Slovenský .....	40
SV Svenska .....	42
TR Türkçe .....	44
ZH 中文 .....	46

## pH 3310

### Overview

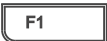
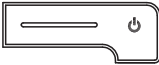
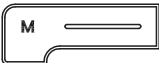
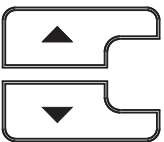
The pH 3310 meter enables you to perform measurements (pH) quickly and reliably.

The pH 3310 meter provides a maximum degree of operating comfort, reliability and measuring certainty.

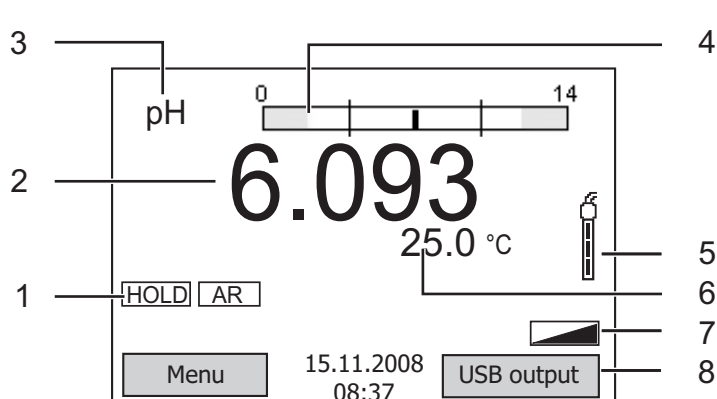
### Keypad

Keys are indicated by brackets <..>.

The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec). A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER\_\_>).

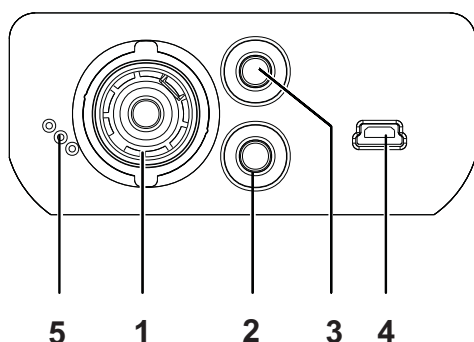
	<F1>: <F1__>:	Situation dependent function (is shown on the display) e.g.: <F1>/[Menu]: Opens the menu for measurement settings <F1__>/[Menu]: Opens the menu for system settings
	<F2>: <F2__>:	
	<On/Off>:	Switches the meter on or off
	<M>:	Selects the measured parameter
	<CAL>: <CAL__>:	Calls up the calibration procedure Displays the calibration data
	<STO>: <STO__> :	Saves a measured value manually Opens the menu for the automatic save function
	<RCL>: <RCL__>:	Displays the manually stored measured values Displays the automatically stored measured values
	<▲><▼>:	Menu control, navigation
	<ENTER>: <ENTER__>:	Opens the menu for measurement settings / confirms entries Opens the menu for system settings
	<AR>:	Freezes the measured value (HOLD function) Switches the AutoRead measurement on or off

## Display

1	Status information
2	Measured value (with unit)
3	Measured parameter
4	Continuous measurement control (CMC function)
5	Sensor symbol (calibration evaluation)
6	Measured temperature (with unit)
7	Further status information
8	Softkeys and date/time

## Jack field /connections

1	pH electrode
2	Reference electrode
3	Temperature sensor
4	USB B interface (device)
5	Service interface

## Initial commissioning

- 1 Insert the supplied batteries. Make sure that the poles are in the correct position.
- 2 Switch on the meter with the <On/Off> key.

### Caution

Follow the safety instructions of the used sensors. The operating manuals of the sensors are available on the supplied CD and on the Internet under [www.WTW.com](http://www.WTW.com).

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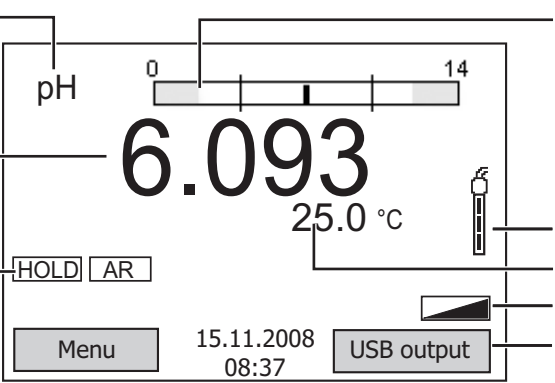
## Visuel

The diagram shows the display screen of the pH 3310 meter. It features a large digital readout of '6.093' with a pH scale above it ranging from 0 to 14. A temperature reading of '25.0 °C' is shown below the main value. At the bottom, there is a date and time '15.11.2008 08:37' and a 'USB output' indicator. On the left side, there are 'HOLD' and 'AR' buttons. On the right side, there is a small icon representing the electrode. The diagram is annotated with numbers 1 through 8 pointing to specific elements.

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# Xylem |'zīləm|

- 1) The tissue in plants that brings water upward from the roots;
- 2) a leading global water technology company.

We're a global team unified in a common purpose: creating advanced technology solutions to the world's water challenges. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. Our products and services move, treat, analyze, monitor and return water to the environment, in public utility, industrial, residential and commercial building services settings. Xylem also provides a leading portfolio of smart metering, network technologies and advanced analytics solutions for water, electric and gas utilities. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise with a strong focus on developing comprehensive, sustainable solutions.

**For more information on how Xylem can help you, go to [www.xylem.com](http://www.xylem.com).**

## **Service and Returns:**

Xylem Analytics Germany  
Sales GmbH & Co. KG  
WTW  
Am Achalaich 11  
82362 Weilheim  
Germany

Tel.: +49 881 183-325  
Fax: +49 881 183-414  
E-Mail [wtw.rma@xylem.com](mailto:wtw.rma@xylem.com)  
Internet: [www.xylemanalytics.com](http://www.xylemanalytics.com)

Xylem Analytics Germany GmbH  
Am Achalaich 11  
82362 Weilheim  
Germany