QUICK START GUIDE

ba77175y01 02/2016

inoLab[®] Multi 9620 IDS

a xylem brand

Contents

EN English	4
DE Deutsch.....	6
CS Česky	8
DA Dansk	10
EL Ελληνικά	12
ES Español	14
FI Suomi	16
FR Français	18
HU Magyar	20
IT Italiano	22
JA 日本語	24
KO 한국어.....	26
NL Nederlands	28
NO Norsk	30
PL Polski.....	32
PT Português	34
RO Română	36
RU Русский	38
SK Slovenský	40
SV Svenska	42
TR Türkçe	44
ZH 中文.....	46

inoLab® Multi 9620 IDS

Overview

The Multi 9620 IDS meter enables you to perform measurements (pH, U, ION, conductivity, D.O.) quickly and reliably.

The Multi 9620 IDS meter provides a maximum degree of operating comfort, reliability and measuring certainty.


Keypad

Keys are indicated by brackets <..>.

The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec). A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER__>).


<F1>	Situation dependent function (is shown on the display)
<F1__>	
<F4>	
<F4__>	
<On/Off>	Switches the meter on or off
<M>	Selects the measured parameter
<CAL>	Calls up the calibration procedure
<CAL__>	Displays the calibration data
<STO>	Saves a measured value manually
<STO__>	Opens the menu for the automatic save function
<RCL>	Displays the manually stored measured values
<RCL__>	Displays the automatically stored measured values
<▲><▼>	Menu control, navigation
<◀><▶>	
<ENTER>	Opens the menu for measurement settings / confirms entries
<ENTER__>	Opens the menu for system settings
<AR>	Freezes the measured value (HOLD function) Switches the AutoRead measurement on or off
<ESC>	Cancels an action
<PRT>	Sends data to the interface (USB)
<PRT__>	Opens the menu for automatic output to USB

Display (example: pH measurement)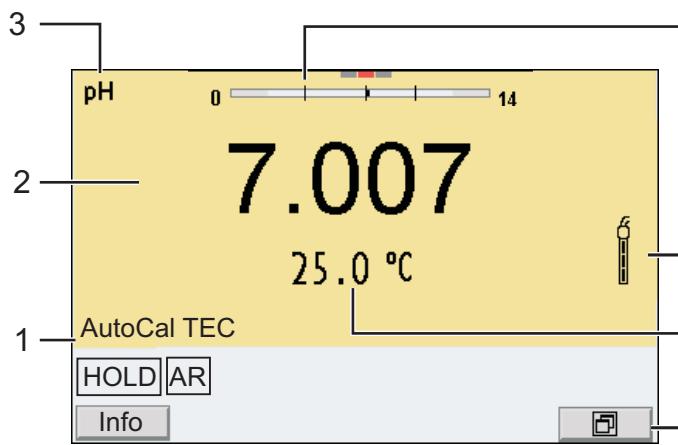

- | | | |
|---|---|---|
| 4 | 1 | Status information |
| 2 | 2 | Measured value (with unit) |
| 3 | 3 | Measured parameter |
| 4 | 4 | Continuous measurement control (CMC function) |
| 5 | 5 | Sensor symbol (calibration evaluation) |
| 6 | 6 | Measured temperature (with unit) |
| 7 | 7 | Softkeys and date/time |

Jack field /connections

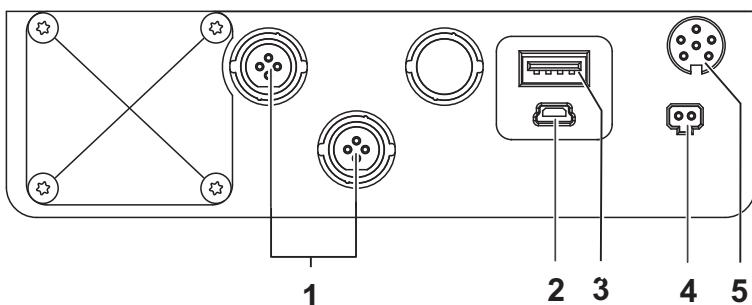- | | | |
|---|---|---|
| 1 | 1 | IDS sensors (pH, U, conductivity, D.O.) |
| 2 | 2 | USB B interface (device) |
| 3 | 3 | USB A interface (host) |
| 4 | 4 | Connection for line adaptor |
| 5 | 5 | Service interface |

Initial commissioning

- | | | |
|---|---|--|
| 1 | 1 | Connect the power pack. |
| 2 | 2 | Switch on the meter with the <On/Off> key. |

Caution

Follow the safety instructions of the used sensors.

The operating manuals of the sensors are available on the supplied CD and on the Internet under www.WTW.com.

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Xylem |'ziləm|

- 1) The tissue in plants that brings water upward from the roots;
- 2) a leading global water technology company.

We're a global team unified in a common purpose: creating advanced technology solutions to the world's water challenges. Developing new technologies that will improve the way water is used, conserved, and reused in the future is central to our work. Our products and services move, treat, analyze, monitor and return water to the environment, in public utility, industrial, residential and commercial building services settings. Xylem also provides a leading portfolio of smart metering, network technologies and advanced analytics solutions for water, electric and gas utilities. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise with a strong focus on developing comprehensive, sustainable solutions.

For more information on how Xylem can help you, go to www.xylem.com.

Service and Returns:

Xylem Analytics Germany

Sales GmbH & Co. KG

WTW

Am Achalaich 11

82362 Weilheim

Germany

Tel.: +49 881 183-325

Fax: +49 881 183-414

E-Mail wtw.rma@xylem.com

Internet: www.xylemanalytics.com

Xylem Analytics Germany GmbH
Am Achalaich 11
82362 Weilheim
Germany