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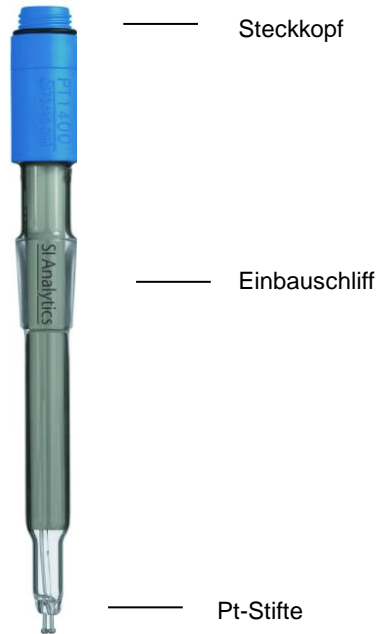
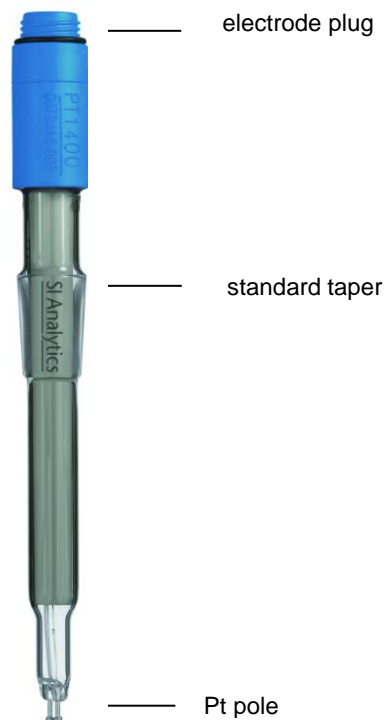
**SI Analytics**  
a xylem brand

### Preparation and general

The double platinum electrodes are delivered ready-to-use. They can be used for dead stop titrations over the whole pH range.

### Storage and maintenance

The electrode should be stored at a temperature between 0 and 40°C.

### Cleaning

Dirt and contamination at the Pt sensor will lead to measurement inaccuracies. They can be removed with the following procedure:

- **Coatings and deposits** with diluted mineral acids (e.g. diss. hydrochloric acid)
- **Organic contaminations** with the aid of suitable solvents
- **Grease** with tenside solution or alcohol
- **Proteins** with hydrochloric pepsin solution (cleaning solution L 510)
- **Persistent films** by carefully rubbing down with fine sandpaper or a diamond file. **Do not bend the Pt pole!**

After the procedure the electrode must be rinsed with distilled water.

### Quality

Every electrode must meet the strict quality requirements of final testing. The working life depends mainly on the usage conditions. Extreme conditions include for example high or frequently fluctuating temperatures, strong acids and caustic solutions. Hydrofluoric acid, sodium hydroxide and hot phosphoric acid corrode glass.

### Technical details

	KF1100	Pt 1200	Pt 1400	KF 1150
Temperature range	-----	-30 ...+135°C	-----	-----
Length (mm)	96	120	103	116
Sensor	-----	2 pole, 1mm ø	-----	-----
Standard taper	NS 7.5	no	NS 14.5	NS 7.5
Connection	fixed cable with 2X4mm plug	plug head (cable e.g. L1NN)	plug head (cable e.g. L1NN)	fixed cable with 2X4mm plug

### Further information

Further information can be found in our product catalogs.

Subject to technical amendment.

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### Xylem Analytics Germany GmbH