**BEDIENUNGSANLEITUNG**  
**OPERATING MANUAL**  
**MODE D'EMPLOI**  
**INSTRUCCIONES DE OPERACIÓN**

ba15350defs03 03/2015

**Ag/S 500 DIN, Ag/S 500 BNC**  
**Ag/S 800 DIN, Ag/S 800 BNC**

a xylem brand

---

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---

# Ag/S 500 DIN, Ag/S 500 BNC Ag/S 800 DIN, Ag/S 800 BNC

SILVER ION AND SULFIDE SENSITIVE ELECTRODE

---

a xylem brand

**Note**

The latest version of the present operating manual can be found on the Internet under [www.WTW.com](http://www.WTW.com). It is also available in other languages.

**Copyright**


© 2014, WTW GmbH

Reprinting - even as excerpts - is only allowed with the explicit written authorization of WTW GmbH.

Printed in Germany.

## Overview

View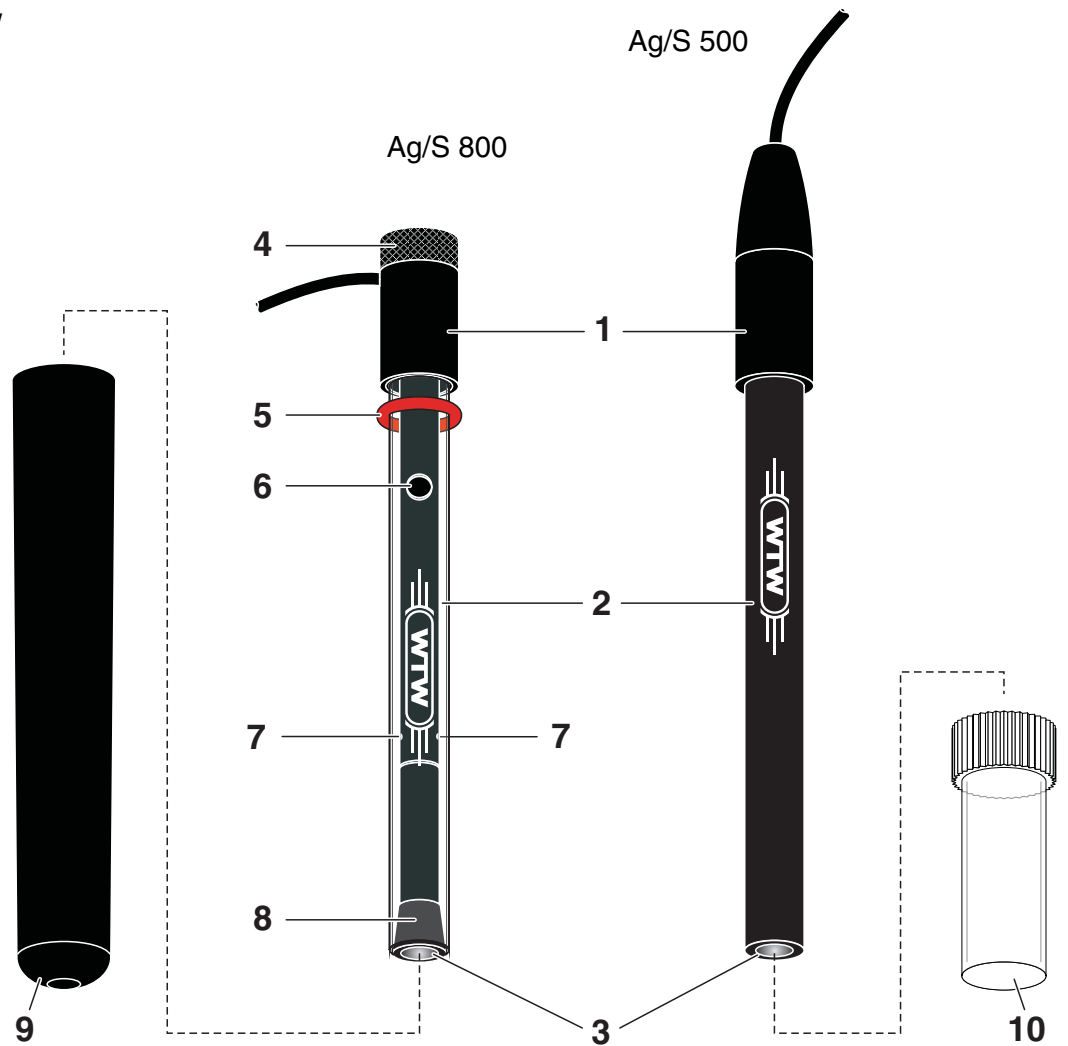

1	Connection head with connection cable
2	Shaft
3	Membrane
4	Knurled nut
5	O-ring
6	Filling opening for bridge electrolyte
7	Inner junctions
8	Ground junction
9	Storage vessel for Ag/S 800
10	Protective cap for Ag/S 500

## Commissioning

### Combination electrode Ag/S 800

1	Remove the storage vessel.
2	Fill the bridge electrolyte ELY/BR/503 into the filling opening.
3	Shortly press the shaft of the combination electrode against the connection head to wet the ground junction with bridge electrolyte.
4	When doing so refill any spilled bridge electrolyte so that the inner junctions are covered with bridge electrolyte.
5	Rinse the combination electrode with deionized water.
6	Wipe the shaft using a clean paper towel.


### Double rod electrode Ag/S 500 + R 503

For measurements with the Ag/S 500 electrode, a reference electrode is required (e.g. R 503). The two electrodes together form a double rod combination electrode.

1	Put the reference electrode into operation (see operating manual of the reference electrode).
2	Remove the protection cap of the electrode.
3	Rinse the electrode with deionized water.
4	Wipe the shaft using a clean paper towel.

## Conditioning, calibration, measurement

### General information

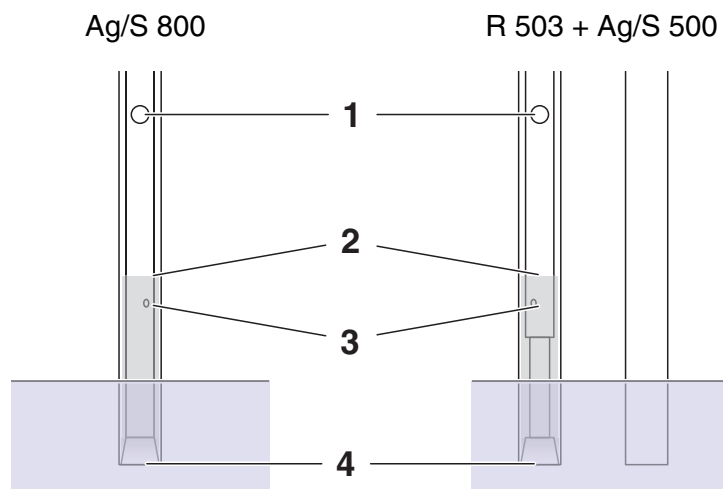When operating the electrode ensure that

- the filling opening (1) for the bridge electrolyte is open
- the inner junctions (3) are covered with bridge electrolyte
- no air bubbles are in the bridge electrolyte
- the depth of immersion is within the optimum range:

Minimum depth of immersion

The ground junction (4) must be covered

Maximum depth of immersion

Approx. 1 cm below the fluid level (2) of the bridge electrolyte

### Before measuring

1	Condition every ion sensitive electrode in standard solution before using it. To achieve a good responding behavior, the concentration of the standard solution should be similar to the concentration expected of the test sample.
2	Remove any air bubbles in the bridge electrolyte by slightly knocking against the shaft.
3	Calibrate according to the operating manual of the meter and the analysis specification.

The conditioning time depends on the previous history of the ion sensitive electrode. Connect the electrode(s) to the meter and switch to the measured value display (see operating manual of the meter). The ion sensitive electrode is sufficiently conditioned when the measured value no longer changes considerably under otherwise constant conditions (temperature, composition of the solution).

## Sample preparation

### Ag<sup>+</sup> measurement

Add 2 % ISA/FK solution. This sample conditioning solution creates optimum conditions for measuring. It provides a constant ionic strength and similar diffusion potentials at the reference electrode in standard solution and test sample.

### S<sup>2-</sup> measurement


Add 50 % SAOB II solution.

Preparation of the SAOB II solution:

- 200 ml/l 10 molar NaOH
- 35 g/l ascorbic acid
- 67 g/l EDTA

A freshly prepared SAOB II solution (Sulfide Anti-Oxidant Buffer) is colorless or slightly yellowish brown. If the solution has already strongly oxidized its color is dark brown. The solution can no longer be used.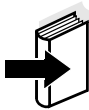

If you would like to have more detailed information concerning sample preparation and measuring procedures, WTW provides a large number of application reports for various applications.

## Response times

The response time depends on the concentration range. It is

- several seconds at high concentrations,
- several minutes near the detection limit.

The measured value is stable if the value does not change by more than 0.1 mV within 30 seconds.

## Interferences

- Mercury
- Protein

## Aging

Please note that every electrode undergoes a natural aging process. The response time increases and the slope decreases with the age of the electrode. The following factors shorten the lifetime considerably:

- Unsuitable storing conditions (see section STORAGE)
- Special measuring conditions (e.g. organic solutions, frequent measuring with mercury or protein concentrations)
- High temperatures
- High changes in temperature

The warranty does not cover damage caused by measuring conditions and mechanical damage.

## Maintenance

- Combination electrodes:  
Refill any used up bridge electrolyte.
- In the case of increased response times of the electrode perform the following activities:
  - Polish the membrane surface using an SF/K polishing strip (see section RECOMMENDED ACCESSORIES). To do so, moisten a polishing strip with deionized water, and with circular movements polish the membrane surface for approx. 30 seconds.
  - Then recondition the electrode (see section CONDITIONING, CALIBRATION, MEASUREMENT).
- Clean the inside of the combination electrode. To do so, open the combination electrode as follows:

### Opening the combination electrode

1	Unscrew the knurled nut from the connection head.
2	Position the connection cable upward in a straight line with the connection head.
3	Unscrew the connection head.
4	Push the connection head and pressure spring over the connection cable.
5	Push the inner body out of the shaft.

### **NOTE**

**Never pull the connection cable of the electrode.  
The cable might be damaged.**

## Storage

### Between two measurements

Condition the electrode in standard solution.

### Overnight

**Ag/S 500:** Condition the electrode in standard solution.

**Ag/S 800:** Condition the combination electrode in standard solution with the filling opening open. To avoid a contamination of the bridge electrolyte with standard solution, the level of the bridge electrolyte must be clearly above the level of the conditioning solution.

For the following measurement, exchange the bridge electrolyte. To do so, press the shaft of the combination electrode against the connection head and let the bridge electrolyte drain off. Then fill new bridge electrolyte into the filling opening.

### Longer than overnight

**Ag/S 500:** Rinse the electrode with deionized water, then dab it dry with a clean paper towel and put on the protective cap. Store the electrode in a dry place.

**Ag/S 800:** Remove the bridge electrolyte, then rinse the combination electrode with deionized water and dab it dry with a clean paper towel. Store the combination electrode in the storage vessel in a dry place. For the following measurement, prepare the combination electrode as described in section COMMISSIONING.

Store the reference electrode according to the instructions in its operating manual.

## Recommended accessories

Description	Model	Order no.
Reference electrode for electrode Ag/S 500	R 503/P*	106570
	R 503/D**	106571
Bridge electrolyte	ELY/BR/503	106 575
ISA sample conditioning solution	ISA/FK	140 110
Polishing strips (24 pieces)	S/FK	180130

\* Pin plug


\*\* Banana plug

### Calibration line of a silver ion / sulfide sensitive electrode


**Ag<sup>+</sup> measurement**

Electromotive force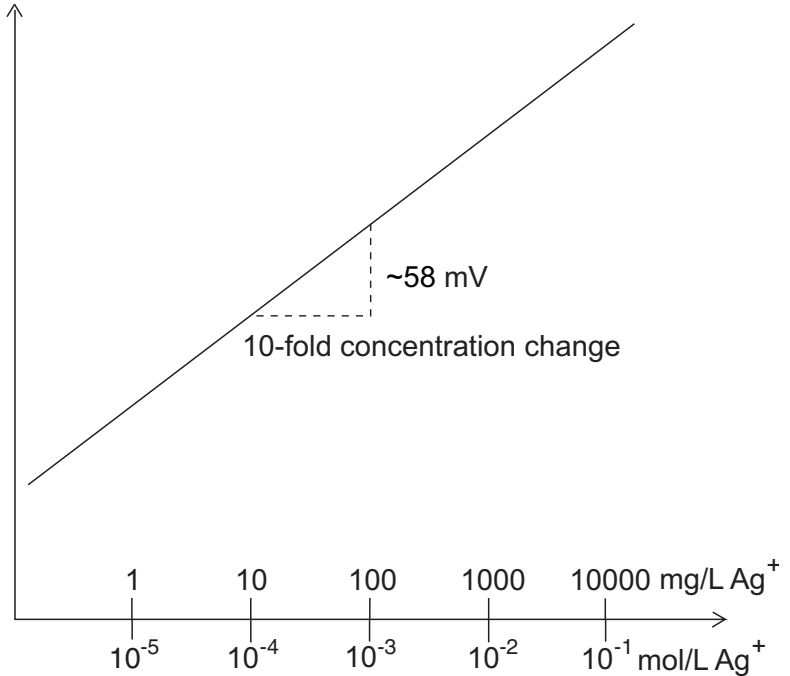

**S<sup>2-</sup> measurement**

Electromotive force

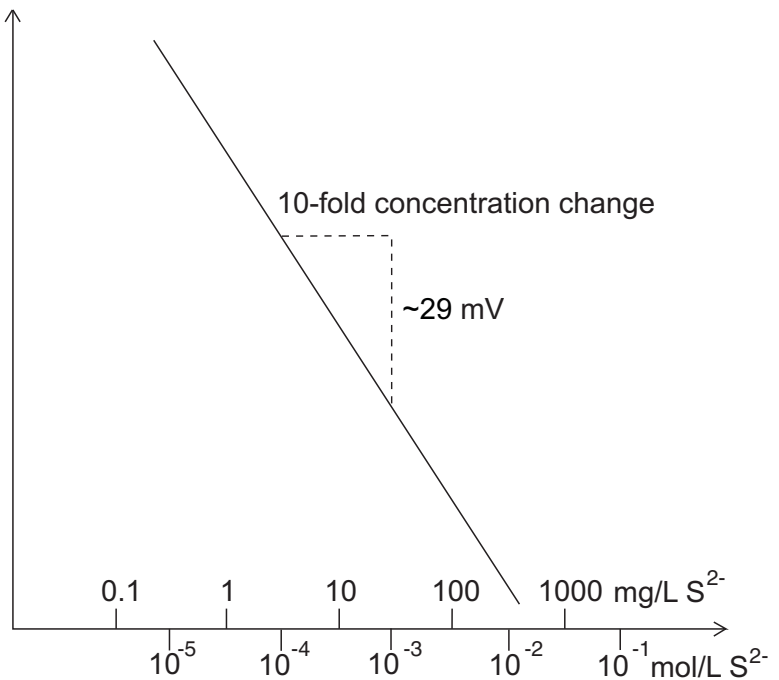## What to do if ...

Measured value unstable	Cause	Remedy
	– Inner junctions not sufficiently wetted with bridge electrolyte (Ag/S 800)	– Fill up bridge electrolyte until the inner junctions are covered with bridge electrolyte
– Inner junctions encrusted (Ag/S 800)	– Leave the bridge electrolyte to react on the inner junctions for some hours until the crusts have dissolved.	
– Ground junction contaminated (Ag/S 800)	– Rinse ground junction with bridge electrolyte	
– Cable broken	– Exchange the electrode	

Slope too low	Cause	Remedy
	– Membrane surface contaminated	– (see MAINTENANCE)
– Conditioning time too short	– Extend the conditioning time	
– Standard solutions too old	– Use new standard solutions	
– Inner junctions encrusted (Ag/S 800)	– Leave the bridge electrolyte to react on the inner junctions for some hours until the crusts have dissolved.	
– Electrode defective	– Exchange the electrode	

## Technical data

<b>Measuring ranges</b>	Ag <sup>+</sup>	0.01 ... 108,000 mg/l	(10 <sup>-7</sup> ... 1 mol/l Ag <sup>+</sup> )
	S <sup>2-</sup>	0.003 ... 32,000 mg/l	(10 <sup>-7</sup> ... 1 mol/l S <sup>2-</sup> )
<b>Reproducibility</b>	± 2 % for silver; ± 4 % for sulfide		
<b>pH range</b>	2 ... 12	(silver determination: pH < 8; sulfide determination with SAOB II solution)	
<b>Temperature range</b>	0 ... 80 °C	(shortly up to 100 °C)	
<b>Resistance</b>	< 1 MΩ		
<b>Length</b>	Ag/S 500:	170 mm (including 50 mm connection head)	
	Ag/S 800:	153 mm (including 33 mm connection head)	
<b>Diameter</b>	Shaft:	12 mm	
	Connection head:	16 mm	
<b>Cable length</b>	1 m		
<b>Plug</b>	DIN plug or BNC plug, depending on the design.		

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# What can Xylem do for you?

We're a global team unified in a common purpose: creating innovative solutions to meet our world's water needs. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. We move, treat, analyze, and return water to the environment, and we help people use water efficiently, in their homes, buildings, factories and farms. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise, backed by a legacy of innovation.

**For more information on how Xylem can help you, go to [xylem.com](http://xylem.com).**

**Service address:**

Xylem Analytics Germany  
Sales GmbH & Co. KG  
WTW  
Dr.-Karl-Slevogt-Str. 1  
82362 Weilheim  
Germany

Tel.: +49 881 183-325  
Fax: +49 881 183-414  
E-Mail [wtw.rma@xylem.com](mailto:wtw.rma@xylem.com)  
Internet: [www.WTW.com](http://www.WTW.com)

Xylem Analytics Germany GmbH  
Dr.-Karl-Slevogt-Str. 1  
82362 Weilheim  
Germany