OPERATING MANUAL

ba75739e02 07/2013

SenTix[®] H
SenTix[®] HW
SenTix[®] HWD

SenTix®

PH COMBIN. ELECTRODES WITH LIQUID REFERENCE SYSTEM AND GROUND JUNCTION





For the most recent version of the manual, please visit www.WTW.com.

Copyright

© Weilheim 2013, WTW GmbH

Reproduction in whole - or even in part - is prohibited without the express written permission of WTW GmbH, Weilheim.

Printed in Germany.

SenTix[®] Technical data

Technical data

General data

| WTW model | Reference electro- lyte | Junction | NTC | Special features |
|-------------------------|-----------------------------------|----------|-----|----------------------------|
| SenTix [®] H | KCI 3 mol/L, Ag ⁺ free | Ground | No | Adjustable ground junction |
| SenTix [®] HW | KCI 3 mol/L, Ag ⁺ free | Ground | No | Adjustable ground junction |
| SenTix [®] HWD | KCI 3 mol/L, Ag ⁺ free | Ground | Yes | |

Measurement and application characteristics

| WTW model | pH measur- ing range | Allowed temperature range | Membrane resistance at 25 °C | Typical application |
|-------------------------|-------------------------|---------------------------|------------------------------|--|
| SenTix [®] H | 0 14 | 0 80 °C | < 2 GOhm | Highly alkaline and highly saltine solutions |
| SenTix [®] HW | 0 14 | 0 60 °C | < 800 MOhm | Solutions with low ion |
| SenTix [®] HWD | 0 14 | -5 100 °C | < 600 MOhm | contents, emulsions, suspensions |

Shaft dimensions, shaft material, electrical connection

| WTW model | Shaft | | | Electrical connection | | |
|-------------------------|----------------|--------|----------|----------------------------------|------------------------|-----------------|
| | Length [mm] | Ø [mm] | Material | Combination electrode connection | Meter connection | Cable length |
| SenTix [®] H | 170 | 12 | Glass | S7 plug-in connector | depending on S7 cable* | |
| SenTix [®] HW | 170 | 12 | Glass | S7 plug-in connector | depending on S7 cable* | |
| SenTix [®] HWD | 170 | 12 | Glass | Fixed cable | DIN** + banana | 1 m |

 $^{^{\}star}$ Connection cable not included in the scope of delivery of the combination electrode (see Wear parts and accessories)

ba75739e02 07/2013 **3**

^{**} Coaxial plug according to DIN 19262

Technical data SenTix[®]


Commissioning, measuring, calibration

Commissioning

The combination electrode is filled with reference electrolyte solution in the factory. Prepare the combination electrode for measuring as follows:

Open the refilling opening for the reference electrolyte solution. Depending on the model, the stopper of the refilling opening is an elastomer stopper or a slider.

The refilling opening must always be open during calibration and measurement!

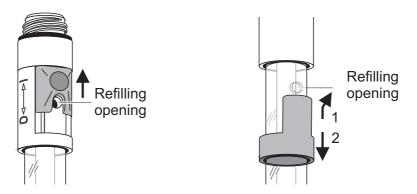

 Remove the watering cap from the electrode tip. Possible salt deposits in the area of the watering cap do not affect the measuring characteristics and can easily be removed with deionized water.



Note

Please keep the watering cap. It is required for the combination electrode to be stored. Always keep the watering cap clean.

- Connect the combination electrode to the meter
- Calibrate the combination electrode according to the operating manual of the meter. When doing so, observe the general rules on the following page.

Ground junction of the SenTix HWD

When the SenTix HWD combination electrode is delivered, the ground sleeve is fixed with a shrinkable tubing. For commissioning, carefully remove the shrinkable tubing with a sharp knife or scalpel.

Caution:

The ground sleeve is made of synthetic material and can easily be damaged!


To moisten the ground surfaces hold the combination electrode in a vertical position with the refill opening open and carefully loosen the ground junction by turning it slightly until reference electrolyte solution comes out of the junction. Then put the ground sleeve on again while turning it slightly.

SenTix[®] Technical data

Ground junction of SenTix H and SenTix HW

With the ground junction of the SenTix H and SenTix HW combination electrodes, the electrode outflow can be adjusted with a thread:

 Hold the combination electrode with the pH glass membrane pointing downward. Using the provided special wrench, open the ground junction until some electrolyte solution can be seen to come out of the junction and all ground surfaces are completely moist.

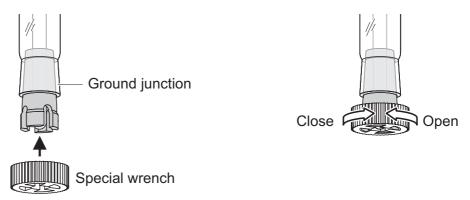

- Carefully and using little power close the junction until the contact between the rough ground surfaces can be noticed as a resistance. The combination electrode is now ready to measure.
 - Caution: Always use very little power to close the junction so the ground surfaces are not damaged!
- The ground junction can be opened further so that more electrolyte solution comes out.

Calibration and measurement: General rules

- Make sure the refilling opening for the reference electrolyte solution is open.
- Avoid the displacement of any solution (sample or buffer solution) from one measurement to the next by taking the following measures:
 - Shortly rinse the calibration and sample beakers with the solution the beakers are to be filled with next.
 - Between measurements, rinse the combination electrode with the solution that follows. Alternatively, you can also rinse the combination electrode with deionized water and then carefully dab it dry.
- Immerse the combination electrode in the solution in a vertical or slightly tilted position.
- Make sure the immersion depth is correct. The ground junction must be completely submersed in the solution. At the same time, the level of the reference electrolyte must be at least 2 cm above the level of the solution.
- Provide approximately the same stirring conditions for measuring as for calibrating.

ba75739e02 07/2013 **5**

Storage SenTix[®]

Subsequent calibrations

The frequency of subsequent calibrations depends on the application. Many meters provide an option where you can enter a calibration interval. After the calibration interval has expired, the meter will automatically remind you of the due calibration.

Storage

During short measuring breaks

With the refilling opening open, immerse the combination electrode in reference electrolyte (KCI 3 mol/L, Ag⁺ free). Observe the minimum depth of immersion. Prior to the next measurement, shortly rinse the combination electrode with the test sample or deionized water.



Note

Prevent contact of the pH membrane to the beaker bottom to avoid scratches on the pH membrane.

Overnight or longer

Put the clean combination electrode in the watering cap that is filled with reference electrolyte (KCl 3 mol/L, Ag⁺ free) and close the refilling opening.



Note

pH combination electrodes must not be stored dry or in deionized water. The combination electrode could be permanently damaged by this. If the liquid in the watering cap has dried up, condition the combination electrode in reference electrolyte (KCI 3 mol/L, Ag+ + free) for at least 24 hours.



Note

During longer storing periods, salt sediments may develop on the watering cap. They do not affect the measuring characteristics and can easily be removed with deionized water when the combination electrode is put into operation again.

Aging

Every pH combination electrode undergoes a natural aging process. With aging, the responding behavior becomes slower and slower and the electrode slope and asymmetry change. Moreover, extreme operating conditions can considerably shorten the lifetime of the combination electrode. These are:

- Strong acids or lyes, hydrofluoric acid, organic solvents, oils, fats, bromides, sulfides, iodides, proteins
- High temperatures
- High changes in pH and temperature.

The warranty does not cover failure caused by measuring conditions and mechanical damage.

Maintenance and cleaning

Refilling the reference electrolyte

During operation, a small amount of reference electrolyte leaks through the junction from the combination electrode into the test sample. If the level of reference electrolyte becomes too low with time, refill it through the refilling opening. Refilling is very easy with the enclosed dropping bottle. Proceed as follows:

- Cut off the tip of the dropping bottle at a right angle until the opening in the tip can be seen
- Open the refilling opening of the combination electrode
- Press the tip of the dropping bottle into the refilling opening while turning it slightly
- Pump the reference electrolyte in the shaft using the dropping bottle
- Pull the dropping bottle out of the refilling opening while turning it slightly if necessary.



Cleaning

Remove water-soluble contamination by rinsing with deionized water. Remove other contamination as follows:

| Contamination | Cleaning procedure |
|-----------------------------|--|
| Fat and oil | Rinse with water containing household washing-up liquid |
| Lime and hydroxide deposits | Rinse with citric acid (10 % by weight) |
| Proteins | Immerse in pepsin cleaning solution PEP/pH for approx. 1 hour. Note: Make sure the level of the reference electrolyte is above that of the cleaning solution. |



Note

Hydrofluoric acid, hot phosphoric acid and strong alkaline solutions destroy the glass membrane.

ba75739e02 07/2013 **7**

If you want to clean and rinse the ground junction, you can open it as follows:

- SenTix H and SenTix HW: completely unscrew the ground stopper (see figure on Seite 5).
- SenTix HWD: slide the ground sleeve from the ground stopper towards the shaft while turning it slightly.

After cleaning

Rinse the combination electrode with deionized water and condition it in reference electrolyte solution for at least 1 hour. Then recalibrate the combination electrode.

Wear parts and accessories

| Description | Model | Order no. |
|---|----------|-----------|
| Reference electrolyte solution 250 mL (KCl 3 mol/L, Ag ⁺ free) | KCI-250 | 109 705 |
| Pepsin cleaning solution 3 x 250 ml | PEP/pH | 109 648 |
| Connection cable S7 plug-in connector/DIN, 1 m | AS/DIN | 108 110 |
| Connection cable S7 plug-in connector/DIN, 3 m | AS/DIN-3 | 108 112 |
| Connection cable S7 plug-in connector/BNC, 1 m | AS/BNC | 108 114 |



Note

Detailed information on our wide range of buffer solutions and more accessories is given in the price list of the WTW catalog "Laboratory and field instrumentation".

Xylem | zīləm

- 1) The tissue in plants that brings water upward from the roots;
- 2) a leading global water technology company.

We're 12,500 people unified in a common purpose: creating innovative solutions to meet our world's water needs. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. We move, treat, analyze, and return water to the environment, and we help people use water efficiently, in their homes, buildings, factories and farms. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise, backed by a legacy of innovation.

For more information on how Xylem can help you, go to www.xyleminc.com



Wissenschaftlich-Technische Werkstätten GmbH

Dr.-Karl-Slevogt-Straße 1 D-82362 Weilheim Germany

Tel: +49 881 183-0 Fax: +49 881 183-420 E-Mail: info@wtw.com Internet: www.wtw.com